

Installing the Analyst Administrator Console 3.0

Topics in this section:

- [Important Information to Know Before Installing on page 1](#)
- [Contents of the Analyst Administrator Console DVD on page 2](#)
- [Upgrading the Administrator Console Server on page 3](#)
- [Installing the Administrator Console Server and Client on page 3](#)
- [Installing the MultiQuant Software Update on the Administrator Console Server on page 4](#)

Important Information to Know Before Installing

To ensure your installation is successful, read the following points carefully before you start any of the procedures in this document.

- The person installing the Administrator Console server must have Administrator rights on the computer.
- The Administrator Console server cannot be installed on the same workstation as the Analyst[®] software.
- If you have a firewall running on the computer running the Administrator Console server, the following three ports, 1633(tcp), 1634(tcp), and 6001(tcp) must be opened on both the client and server computers in order for the Administrator Console to function correctly.

If you already have the Administrator Console 1.0 or 2.0 installed and you want to upgrade the server, then see [Upgrading the Administrator Console Server on page 3](#).

Operating System Requirements

The following are the operating system requirements for the Administrator Console server:

- Microsoft Windows XP SP2
- Microsoft Windows XP SP3
- Windows 7 SP1 (only tested on SP1)

Network Environment Requirements

The Analyst 1.6 software is supported in the following network environment:

- Microsoft Windows Server 2003 and 2008 (for Active Directory networks).

Compatible Software

- Analyst software version 1.4.1 and corresponding software components and HotFix releases.
- Analyst software version 1.4.2 and corresponding HotFix releases.
- Analyst software version 1.5 and corresponding software components and HotFix releases.
- Analyst software version 1.5.1 and corresponding HotFix releases.
- Analyst software version 1.5.2 and corresponding HotFix releases.
- Analyst software version 1.6.
- The Administrator Console server is compatible with the Administrator Console client versions 1.0, 2.0, and 3.0.
- MultiQuant™ software versions 2.0.x and 2.1

Minimum Computer Requirements

For optimal performance, we recommend a computer with a minimum 1.7 GHz processor.

Contents of the Analyst Administrator Console DVD

The following software applications are included on the Analyst Administrator Console software DVD:

- AdminConsole.exe: Installs the Administrator Console client only
- ACServer.exe: Installs both the Administrator Console client and server
- AAC Release Notes.pdf
- Analyst Administrator Console Installation Instructions.pdf

Administrator Console Components

The Administrator Console is composed of the following:

- Administrator Console server
- Administrator Console client

Administrator Console Server

The Administrator Console server is installed on a computer from the Administrator Console installation DVD. The Administrator Console client and the FastObjects server are automatically installed during server installation.

All security information is stored in the database on the server. During installation the security database is automatically populated with the default Analyst software roles and users.

Additionally, each time the Analyst software is started, the backup copy of the database on each registered workstation is updated to reflect the master copy on the Administrator Console server.

Administrator Console Client

The Administrator Console client is a Microsoft Management Console (MMC) snap-in and is installed on a separate computer. When the Administrator Console client is installed on a workstation, the console administrators can use it to access the server remotely.



Note: The Administrator Console client is also automatically installed on a workstation as part of the Analyst software installation.

Upgrading the Administrator Console Server

(Optional) If you are not sure of the version installed, click the **Click Here for support information** link in the Add/Remove Programs. The previously released version is 2.0.

1. Close all Administrator Console clients, local and remote.
2. Close the Administrator Console client on the Administrator Console Server.
3. Back up the security database; make a copy of the file C:\Program Files\Administrator Console Server\SecurityDB.odc
4. Open the Control Panel and in the Add/Remove programs uninstall the Administrator Console Server.



Note: If you want to keep the existing database, then during installation do not overwrite the security database. Select No when you receive the message “Old security database is found. Would you like to overwrite it?” on the installation wizard page. This message appears after you have accepted the licence agreement.

5. Install the Administrator Console Server 3.0. See [Installing the Administrator Console Server and Client on page 3](#).
6. Restart the computer.
7. Open the Administrator Console client on the Administrator Console Server computer. Initial startup will take several minutes to upgrade the database to the new format. The upgrade will not change the contents of the database.
8. Start and connect the Administrator Console clients.

Installing the Administrator Console Server and Client

If you already have the Administrator Console 3.0 installed and you want to upgrade the server, then see [Upgrading the Administrator Console Server on page 3](#).

The Administrator Console server can be installed on any computer that does not have the Analyst software installed.

Installing the Administrator Console Server

1. Insert the DVD, navigate to the DVD drive, and then double-click **ACServer.exe**.
The Welcome page appears.
2. Click **Next**.
The License Agreement page appears.
3. Read the agreement, and then click **I accept the terms in the license agreement**.
4. Click **Next**.
The Ready to Install the Program page appears
5. Click **Install**.
The InstallShield Wizard Completed page appears.
6. Click **Finish**.
The Administrator Console Server Installer Information dialog box appears.
7. Click **Yes** to restart the computer.
8. Remove the DVD from the drive.
The Administrator Console server files are installed in the Program Files\Administrator Console Server folder. The FastObjects server runs as a service and it is found in Administrative Tools, Services.

Installing the MultiQuant Software Update on the Administrator Console Server

If your organization is using the MultiQuant software with Audit Trail and Security Features, then you must update the AAC Server to add the MultiQuant software security items. The MultiQuant software update file (AACUpdateSetup.exe) is located on the MultiQuant software installation DVD in the InstallerAdminConsole folder.

- Double-click the AACUpdateSetup.exe and then follow the on-screen instructions.

Installing the Administrator Console Client

1. Insert the DVD, navigate to the DVD drive and then double-click **AdminConsole.exe**.
The Welcome page appears.
2. Click **Next**.
The Ready to Install the Program page appears.
3. Click **Install**.
The Setup Complete page appears.
4. Click **Finish**.
5. Remove the DVD from the drive.
The Administrator Console client files are installed in the Program Files\Administrator Console folder.

This document is provided to customers who have purchased AB SCIEX equipment to use in the operation of such AB SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as AB SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such usage is intended only to designate those manufacturers' products as supplied by AB SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

AB SCIEX makes no warranties or representations as to the fitness of this equipment for any particular purpose and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use to which the purchaser may put the equipment described herein, or for any adverse circumstances arising therefrom.

For research use only. Not for use in diagnostic procedures.

The trademarks mentioned herein are the property of AB Sciex Pte. Ltd. or their respective owners.
AB SCIEX™ is being used under license.



AB SCIEX

71 Four Valley Dr., Concord, Ontario, Canada. L4K 4V8.

AB SCIEX LP is ISO 9001 registered.

© 2011 AB SCIEX.

Printed in Canada.
