

# Forensic Method

## Forensic drug screening analysis

### Elevate your forensic testing with the X500R QTOF System

Method details and access to HR-MS/MS libraries to screen for forensic drugs in blood extracts using HPLC coupled with the X500R QTOF system, powered by SCIEX OS Software.



### Sample Prep

- Step 1** • Add 0.2ml of blood into centrifuge tube
- Step 2** • Add 0.8mL of cold acetonitrile dropwise while vortexing
- Step 3** • Centrifuge for 10 minutes at 5000rpm
- Step 4** • Transfer supernatant to clean glass vials and dry down.
- Step 5** • Reconstitute in 200µL of 20% methanol in water.
- Step 6** • Vortex for 1 minute and centrifuge for 10 minutes at 10,000xg
- Step 7** • Transfer supernatant to clean glass HPLC vials

### LC Method

<i>Column</i>	Phenomenex Kinetex Biphenyl, 100 x 3.0 mm, 2.6 um column	
<i>Mobile Phase A</i>	0.1% formic acid in water	
<i>Mobile Phase B</i>	0.1% formic acid in methanol	
<i>Flow rate</i>	0.6 mL/min	
<i>Column temperature</i>	30°C	
<i>Injection volume</i>	10 µL	
<i>Gradient profile</i>	<b>Time (min)</b>	<b>% B</b>
	0	2
	1	2
	7	65
	7.1	100
	9	100
	9.1	2
	12	2



SCIEX OS can deliver faster method set-up

# MS Method

**Forensic Screen IDA**

Method duration: 6.5 min | Total scan time: 0.41118 sec  
 Estimated cycles: 948

**Source and Gas Parameters**

Ion source gas 1: 50 psi | Curtain gas: 30 | Temperature: 600 °C  
 Ion source gas 2: 50 psi | CAD gas: 7

**Experiment IDA**

Polarity: Positive V | Spray voltage: 5500 V

**TOF MS**

TOF start mass: 100 Da | Declustering potential: 80 V | Collision energy: 10 V  
 TOF stop mass: 1000 Da | DP spread: 0 V | CE spread: 0 V  
 Accumulation time: 0.1 sec

**IDA Criteria** (Small molecule)

Maximum candidate ions: 10 |  Dynamic background subtraction  
 Intensity threshold exceeds: 10 cps |  Exclude former candidate ions  
 For: 6 sec | After: 1 occurrences

**Advanced Criteria**

**TOF MSMS**

Presursor ion: 6.10 Da | Declustering potential: 80 V | Collision energy: 35 V  
 TOF start mass: 40 Da | DP spread: 0 V | CE spread: 15 V  
 TOF stop mass: 1000 Da | Accumulation time: 0.025 sec

Buttons: Start, Stop, Save...

Suggested IDA (Information Dependent Acquisition) conditions for routine forensic drug screening as displayed in SCIEX OS

## Data Processing

Analytics interface showing a chromatogram with a significant peak at approximately 4.5 minutes. The mass spectrum plot shows relative intensity versus m/z, with a base peak at m/z 200. The interface includes various toolbars and data tables for sample and compound information.

Review your results with utmost efficiency using SCIEX OS for simultaneous quantitation and MS/MS library confirmation.

## X500R High Resolution Libraries

Download the XIC List

Download a FREE Trial of the MS/MS Library

Download a free XIC compound list detailing a full list of forensic drug compounds including molecular formula and accurate mass.

Download a free trial of the forensic high resolution MS/MS library, containing 1703 compounds.

[Learn more at sciex.com/X-Forensics.](http://sciex.com/X-Forensics)

AB Sciex is doing business as SCIEX.

© 2015 AB Sciex. For Research Use Only. Not for use in diagnostic procedures. The trademarks mentioned herein are the property of AB Sciex Pte. Ltd. or their respective owners. AB SCIEX™ is being used under license.

Document number: RUO-MKT-07-2865-A