

A novel fragmentation technology allows for in-depth glycopeptide characterization in glycoproteins

Jenny Albanese and Zoe Zhang
SCIEX, USA

INTRODUCTION

A newly developed electron activated dissociation (EAD) approach allows to tune electron energy to produce different fragmentation patterns of the electron-based dissociation (ExD) family. The data presented here demonstrate the advantage for glycoproteins using this novel fragmentation cell over traditional collision-induced dissociation (CID) used for peptide mapping, which does not allow for consistent identification and localization of glycans on peptides. Combining increased detection of fragments using a mechanism for increased duty cycle on a Q-TOF instrument (the Zeno trap) enables higher confidence in data assignment, making the Zeno EAD combination ideal for in-depth analysis of glycopeptides.

Key features of the SCIEX solution

- EAD allows in combination with fast data-dependent acquisition (DDA) greater in-depth analyses of next generation glycoprotein therapeutics and standard mAbs by offering an alternative fragmentation
- The tunable electron energy provides a higher level of structural information for glycopeptide characterization. It resulted in rich fragmentation with predominantly peptide backbone fragments leaving the glycan intact on the glycosylation site, enabling localization information.
- Increased detection of fragments (5 to 10-fold) was achieved using the Zeno trap. Enhanced sensitivity is required with alternative fragmentation compared to CID, due to the greater number of diagnostic fragments created.
- Reproducible fragmentation with EAD for singly, doubly and multiply charged ions. EAD is capable of fragmenting peptides with a variety of charge states.
- Fully automated DDA using EAD with SCIEX OS software and automated data interpretation with the new Biologics Explorer software from SCIEX offer a streamlined characterization workflow

MATERIALS AND METHODS

Sample preparation:

Therapeutic monoclonal antibodies adalimumab, rituximab as well as an antibody-drug-conjugated sample (ado- trastuzumab emtansine) were denatured with 7.2 M guanidine hydrochloride, 100mM Tris buffer (pH 7.2), followed by reduction with 10mM DL-dithiothreitol and alkylation with 30mM iodoacetamide. Digestion was performed with trypsin/Lys-C enzyme at 37°C for 16 h.

Time [min]	Mobile phase A [%]	Mobile phase B [%]
Initial	98	2
5	98	2
6	90	10
40	55	45
44	10	90
46	10	90
47	98	2
50	98	2
51	10	90
54	10	90
55	98	2
60	98	2

HPLC conditions:

An ExionLC system equipped with a CSH C18 column (1.7 μ m particle size, 130 \AA , 2.1x100 mm, Waters) was used to separate 10 μ L (4 μ g) of the trypsin/Lys-C digest. The mobile phase A consisted of water with 0.1% formic acid, while the organic phase B was acetonitrile 0.1% formic acid. A gradient profile was used at a flow rate of 300 μ L/min. The column temperature was maintained at 50°C.

Parameter	CID	EAD
Collision energy	rolling	12 V
Electron KE	NA	7 eV
Electron beam current	NA	5500 nA
ETC	NA	100
Zeno trap	ON	ON
Accumulation time	0.05 s	0.09 s

MS/MS conditions: Data were acquired using DDA on a QTOF system (ZenoTOF 7600 system, SCIEX) with either CID or EAD, a novel type of ExD.¹⁻⁴

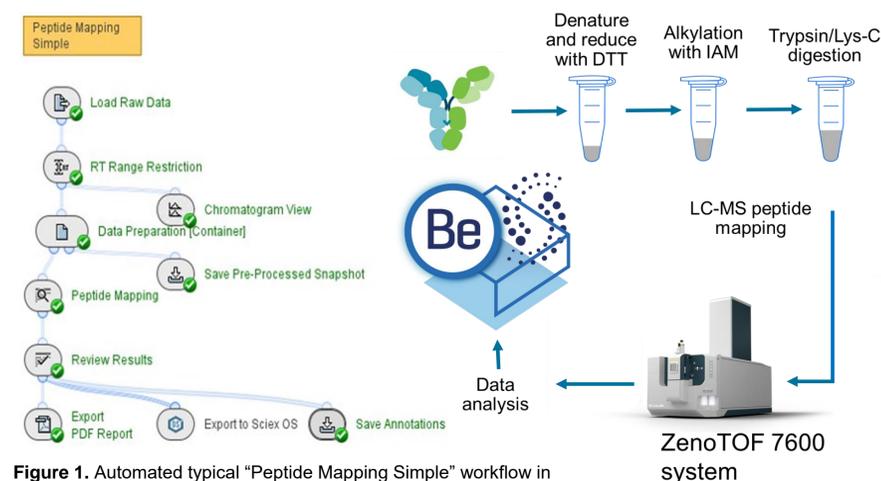


Figure 1. Automated typical "Peptide Mapping Simple" workflow in Biologics Explorer software used for N-glycopeptide analysis of CID and EAD data sets.

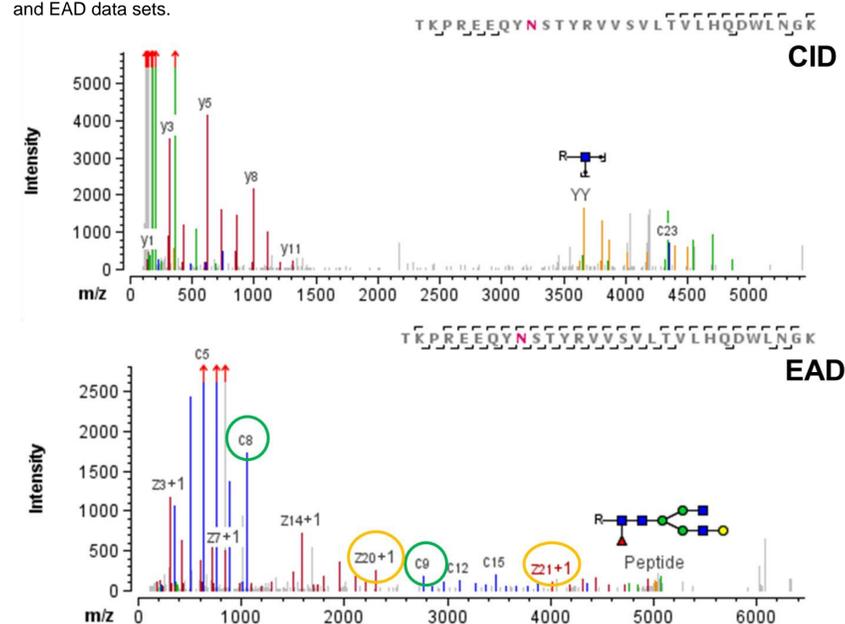


Figure 2. Comparison of MS/MS data from a glycopeptide containing G1F using CID and EAD. Top panel shows CID data, while bottom shows EAD data. Circled in green are indicative c-ions (c8, c9) for the glycan while circled in orange are the indicative z-ions (z20, z21). Blue (c- and b-ions) and red (z- and y- ions) deisotoped, deconvoluted and centroided peaks indicate the ions identified. In addition, diagnostic fragment ions confirm the localization of the glycosylation (encircled ions) in the case of EAD; whereas CID did not provide this information. In the CID spectrum red peaks show peptide fragments (y-ions) and orange peaks show oxonium ions of glycan portion.

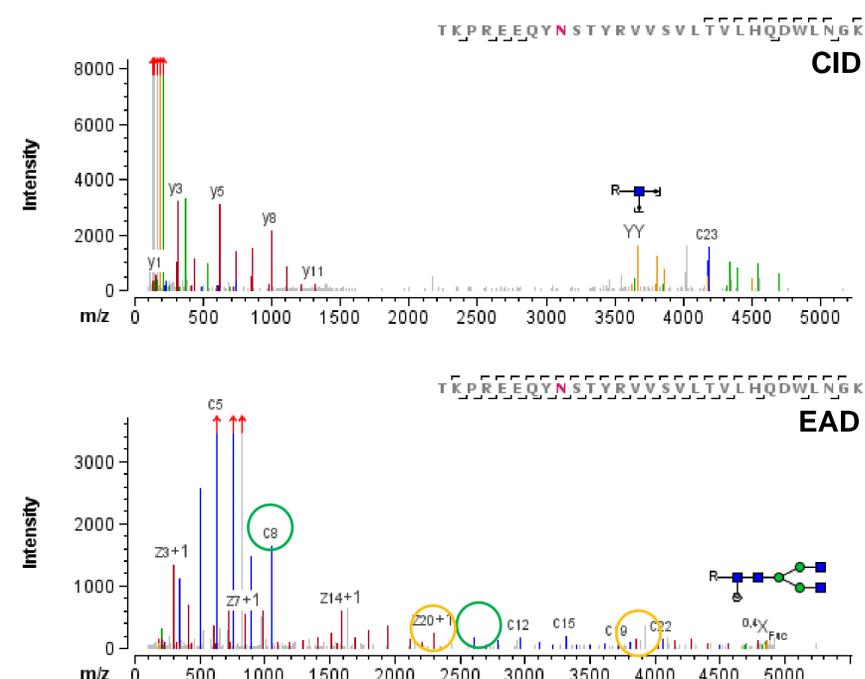


Figure 3. Comparison of CID and EAD MS/MS data for glycopeptide with G0F. Top: CID did not provide descriptive fragment information, neither positional information for the glycosylation. Bottom: close to 100% fragment coverage was achieved with EAD and diagnostic fragment ions confirm the localization of the glycosylation (encircled ions c8/c9 and z20/z21).

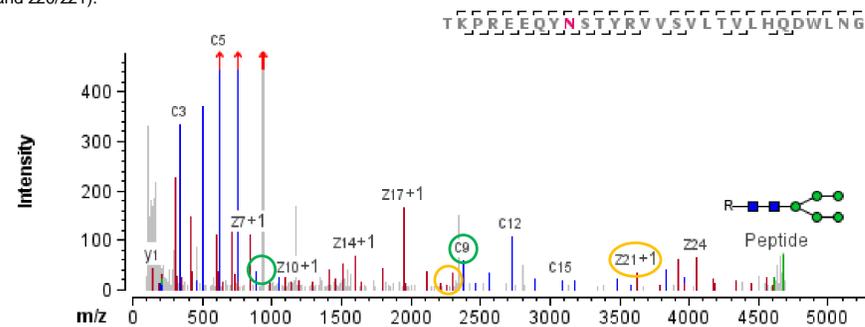


Figure 4. EAD MS/MS spectrum for glycopeptide with Man5. MS/MS spectrum is shown for a low abundance glycopeptide (z =+5) of ~3% relative abundance. EAD achieved extensive fragment coverage and localization of the glycosylation site. Diagnostic fragments c8/c9 and z20/z21 are circled in the spectrum.

Modifi...	20210130 Herceptin IDA ECD ...	20210130 Herceptin IDA ECD ...	20210130 Herceptin IDA ECD Most intense _3 [%]
G0	3.23	3.20	3.13
G0-GlcNAc	0.57	0.58	0.60
G0F	45.08	46.09	45.48
G0F-GlcNAc	6.03	6.26	6.42
G1	1.08	1.07	1.03
G1F	34.88	34.60	35.17
G2F	4.36	4.32	4.41
Man5	3.27	3.35	3.24
Man6	0.42	0.43	0.40

Figure 5. Identification summary of N-linked glycosylations in trastuzumab at N300. The table from the Biologics Explorer software summarizes the identified glycan species based on MS/MS with EAD and the relative abundance based on the XIC of the MS1 for 3 replicate injections. G0F and G1F are the most abundant glycoforms observed. Trace level of high mannose were also detected. The table confirms high reproducibility from high-abundance modified peptides to very low-abundance forms.

CONCLUSIONS

- The robust, reproducible and easy-to-use alternative fragmentation mechanism of EAD enables users to identify, fully characterize and relatively quantify glycopeptides along with a general peptide mapping analysis in 1 injection
- Excellent fragment coverage and localization of fragile modifications can be achieved with Zeno EAD with very high reproducibility, allowing for full confidence in peptide ID
- MS/MS fragment detection can be significantly enhanced compared to traditional MS/MS analyses, enabling collection of high-quality data for confident fragment assignment utilizing Zeno EAD⁴⁻⁵
- Automatic data processing enables the routine and advanced characterization of complex biotherapeutics and standard mAbs in a reproducible manner using the new Biologics Explorer software from SCIEX⁶

REFERENCES

- Baba T et al. (2015) Electron capture dissociation in a branched radio-frequency ion trap. *Anal Chem.* **87**, 785–792.
- Baba T et al. (2021) Dissociation of biomolecules by an intense low-energy electron beam in a high sensitivity time-of-flight mass spectrometry. *J. Am. Soc. Mass Spectrom.* **32**(8), 1964–1975.
- Comprehensive peptide mapping of biopharmaceuticals utilizing electron activated dissociation (EAD). [SCIEX technical note, RUO-MKT-02-12639-B](#).
- A comprehensive peptide level characterization of a multispecific monoclonal antibody (mAb). [SCIEX technical note, RUO-MKT-02-12975-B](#).
- Evaluation of biotherapeutic sequence coverage using electron activated dissociation (EAD). [SCIEX technical note, RUO-MKT-02-1290-A](#).
- Peptide mapping analysis of etanercept – a highly glycosylated dimeric fusion protein. [SCIEX technical note, RUO-MKT-02-12653-A](#).

TRADEMARKS/LICENSING

The SCIEX clinical diagnostic portfolio is For In Vitro Diagnostic Use. Rx Only. Product(s) not available in all countries. For information on availability, please contact your local sales representative or refer to www.sciex.com/diagnostics. All other products are For Research Use Only. Not for use in Diagnostic Procedures. Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd. or their respective owners in the United States and/or certain other countries (see www.sciex.com/trademarks).

© 2022 DH Tech. Dev. Pte. Ltd. RUO-MKT-10-14691-A.

To receive a copy of this poster:

- Scan the code with your phone camera
- Complete the form

