



Characterization of sialylated N-glycopeptide isomers on RP LC-EAD mass spectrometry



Suya Liu, Zoe Zhang, Pavel Ryumin and Takashi Baba

2023 June 8

DOC-MKT-28075-A © 2023 DH Tech. Dev. Pte. Ltd.

Biological significance of sialic acid

- Sialic acid has important biological functions
 - Cell–cell interaction
 - Immune reactions
 - Development processes
 - Virus host cell interaction
- Sialic acid linkage is also very important

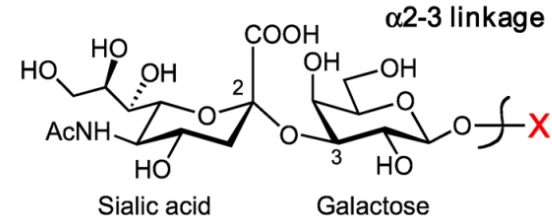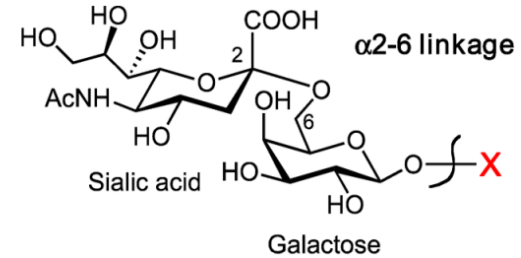


Cited from Zhu, X.Y. et al, [dx.doi.org/10.1021/cb5004114](https://doi.org/10.1021/cb5004114) | ACS Chem. Biol. 2014, 9, 1877–1884

LC separation of sialylated N-glycopeptide isomers

MODEL GLYCOPEPTIDE FROM FETUIN DIGEST


- Current LC methods
 - HILIC, PGC, RP HPLC with high column temperature
- Sample: fetuin digest, LCPDCPLLAPLNDSR

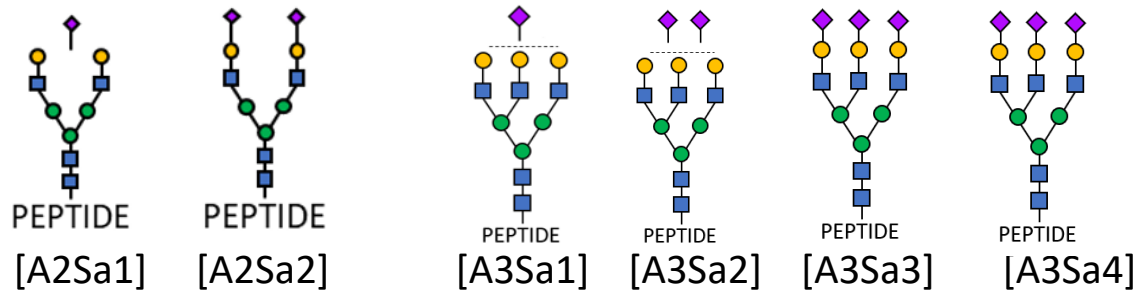

— Sialic acid  Gal  GLcNac  Man 

Optimization to achieve superior separation of sialylated N-glycopeptide isomers

LCPDCPLLAPLN[A3SA3]DSR, M/Z = 1151.215, +4

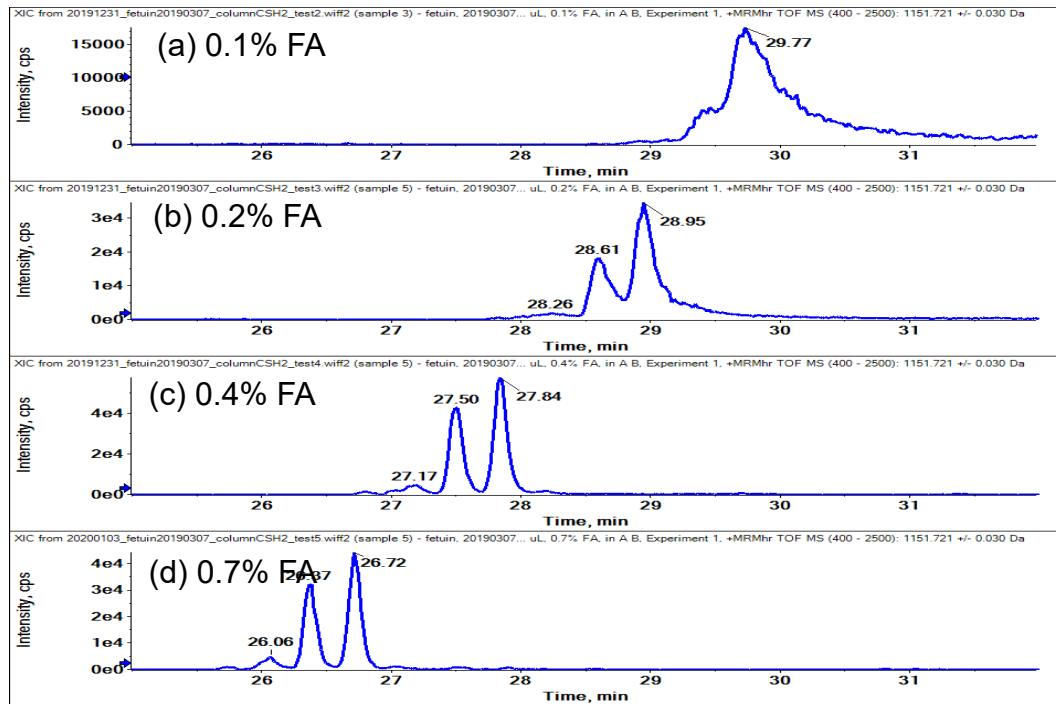

Shimadzu Prominence LC

ACQUITY UPLC

CSH C18 column (130 Å,
1.7 µm, 2.1 mm × 100
mm, 148 Waters)

A: 0.1 % FA water





B: 0.1% FA acetonitrile

Flow rate 0.15 mL/min

Zeno TOF 7600 system

Optimization to achieve superior separation of sialylated N-glycopeptide isomers

LCPDCPLLAPLN[A3SA4]DSR, $M/Z = 1223.99$, +4

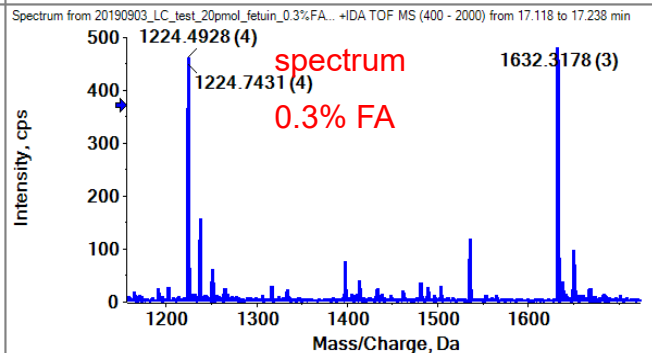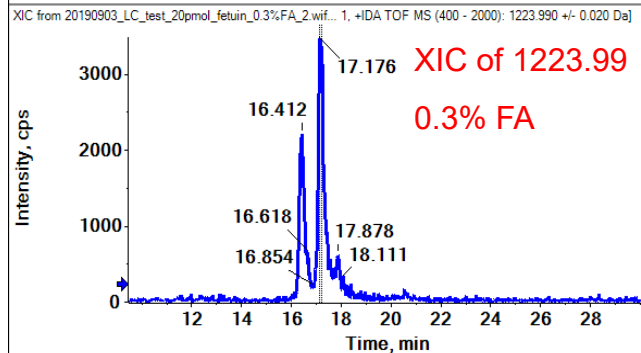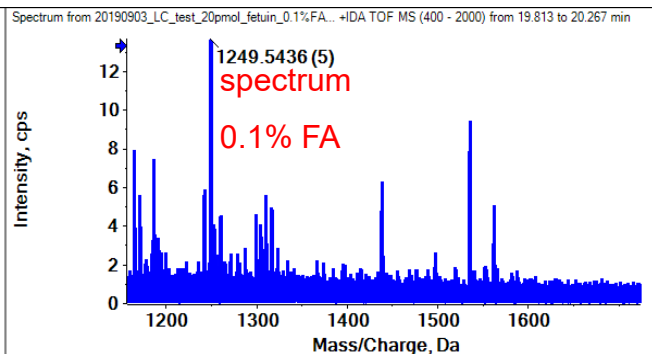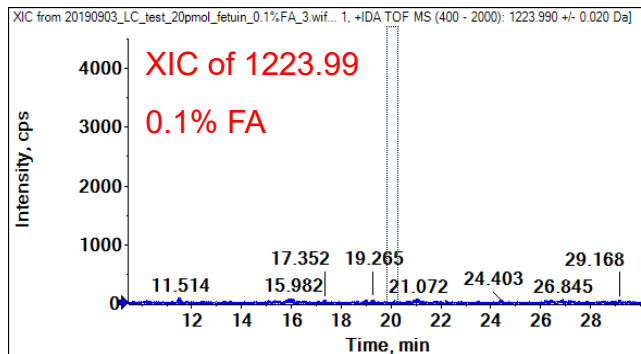

The glycopeptide was only detected when >0.3% FA in the mobile phase


EAD of glycopeptide characterization

LCPDCPLLAPLN[A3SA2]DSR, M/Z = 987.17, +4

4585.8	4472.7	4312.7	4215.7	4100.6	3940.6	3843.5	3730.5	3617.4	3546.3	3449.3	3336.2	361.2	246.1	159.1
Z15	Z14	Z13	Z12	Z11	Z10	Z9	Z8	Z7	Z6	Z5	Z4	Z3	Z2	Z1
L	C	P	D	C	P	L	L	A	P	L	N-glycan	D	S	R
C1	C2	C3	C4	C5	C6	C7	C8	C9	C10	C11	C12	C13	C14	C15
131.1	291.1	388.2	503.2	663.3	760.3	873.4	986.5	1057.5	1154.6	1267.7	3586.5	3701.5	3788.5	3944.5

LC-EAD on Zeno TOF 7600 KE 7,



Spectrum from 20190222_s1MRM_fetuin_67_100ms_CE10_ETC0_087_7_Z_60min_20.pnd.vi12 (sample 1) - fetuin 20.pnd, Experiment 25, 987.17, s1MRM TOF MS/MS of 9872 (200 - 2000) from 26.235 to 26.431 min, Gaussian smoothed (1.0 points)

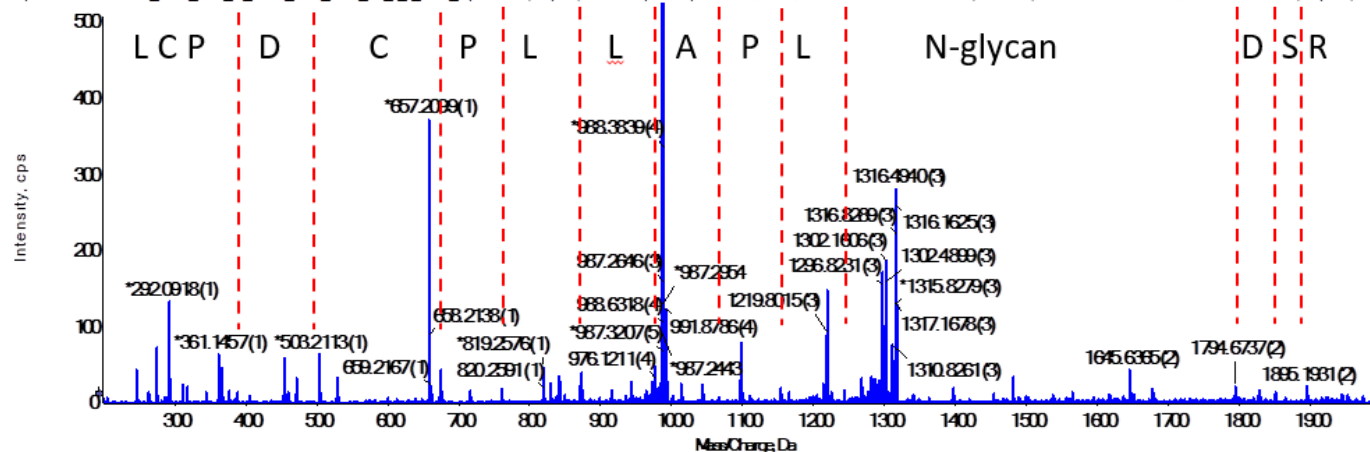

Identification of sialic acid linkage isomers

LCPDCPLLAPLN[A3SA]DSR, $M/Z = 914.38$, +4


TIC from 20190927_Fetuin_20pmol_MRM_91...sMRMhr TOF MSMS of 914.4 (100 - 2500)

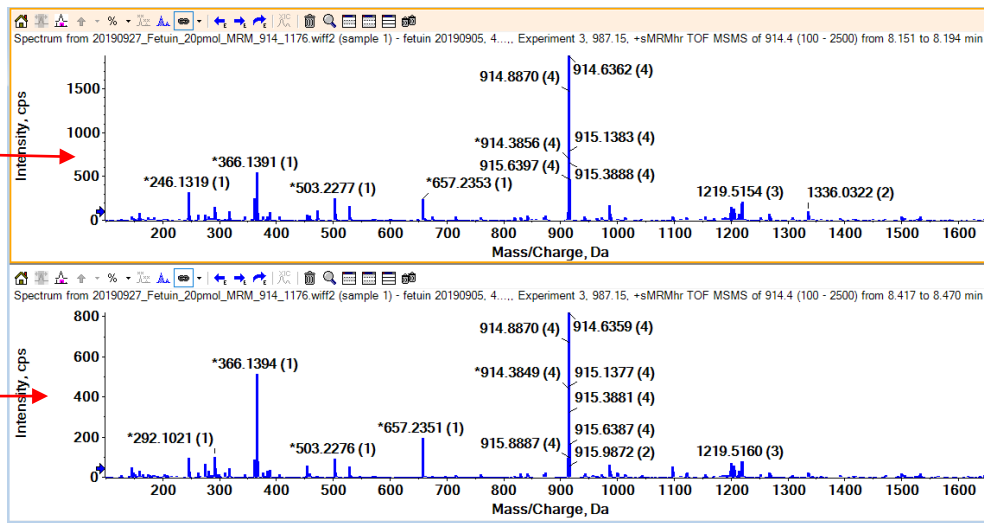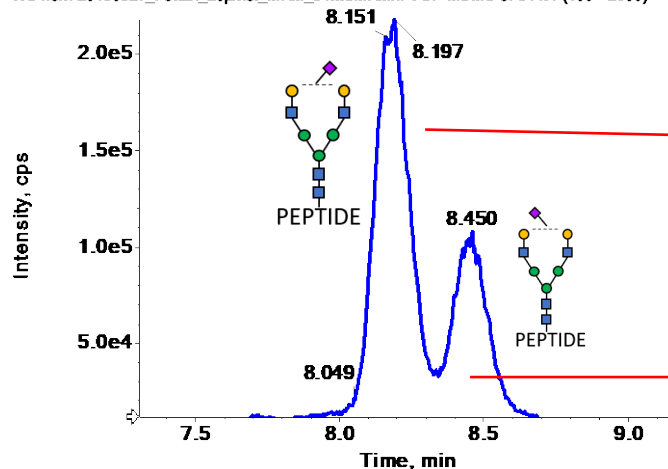



The two isomers actually generated the same m/z fragment ions

Identification of sialic acid linkage isomers

FRAGMENTATION OF SIALIC ACID LINKED GLYCANS

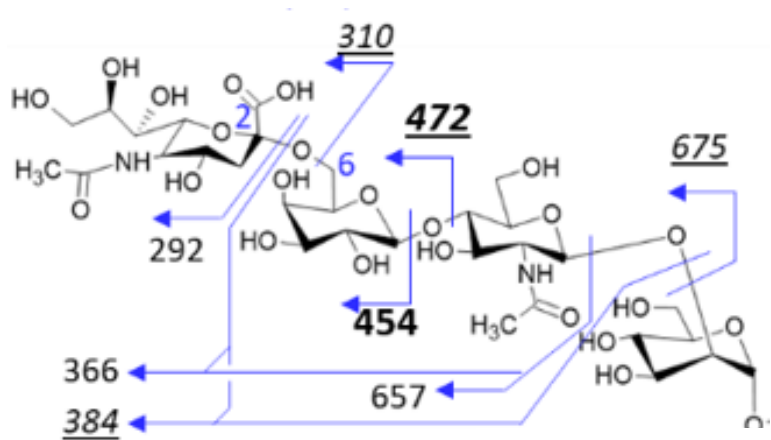


Neu5Ac - $\alpha(2-6)$ - Gal

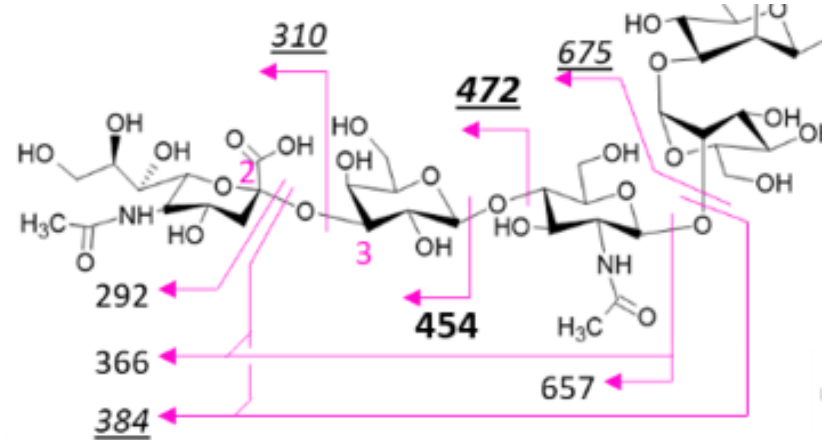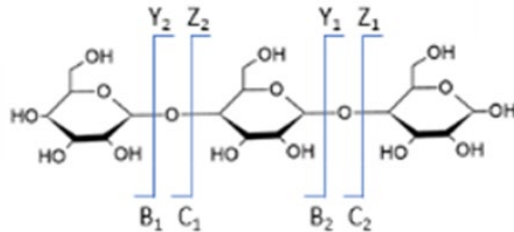

Neu5Ac - $\alpha(2-3)$ - Gal

Identification of sialic acid linkage isomers

LCPDCPLLAPLN[A2SA]DSR M/Z = 914.38, +4

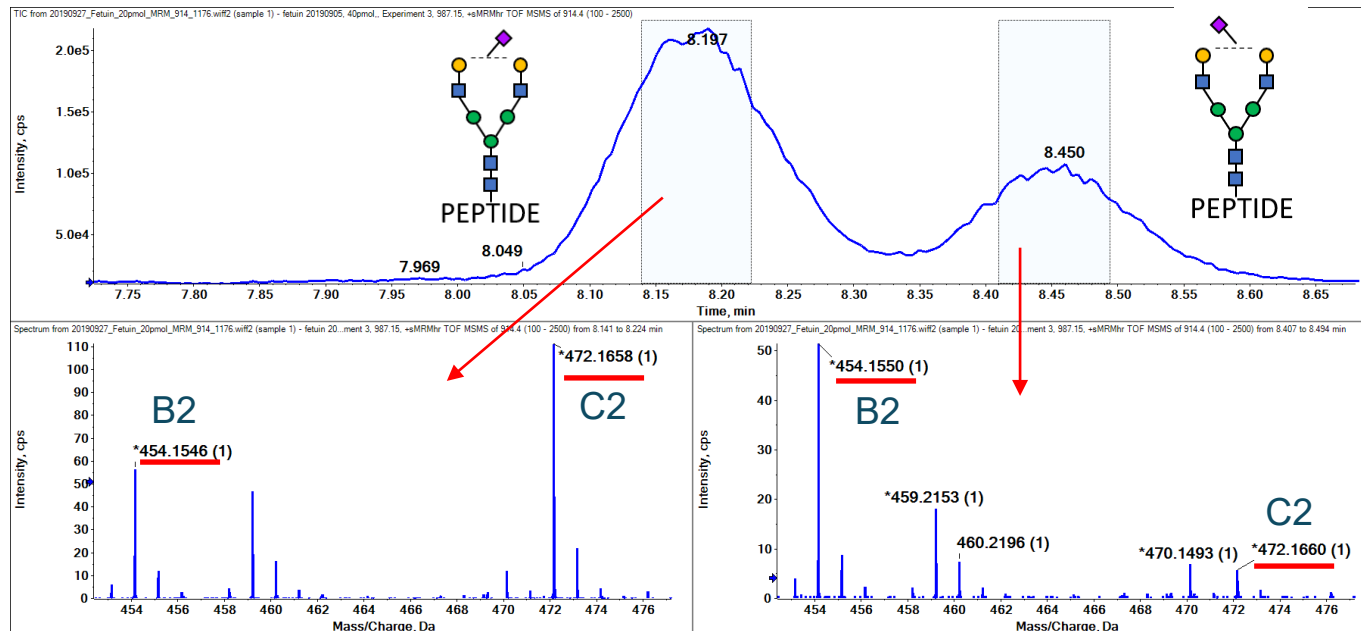


However, the fragment ion ratios showed difference between isomers

$$\begin{aligned} 472/454 &= 2 \quad \text{for } \alpha 2-6 \\ 472/454 &= 0.2 \quad \text{for } \alpha 2-3 \end{aligned}$$

Identification of sialic acid linkage isomers

B AND C ION INTENSITY VS. KE

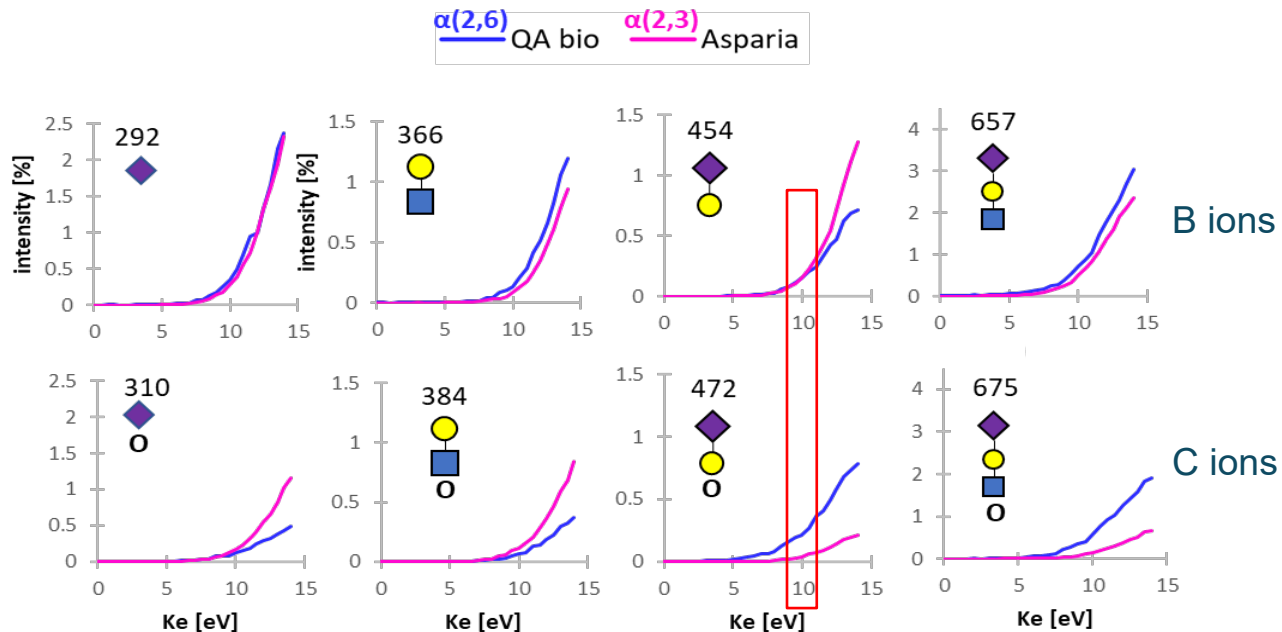

KVAN[A2Sa2]KT

QA Bio $\alpha 2-6$
Asparia $\alpha 2-3$

C and B ion intensity normalized to charge reduced precursor

Identification of sialic acid linkage isomers


OPTIMIZATION OF ECD KE FOR SIALIC ACID LINKAGE ISOMERS

- Diagnostic parameter (DP) for linkage differentiation
 - $$DP = \frac{\text{intensity of B2 ion at 454}}{\text{intensity of C2 ion at 472}}$$
- Ke 10 V is optimal for the linkage differentiation of the glycopeptides studied
- Other glycan ions are helpful to assign the Sa linkage and position
 - Glycan specific fragment ions, precursor - sugar ion



Identification of sialic acid linkage isomers

LCPDCPLLAPLN[A2SA]DSR, M/Z = 914,38, +4

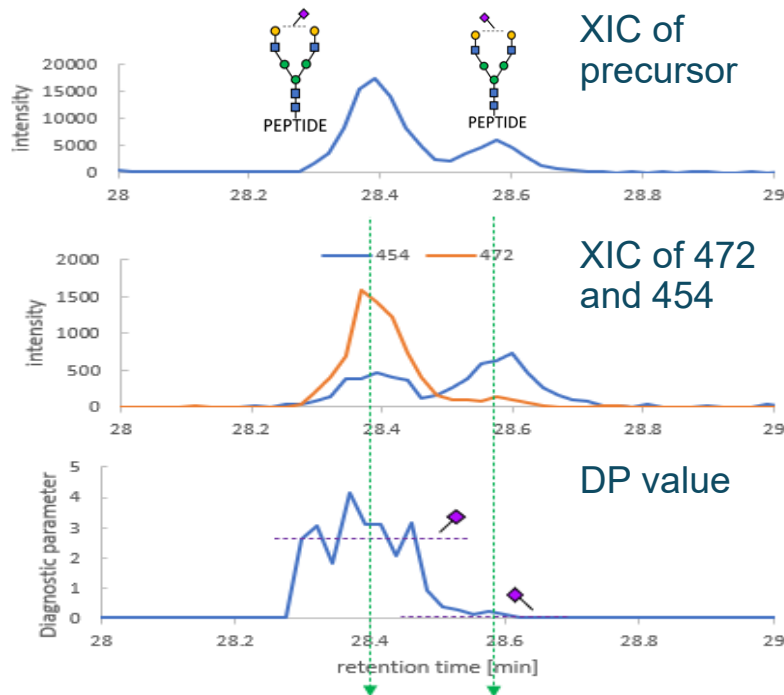


For this single sialylated glycopeptide
the DP clearly assigned the α 2-3 and
 α 2-6 isomers

$$DP = \text{intensity } 472 / \text{intensity } 454$$

Identification of sialic acid linkage isomers

LCPDCPLLAPLN[A2SA2]DSR, M/Z = 987, 17, +4

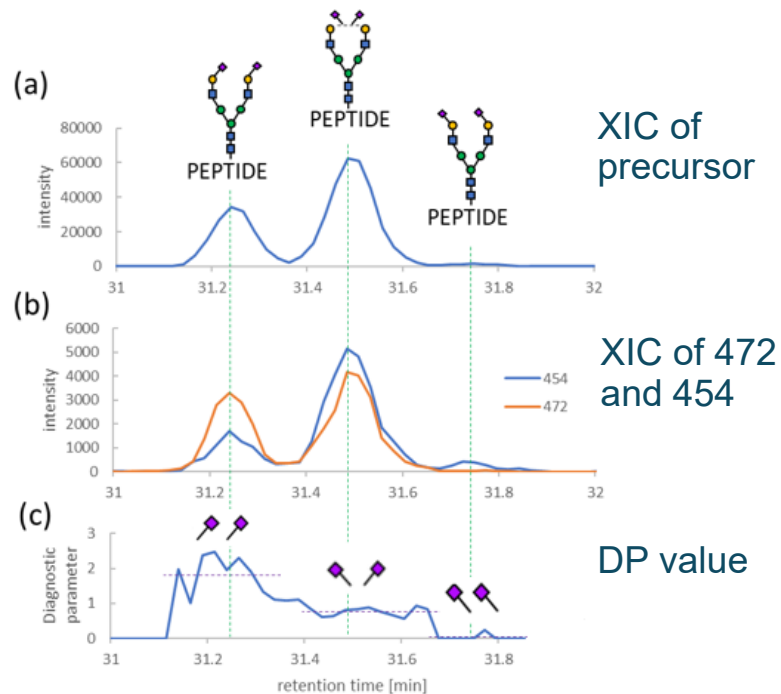


For this double sialylated glycopeptide the DP clearly assigned the α 2-3 and α 2-6 isomers

The DP value for the middle peak depends on the ratio of α 2-3 and α 2-6 linkage


Identification of sialic acid linkage isomers

LCPDCPLLAPLN[A2SA3]DSR M/Z = 1151.21, +4

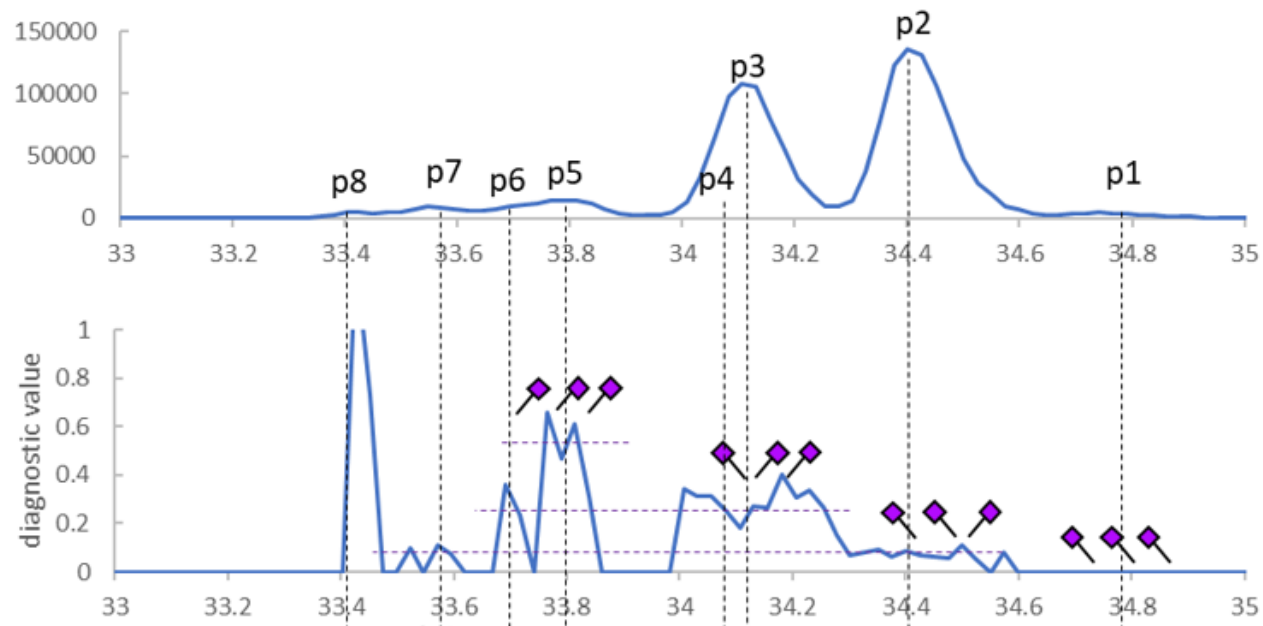

Identification of sialic acid linkage isomers

LCPDCPLLAPLN[A2SA3]DSR M/Z = 1151.21, +4

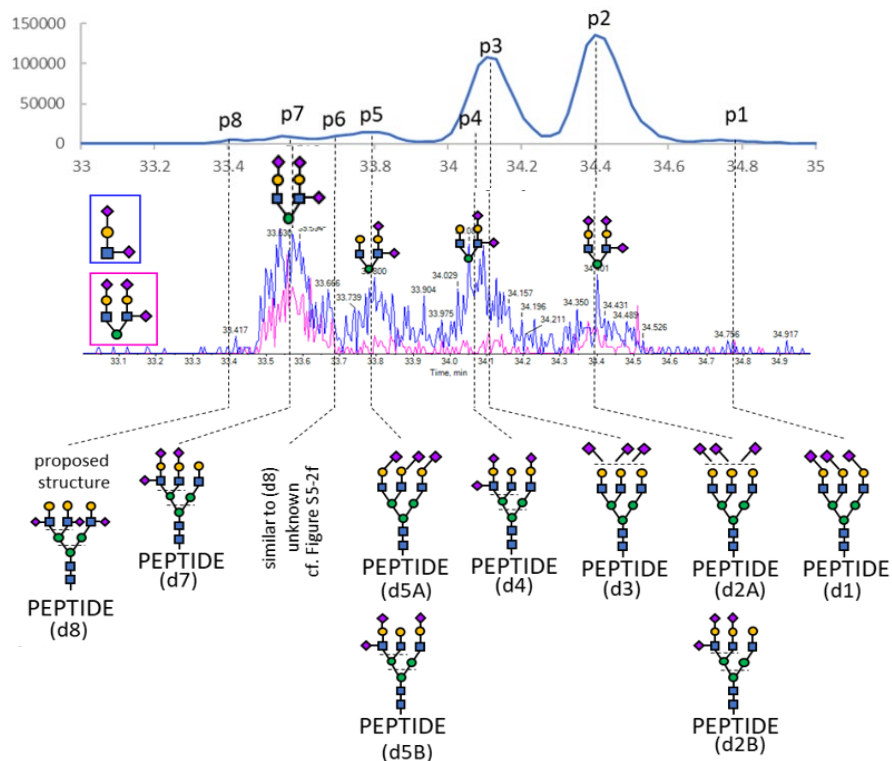

Identification of sialic acid linkage isomers

CONCLUSION

- Single RP LC MS method for sialylated N-glycopeptide
 - 0.3-0.4 % FA in mobile phase
 - High sensitivity
 - Better separation of sialic acid position and linkage isomer
- EAD generated B and C glycan fragment ions, ratio can be used to differentiate Sa linkage and position isomer

Ref: Liu, S.; Simmons, D.; Ryumin, P.; Baba, T. 69th ASMS Conference on Mass Spectrometry and Allied Topics; ASMS, 2021; p 265

Ref: Liu, S.; Ryumin, P.; Albanese, J.; Zhang, Z.; Baba, T. *Analytical Chemistry* **Article ASAP**, DOI: 10.1021/acs.analchem.2c04581

Thank you!



Trademarks / Licensing

The SCIEX clinical diagnostic portfolio is For In Vitro Diagnostic Use. Rx Only. Product(s) not available in all countries. For information on availability, please contact your local sales representative or refer to www.sciex.com/diagnostics. All other products are For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd. or their respective owners in the United States and/or certain other countries (see www.sciex.com/trademarks).

© 2023 DH Tech. Dev. Pte. Ltd. **DOC-MKT- 28075-A**

