

LC-MS/MS Screening and Quantification Methods for the Analysis of 41 Common Pain Drugs in Oral Fluid

Using SCIEX QTRAP® LC-MS/MS systems and technology

Introduction

Liquid chromatography coupled to tandem mass spectrometry (LC-MS/MS) is a widely used analytical tool for screening and confirmation of drugs of abuse in forensic samples. The detection of several compound classes, such as common pain panel and drugs of abuse is of much importance. While this analysis has generally been performed utilizing urine sample matrices, many research laboratories are investigating the analysis of these compounds in oral fluid matrices. Using oral fluid as a matrix provides several advantages over urine analysis. It provides a simpler collection procedure that can be witnessed which means there is less opportunity to adulterate the sample. Typical collection of oral fluid matrix involves a collection swab that retains the oral fluid matrix. After collection, the swab is subjected to extraction buffer. The resulting matrix is then analyzed for the drug compounds. The outcome for these investigations into the use of oral fluid needs to provide a means to accurately and precisely quantify compounds in a timely fashion, while having a simplified workflow that is conducive to their setting. This current effort has shown to accomplish just that. In this study we present a fast, robust, and reliable method that has afforded the detection of 41 compounds in an oral fluid matrix. The compounds included 6-MAM, alprazolam, amphetamine, benzoylecgonine, buprenorphine, carisoprodol, clonazepam, codeine, diazepam, EDDP, fentanyl, flunitrazepam, flurazepam, hydrocodone, hydromorphone, hydroxalprazolam, lorazepam, MDA, MDEA, MDMA, meperidine, meprobamate, methadone, methamphetamine, midazolam, morphine, naloxone, naltrexone, norbuprenorphine, nordiazepam, norfentanyl, normeperidine, norpropoxyphene, oxazepam, oxycodone, oxymorphone, PCP, propoxyphene, sufentanil, temazepam and tramadol.

Key Features of SCIEX QTRAP® Multiplexed LC-MS/MS Systems for the identification and quantification of designer drugs.

- The QTRAP system offers high sensitivity data acquisition strategies for both quantitative and qualitative experiments
- System sensitivity allows small sample volumes and simplified sample processing procedures



QTRAP® system for quantification, identification and confirmation of analytes

- High-throughput quantitative analysis is achieved using MRM (Multiple Reaction Monitoring) and the Scheduled MRM™ Algorithm
- MRM triggered EPI (Enhanced Product Ion) scans generate high sensitivity MS/MS spectra to enhance confidence of identification of analytes, even at low concentrations.
- MPX™ 2.0 High Throughput System allows significant throughput increases by maximizing utility of available MS analytical windows focusing only on areas of the chromatogram showing peak elution.

Experimental

Calibrator Preparation: An oral fluid matrix was spiked to various levels of the relevant compounds. The calibration curve extended above and below the industry accepted cutoff and confirmation levels.

Sample Preparation:

- Add 150 µL preserved buffer to 50 µL oral fluid matrix
- Add 200 µL MeOH:H₂O:Formic Acid containing internal standards
- Vortex, Centrifuge and transfer to an autosampler vial for injection

HPLC Conditions: Chromatography was performed using a Phenomenex Kinetex C18 2.6 μm 50 x 3.0 mm column using ammonium formate in H_2O as mobile phase A and formic acid in methanol as mobile phase B. A gradient was applied during chromatographic run and the total run time (not multiplexed) per stream was 6 minutes. The acquisition window was set to 3.5 minutes per stream (Figure 1) providing an increase in sample throughput of 171% during multiplexed analysis.



Figure 1. MPX chromatographic method. By applying MPX, the injection-to-injection time can be shortened to 3.5 min

MS/MS Conditions: MS/MS analysis was carried out using a QTRAP® system operating in positive electrospray ionization mode. Analytes were analyzed using the Scheduled MRM Algorithm (Figure 2) to accommodate for quantification of all components listed above. In addition to the Scheduled MRM™ Algorithm (sMRM) quantification, an information dependent acquisition (IDA) method was set up for targeted drug screening which involved an sMRM survey scan followed by enhanced product ion scans (EPI), the details of which are shown in Figure 3.

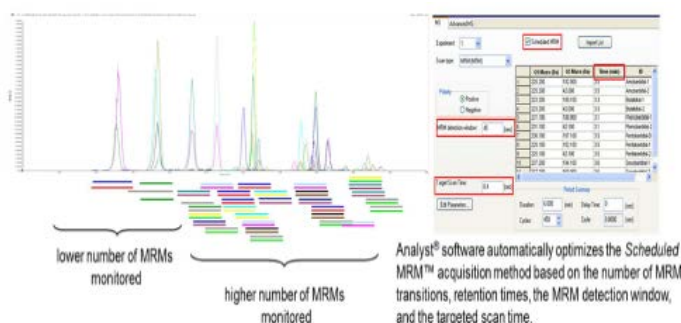


Figure 2. The Scheduled MRM™ Algorithm detects MRM transitions only during a specified time window. This allows many more MRM transitions to be monitored, while maximizing dwell time cycle times.



Figure 3. The sMRM quantitative method for all 41 pain drugs and internal standards (left) and the sMRM-IDA-EPI drug screening method for all 41 pain drugs (right). All experiments were carried out on a SCIEX QTRAP® LC-MS/MS system.

Results and Discussion

Figures 4 and 5 show typical chromatograms achieved using the described approach and quantitative results for the following analytes: codeine, methamphetamine, midazolam and temazepam. Excellent linearity and reproducibility were observed throughout the dynamic range assessed in this effort. Table 1 shows a summary of the calibration range for all 41 pain drugs. Within the listed concentration range, excellent linearity and reproducibility were achieved.

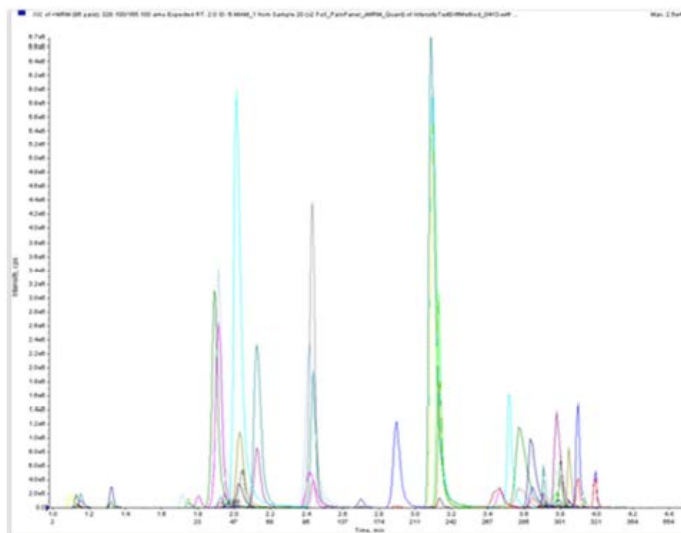


Figure 4. A typical chromatogram of all 41 analytes and internal standards

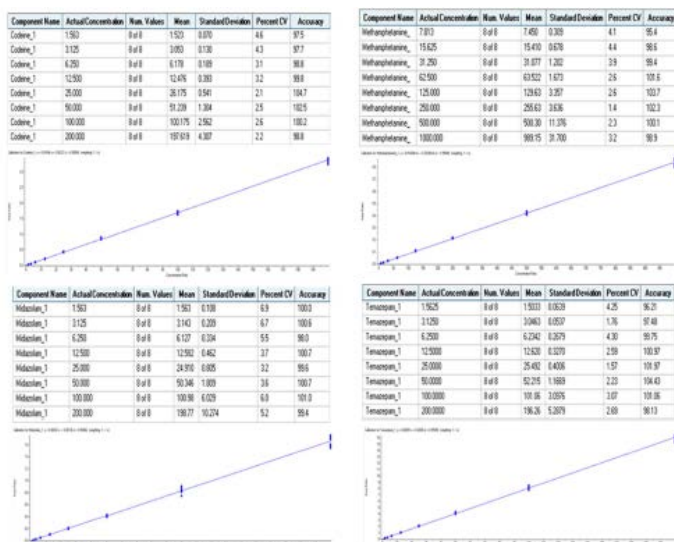


Figure 5. Quantitative analysis examples, codeine (top left), methamphetamine (top right), midazolam (bottom left) and temazepam (bottom right).

Figure 6 shows an excerpt from a report generated for the results of the analysis performed using a QTRAP® LC-MS/MS system in which both the triple quadrupole and linear ion trap functionalities have been combined in one MS method to perform both quantification of detected compounds and confident identification of the compounds through MS/MS spectral library matching.

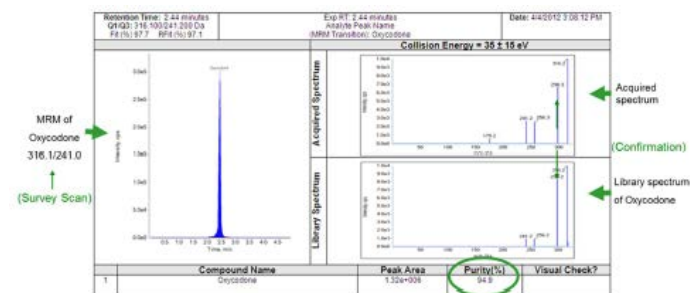


Figure 6. Example of the targeted screening of all 41 common pain drugs. Shown on the left is the chromatography of oxycodeone. Shown on the right is the comparison of acquired EPI spectrum and library spectrum of oxycodeone.

Conclusions

A fast, robust and simple method for the detection of 41 common pain drugs in an oral fluid matrix was developed on a SCIEX QTRAP® LC-MS/MS system. This method utilized a dilute-and-shoot procedure with minimum sample prep needed. Excellent linearity and precision were observed for all 41 compounds in the listed calibration range. Due to the fast scanning speed and trapping capability of the QTRAP® system, a targeted screening method was also developed to accomplish a screen for these 41 common pain drugs.

The application of the MPX™ 2.0 High Throughput System greatly reduced the sample run time and increased sample throughput by 171%. The QTRAP® LC-MS/MS system is a powerful tool for ultra-fast multi-analyte quantification and identification in forensic samples. The high-speed scanning capabilities permit the use of very short dwell times and pause times, resulting in the ideal detector for the ultra-fast LC methods being employed in today's modern laboratories.

Analysis times can be reduced, therefore sample throughput increased, while maintaining data quality (sensitivity, reproducibility, and linear dynamic range, with no crosstalk). This unique LC-MS/MS method utilizing the Scheduled MRM™ Algorithm in combination with acquisition of MS/MS spectra for compound identification through MS/MS library spectral searching allows for the analysis of the ever increasing number of compounds that are required to be quantified in a single method.

The SCIEX clinical diagnostic portfolio is For In Vitro Diagnostic Use. Rx Only. Product(s) not available in all countries. For information on availability, please contact your local sales representative or refer to <https://sciex.com/diagnostics>. All other products are For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein are the property of AB Sciex Pte. Ltd. or their respective owners in the United States and/or certain other countries.

© 2020 DH Tech. Dev. Pte. Ltd. RUO-MKT-02-10918-B. AB SCIEX™ is being used under license.