

Sequencing Analysis as A Tool for Fast and Accurate Microbial Contaminants Identification

• High Specificity • Rapid Results • Superior Over Conventional Microbial Methods

J. Luo¹, K. Uchiumi¹, M. Souquet² and H.Yowanto¹

¹ SCIEX, Brea, California, USA

² SCIEX Darmstadt, Germany

Introduction and Overview

During the development and manufacturing processes of biologics, it is critical to monitor the sterility of pharmaceutical ingredients, water for pharmaceutical use, the manufacturing environment, intermediates and finished products. In addition, Master Cell Bank, each seed lot and cells used in each production run should be tested for adventitious agents including mycoplasma, bacteria, fungi, viruses, and virus-like particles. When microbial growth is observed, an investigation needs to be conducted and FDA guideline requires sterility test isolates be identified to the species level. It has been shown that DNA sequencing is superior in identifying microbes to the species level, when compared against conventional microbiology staining and culture methods. In this poster, we describe a process for sequencing ribosomal RNA genes of microorganisms that offers fast turnaround time (8 hrs) and high microbial identification accuracy. Upon isolation of nucleic acid from each microbial sample, the gene target is amplified using a polymerase chain reaction (PCR), followed by sequencing reaction using the Dye Terminator Cycle Sequencing (DTCS) on the GenomeLab GeXP Genetic Analysis System. Results are compared against the NCBI microbial library to identify the contaminants.

Methods

Genomic DNA Isolation: One loopful of bacteria sample from a single colony was re-suspended in 100 µl of sterile water in a microcentrifuge tube. The tube was placed in a boiling water bath for 10 minutes. After a minute spin at 10,000 g, 5 µl of the supernatant containing microbial DNA was used as the PCR template. For yeast sample, Genomic DNA was extracted using Beckman Coulter Agencourt Genfind V2 kit with a lysis buffer that contained lysozyme and Lyticase from Sigma.

Target Gene Amplification: For Clostridium species and yeast contaminants, each PCR reaction (50 µl total volume) contained 5 µl of sample genomic DNA, 1.5 mM MgCl₂, 200 nM dNTPs, 200 nM Primers and 0.04 units of AmpliTaq Gold. Cycling conditions were: 95° C for 5 min followed by 30 cycles of 95° C for 30 sec, 50° C for 30 sec and 72° C for 1 min; followed by a final extension at 72° C for 10 min and 4° C forever. For non-pathogenic bacteria contaminant, the Takara RR180A kit was used following the manufacturer's instructions. Thermal cycling conditions were: 94° C, 1 min followed by 30 cycles of 94° C for 30 sec, 55° C for 30 sec and 72° C for 1 min; followed by a final extension at 72° C for 3 min and 4° C forever.

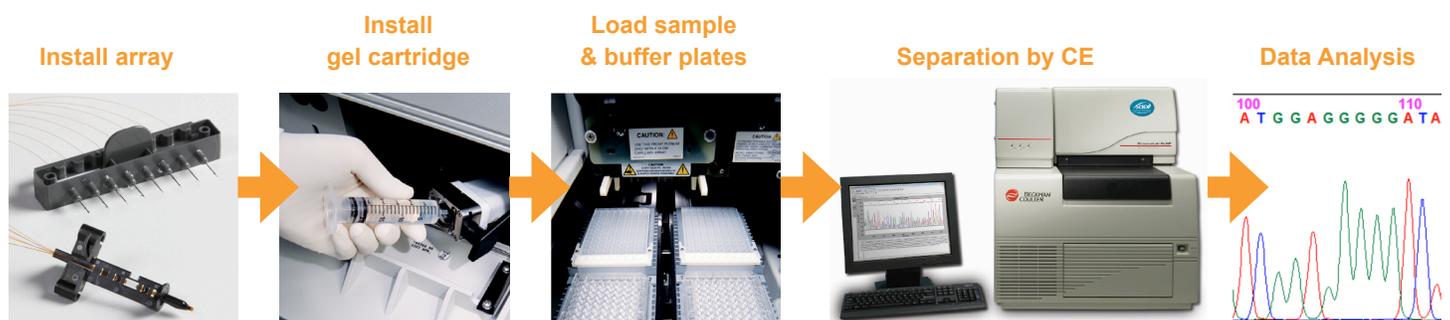


Figure 1. Basic operator steps for using the GenomeLab GeXP Genetic Analysis System.

Post PCR Clean-up: Agencourt AMPure XP from Beckman Coulter was used for post PCR clean-up. The process of Agencourt AMPure XP (Figure 2) takes less than 15 minutes to complete.

Sequencing and Data Analysis: Sequencing reactions were set up per instructions for the DTCS Quick Start Kit from SCIEX. Separation and data analysis are fully automated on the GenomeLab GeXP Genetic Analysis System. Sequences were submitted to NCBI Database for BLAST search to identify the contaminant.

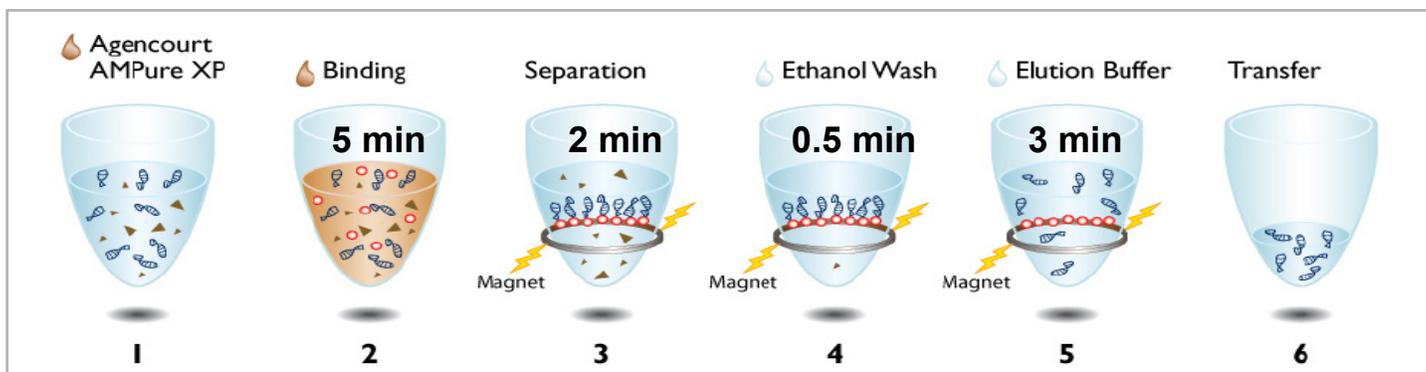


Figure 2. Post PCR clean-up process using Agencourt AMPure XP.

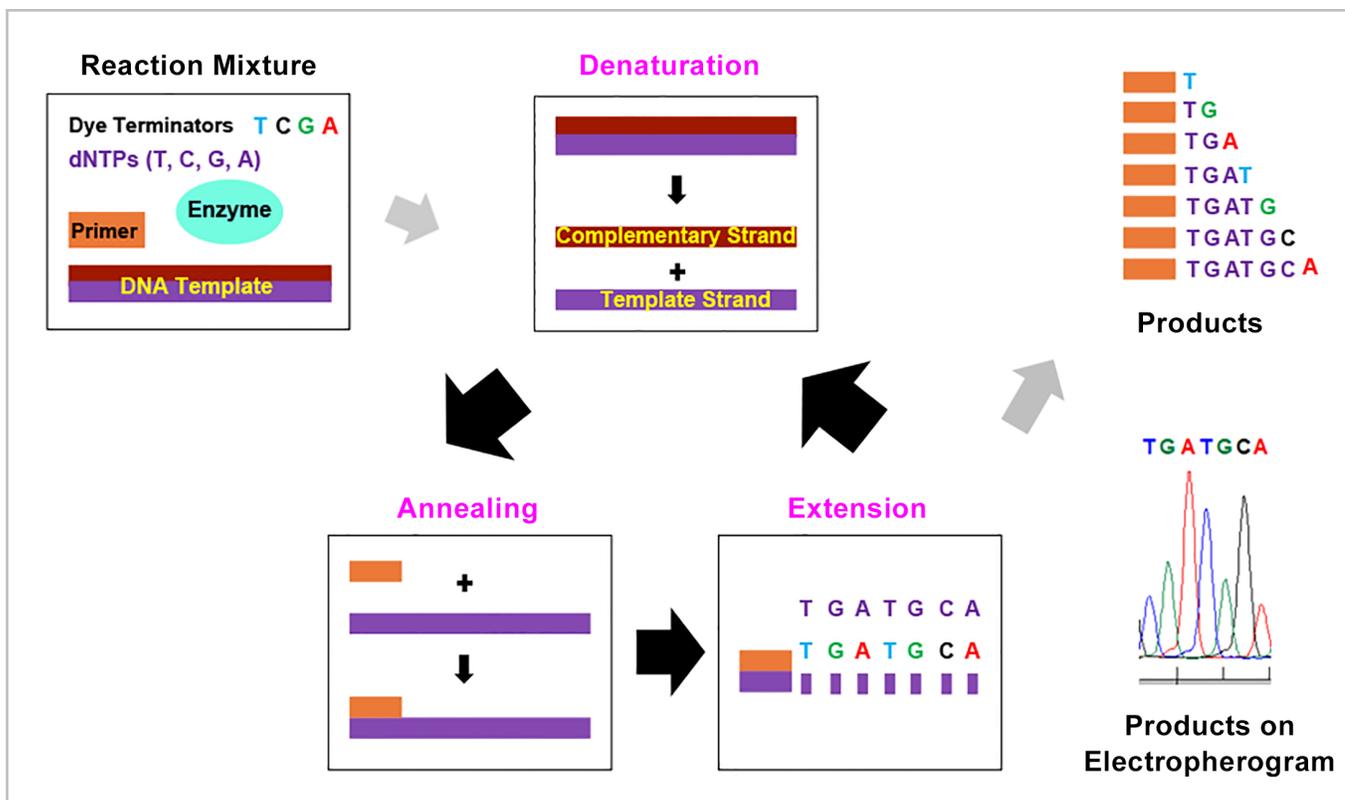


Figure 3. Sanger Dye Terminator Cycle Sequencing.

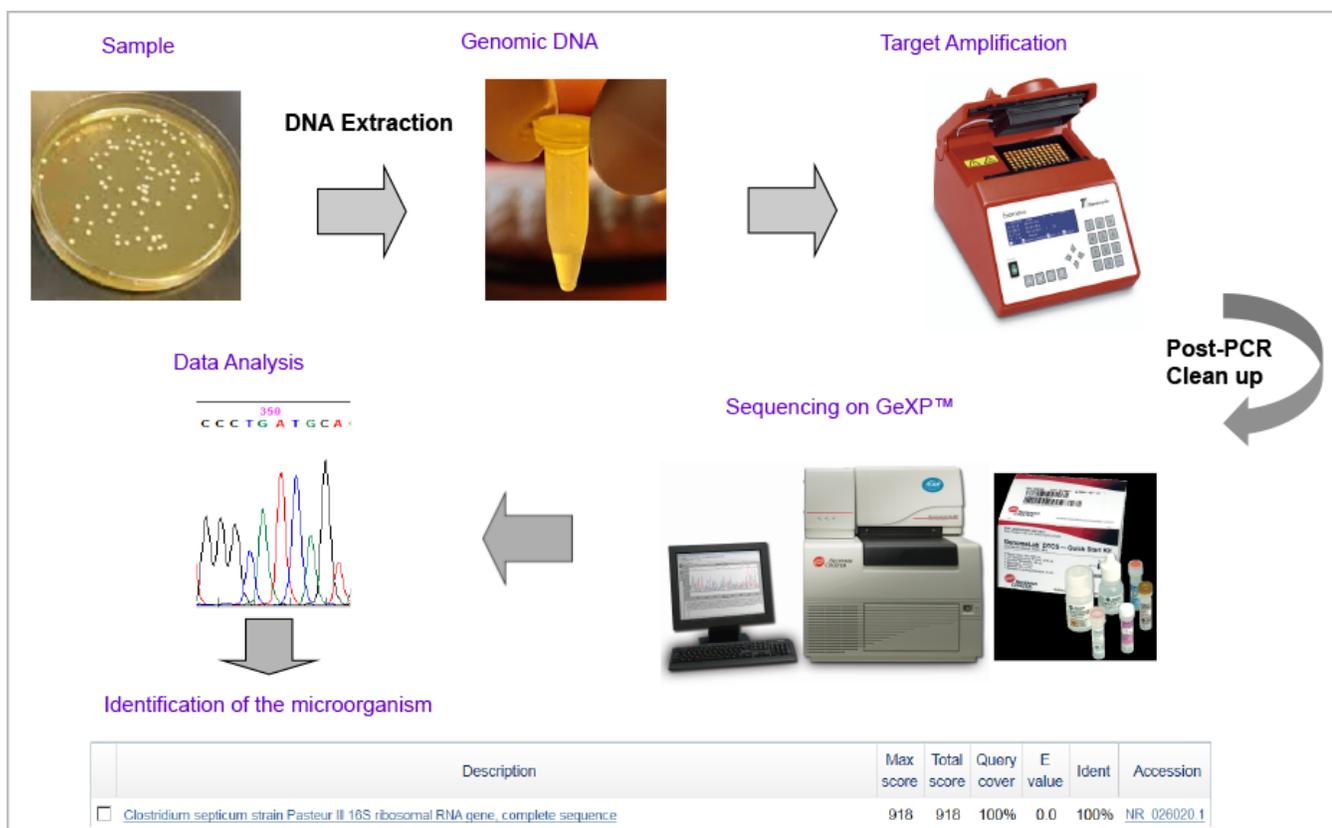


Figure 4. Work Flow for Microbial Identification using GenomeLab GeXP Genetic Analysis System.

Results

Identification of *Clostridium septicum* in a sample. The 16s rDNA PCR product was sequenced using DTCS Quick Start Kit. Sequencing fragments were purified by ethanol precipitation and separated on a GeXP instrument. Figure 5 shows the sequencing

results. After automated quality trimming, a 497bp sequence was submitted for BLAST search against the 16s rRNA database at the NCBI website. Alignment results (Figure 6) indicated that the query sequence was 100% identical to 16s rDNA from *Clostridium septicum*.

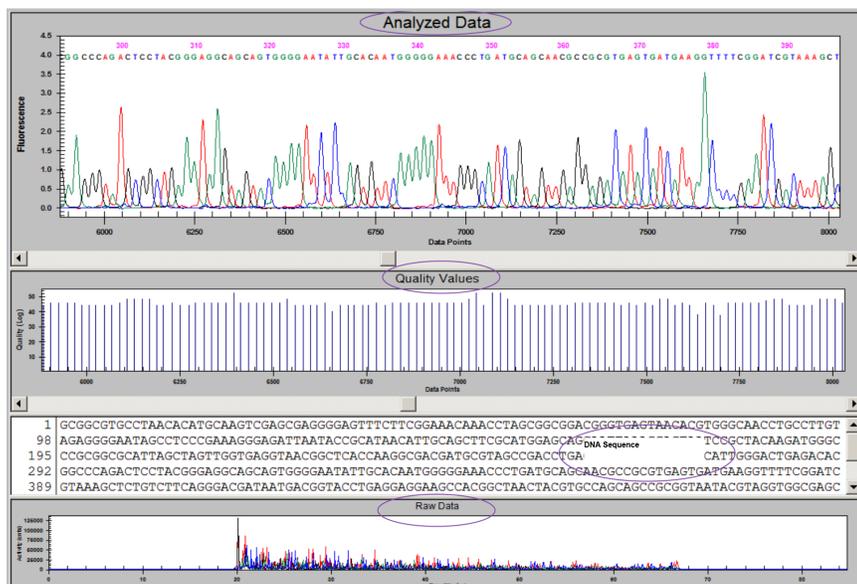


Figure 5. Sequencing results showing raw and analyzed data as well as quality values and DNA sequence.

Alignments						
Description	Max score	Total score	Query cover	E value	Ident	Accession
<input type="checkbox"/> Clostridium septicum strain Pasteur III 16S ribosomal RNA gene, complete sequence	918	918	100%	0.0	100%	NR_026020.1
<input type="checkbox"/> Clostridium chauvoei strain 2585 16S ribosomal RNA gene, complete sequence	874	874	100%	0.0	98%	NR_026013.1

Figure 6. BLAST search results obtained with a 497 bp 16s rDNA sequence in NCBI website.

Identification of *E. coli* in a sample. A non-pathogenic bacteria sample was obtained from Takara, Japan. Genomic DNA was extracted by a quick cell lysis in a boiling water bath. Full length 16s rDNA was amplified and sequenced using DTCS Quick Start Kit and GeXP System. Consensus sequence was created in

Gene Studio (Figure 7) and submitted to NCBI for BLAST search. Search results (Figure 8) showed the query sequence had high homology to pathogenic bacteria and non-pathogenic *E. coli*. Since the sample was known to be non-pathogenic, the sample was identified as an *E. coli* species.

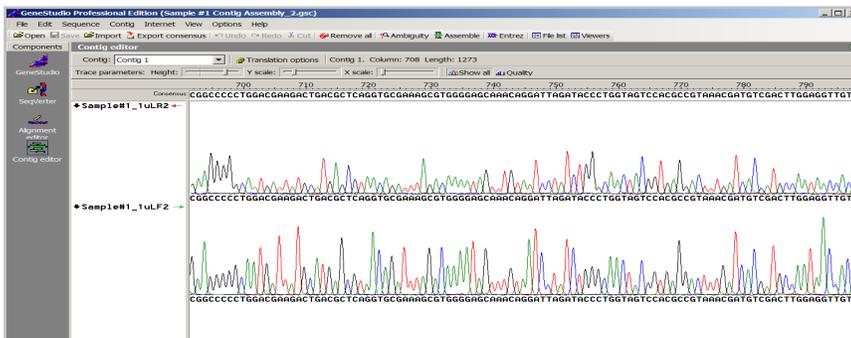


Figure 7. Contig Assembly in Gene Studio to create consensus sequence from forward and reverse sequencing results.

Sequences producing significant alignments:						
Description	Max score	Total score	Query cover	E value	Ident	Accession
<input type="checkbox"/> Escherichia fergusonii strain ATCC 35469 16S ribosomal RNA gene, complete sequence	2649	2649	100%	0.0	99%	NR_074902.1
<input type="checkbox"/> Shigella flexneri strain ATCC 29903 16S ribosomal RNA gene, partial sequence	2649	2649	100%	0.0	99%	NR_026331.1
<input type="checkbox"/> Escherichia coli strain NBRC 102203 16S ribosomal RNA gene, partial sequence	2643	2643	100%	0.0	99%	NR_114042.1
<input type="checkbox"/> Escherichia fergusonii strain ATCC 35469 16S ribosomal RNA gene, partial sequence	2639	2639	100%	0.0	99%	NR_027549.1
<input type="checkbox"/> Escherichia fergusonii strain NBRC 102419 16S ribosomal RNA gene, partial sequence	2638	2638	100%	0.0	99%	NR_114079.1
<input type="checkbox"/> Shigella sonnei strain CECT 4887 16S ribosomal RNA gene, partial sequence	2638	2638	100%	0.0	99%	NR_104826.1
<input type="checkbox"/> Shigella flexneri 2a str. 301 strain 301 16S ribosomal RNA, complete sequence	2625	2625	100%	0.0	99%	NR_074882.1
<input type="checkbox"/> Escherichia coli O157:H7 str. Sakai strain Sakai 16S ribosomal RNA, complete sequence	2615	2615	100%	0.0	99%	NR_074891.1
<input type="checkbox"/> Shigella boydii strain P288 16S ribosomal RNA gene, partial sequence	2615	2615	100%	0.0	99%	NR_104901.1
<input type="checkbox"/> Escherichia coli str. K-12 substr. MG1655 strain K-12 16S ribosomal RNA, complete sequence	2615	2615	100%	0.0	99%	NR_102804.1
<input type="checkbox"/> Shigella boydii Sb227 strain Sb227 16S ribosomal RNA, complete sequence	2610	2610	100%	0.0	99%	NR_074893.1
<input type="checkbox"/> Shigella sonnei Ss046 strain Ss046 16S ribosomal RNA, complete sequence	2604	2604	100%	0.0	99%	NR_074894.1
<input type="checkbox"/> Shigella dysenteriae Sd197 strain Sd197 16S ribosomal RNA, complete sequence	2588	2588	100%	0.0	99%	NR_074892.1
<input type="checkbox"/> Escherichia albertii strain Albert 19982 16S ribosomal RNA gene, partial sequence	2588	2588	100%	0.0	99%	NR_025569.1
<input type="checkbox"/> Shigella dysenteriae strain ATCC 13313 16S ribosomal RNA gene, partial sequence	2580	2580	100%	0.0	99%	NR_026332.1
<input type="checkbox"/> Escherichia coli strain U 5/41 16S ribosomal RNA gene, partial sequence	2567	2567	98%	0.0	99%	NR_024570.1

Figure 8. BLAST search results with a 1443bp 16s rDNA sequence from Sample#1. Sequences with the highest homology to the query sequence are 16s rDNA from pathogenic bacteria (*Escherichia fergusonii*; *Escherichia Albertii*; *Escherichia coli O157* and *Shigella* species) and non-pathogenic bacteria (*Escherichia coli NBRC 102203*, *Escherichia coli* strain K12 and *Escherichia coli* Strain U5/41). Since the sample is non-pathogenic, the sample was identified as an *E. coli* species.

Identification of a Wild Yeast 2 contaminant in a beverage sample. The microbial contaminant was initially typed as *Candida krusei* (Wild Yeast 1) by culture method by a commercial vendor. The 18s rDNA was sequenced using DTCS Quick Start

Kit and GeXP System. After quality based trimming, the 18s rDNA sequence (Figure 9) was submitted to NCBI for BLAST search. The contaminant was identified as *Candida tropicalis* (Wild Yeast 2).

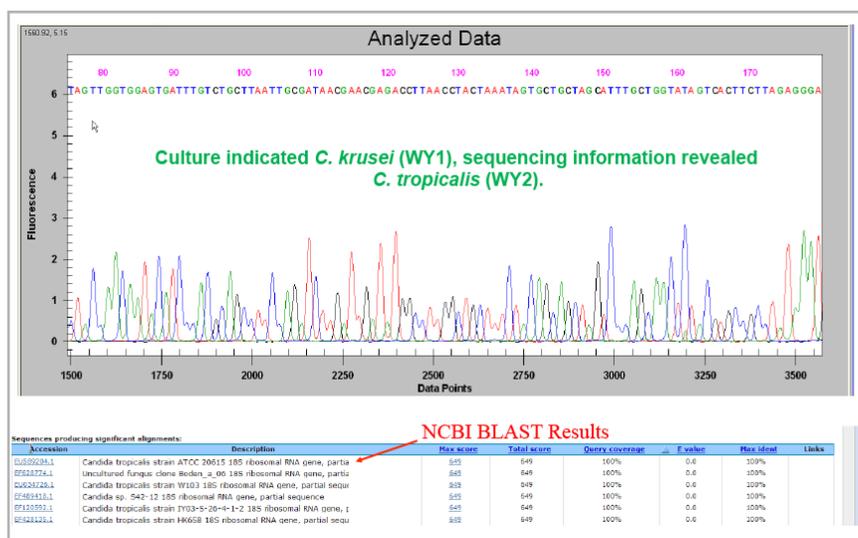


Figure 9. The 18s rDNA sequence of the beverage contaminant and its alignment results in NCBI.

Conclusion

- Sanger Sequencing is more specific and reliable than traditional microbial identification methods such as culture typing.
- Analysis of microbial ribosomal RNA gene sequence on GenomeLab GeXP Genetic Analysis System is simple, fast and accurate.

AB Sciex is doing business as SCIEX.

© 2016 AB Sciex. For research use only. Not for use in diagnostic procedures. The trademarks mentioned herein are the property of the AB Sciex Pte. Ltd. or their respective owners. AB SCIEX™ is being used under license.

RUO-MKT-02-3908 04/2016