

Site-specific differentiation of hydroxyproline isomers using electron activated dissociation (EAD)

Rashmi Madda,¹ Shuai Wu,² John Hui,² Haichuan Liu¹ and Zoe Zhang¹

¹SCIEIX, USA; ²Amgen Inc., USA

This technical note highlights the power of electron activated dissociation (EAD) for clear differentiation of 3- and 4-hydroxyproline (Hyp) isomers to meet the analytical need for biotherapeutic characterization in a case study. The elucidation of Hyp isomers poses an analytical challenge to traditional chromatographic and tandem mass spectrometry (MS/MS) approaches. By contrast, EAD generates signature z -27 and z -43 fragments for 3- and 4-Hyp isomers, respectively, for the unambiguous differentiation of these 2 Hyp isomers.

Natural linker peptides containing proline are commonly used in biotherapeutics.¹ These peptides can undergo structural changes, such as Hyp formation and isomerization, during the manufacturing process.^{2,3} Therefore, there is a need to identify and differentiate Hyp isomers in the development of biotherapeutics to ensure their quality. However, the differentiation of Hyp isomers is challenging using traditional collision-based MS/MS approaches.³ Although conventional electron-based MS/MS techniques can generate signature fragments for the differentiation of Hyp isomers,³ these methods are not suitable for routine biotherapeutic characterization due to slow scan rate and low sensitivity. EAD has proven to be an indispensable tool for rapid differentiation of amino acid isomers, primarily when baseline chromatographic separation cannot be achieved.⁴⁻⁶

In this technical note, optimized EAD workflows were employed to differentiate 3- and 4-Hyp isomers of proline-containing peptides using the ZenoTOF 7600 system equipped with the EAD capability. The detection of diagnostic fragments z -27 for 3-Hyp and z -43 for 4-Hyp enabled confident elucidation of synthetic Hyp-containing peptides and a Hyp isomer identified in real-world samples from the biotherapeutic industry.

Key features of EAD for isomer differentiation and biotherapeutic characterization

- **Unambiguous differentiation of isomers:** EAD generates signature fragments to distinguish amino acid isomers confidently.
- **High-quality MS/MS data:** Gain a 5- to 10-fold increase in MS/MS sensitivity or confident isomer differentiation.
- **Accurate localization:** EAD preserves labile post-translational modifications (PTMs) for their accurate localization.
- **Fast and flexible:** EAD operates in fast and flexible modes of data-dependent acquisition (DDA) or MRM^{HR} with a rapid scanning rate (~20 Hz in DDA mode) and the ability to tune the electron kinetic energy (KE).
- **Single-injection method:** The streamlined EAD DDA workflow provides comprehensive characterization of biotherapeutics in a single injection.

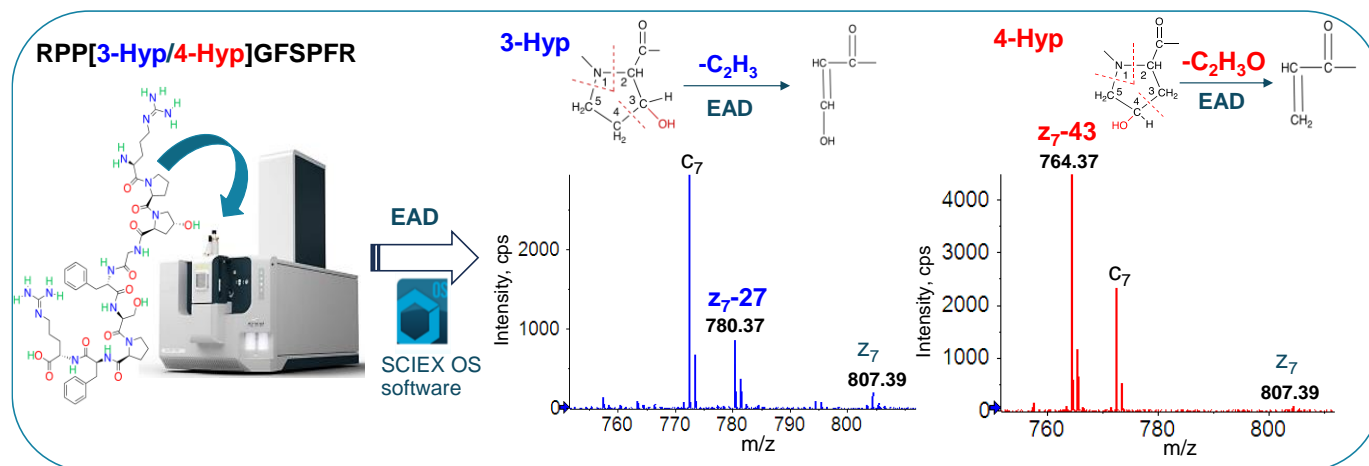


Figure 1. Site-specific differentiation of Hyp isomers using EAD with the ZenoTOF 7600 system. EAD offered by the ZenoTOF 7600 system provides unambiguous differentiation of Hyp isomers by generating signature fragments z -27 (loss of C_2H_3) and z -43 (loss of C_2H_3O) for 3- and 4-Hyp isomers, respectively.

Introduction

The isomerization of amino acid residues, such as aspartic acid (Asp) vs. isoaspartic acid (isoAsp), leucine (Leu) vs. isoleucine (Ile) and 3-Hyp vs. 4-Hyp, can significantly impact the structure and function of biotherapeutics.⁷⁻⁹ Therefore, the differentiation of amino acid isomers plays a role in achieving full characterization of a therapeutic molecule at various stages of its development. It is challenging to differentiate amino acid isomers using conventional collision-based MS/MS. In this study, streamlined EAD workflows were employed to differentiate synthetic 3- and 4-Hyp isomers and determine a Hyp isomer identified in tryptic digests of a reagent protein.

Methods

Samples preparation: Two synthetic bradykinin peptides RPXGFSPFR (X = 3-Hyp or 4-Hyp) were obtained from AnaSpec Inc. The G4P linker peptides containing 3- or 4-Hyp were synthesized and purified at an Amgen facility. Synthetic peptide samples were reconstituted to a final concentration of 10 μ M. The digest of a reagent protein was prepared using trypsin. 1-5 μ l of each sample was injected for LC-MS analysis.

Chromatography: Synthetic bradykinin and G4P linker peptides were separated using the LC gradients 1A and 1B, respectively (Table 1), while the reagent protein digest was analyzed with the LC gradient 2 (Table 1). Peptide separation was performed using a Waters ACQUITY CSH C18 column (2.1 mm x 50 mm, 1.7 μ m, 130 Å, Waters). The flow rate was set to 0.25 mL/min for synthetic Hyp-containing peptides and 0.45 mL/min for the reagent protein digest. The column was kept at 60°C in the column oven of an

Table 1. LC gradients for the separation of Hyp peptides.

Gradient 1A (1B)			Gradient 2		
Time (min)	Mobile phase A (%)	Mobile phase B (%)	Time (min)	Mobile phase A (%)	Mobile phase B (%)
Initial	92	8 (2)	Initial	99	1
1	92	8 (2)	5	99	1
8	75	25 (35)	45	70	30
9	10	90	50	55	45
11	10	90	53	30	70
12	98	2	53.1	30	70
15	98	2	58	30	70
			65	99	1
			70	99	1

ExionLC AD system (SCIEX). The mobile phases A and B consisted of 0.1% formic acid (FA) in water and 0.1% FA in acetonitrile, respectively.

Mass spectrometry: EAD MRM^{HR} data of synthetic Hyp peptides were acquired in SCIEX OS software using the ZenoTOF 7600 system. An EAD data-dependent acquisition (DDA) method was employed to analyze the tryptic digest of the reagent protein. The key TOF MS and MS/MS settings are listed in Tables 2 and 3, respectively.

Table 2. TOF MS parameters.

Parameter	Value
Spray voltage	5500 V
TOF Start mass	300 m/z
TOF Stop mass	1500 m/z
Accumulation time	0.2 s
Source temperature	400°C
Declustering potential	80 V
Collision energy	10 V
Time bins to sum	8

Table 3. EAD MS/MS parameters.

Parameter	MRM ^{HR}	DDA
Q1 resolution		Unit
Electron beam current		5000 nA
Zeno trap		ON
Zeno threshold		100,000 cps
Start mass		100 m/z
Stop mass	3000 m/z	2000 m/z
Accumulation time	1 s	90 ms
Electron KE		7 eV
ETC	100%	Dynamic
EAD RF	100 Da	120 Da
EAD reaction time	10 ms	20 ms
Time bins to sum	10	8

Data processing: The EAD MRM^{HR} and DDA data were analyzed using the Explorer function within SCIEX OS software.

Differentiation of Hyp isomers by EAD

Hydroxylation of a proline residue in CHO-expressed proteins can lead to the formation of 3- and 4-Hyp isomers. The differentiation of these 2 Hyp isomers provides insight into the effect of this modification on the therapeutic molecule. Traditional collision-based MS/MS approaches do not generate diagnostic fragments to differentiate 3- and 4-Hyp isomers.³ By contrast, EAD provides a clear differentiation of Hyp isomers by producing signature z -27 and z -43 fragments for 3- and 4-Hyp, respectively.

In this work, EAD was first employed to characterize 2 synthetic Hyp-containing peptides RPXGSPSPFR (X = 3- or 4-Hyp). These 2 peptides were eluting at similar retention times (RTs) using the gradient 1A shown in Table 1 (Figure 2A and 2D). EAD led to rich MS/MS spectra containing Hyp-signature fragments for confident sequence confirmation and unambiguous differentiation of Hyp isomers (Figure 2B and 2E). The zoomed-in views of EAD spectra showed the diagnostic ions of z_7 -27 for 3-Hyp (Figure 2C) and z_7 -43 for 4-Hyp (Figure 2F). The neutral losses of 27 Da and 43 Da correspond to the cleavage of C_2H_3 and C_2H_3O from the pyrrolidine rings of 3- and 4-Hyp, respectively (Figure 1). The detection of these diagnostic fragments enabled confident differentiation of 2 Hyp isomers.

Differentiation of Hyp isomers of the G4P linker peptides by EAD

The proline residue in the linker peptide—GGGPP—might be hydroxylated and converted into 3- or 4-Hyp during the manufacturing process. To further highlight the power of EAD for Hyp isomer differentiation, 2 G4P linker peptides (SLSLSPGGGGGXAR where X = 3- or 4-Hyp) were synthesized and characterized by EAD. Similar to the example described above (Figure 2), the 2 Hyp-containing G4P linker peptides were eluting at close RTs (Figure 3A and 3D). Excellent EAD fragmentation of 2 isomers led to the detection of a nearly complete series of c/z fragments (Figure 3B and 3E). The expanded views of EAD spectra showed the detection of a z_3 -27 fragment at m/z 316.16 for 3-Hyp (Figure 3C) and a z_3 -43 ion at m/z 300.17 for 4-Hyp (Figure 3F). These results further demonstrate the power of EAD for precise and rapid differentiation of 3- and 4-Hyp isomers of the proline-containing linker peptides.

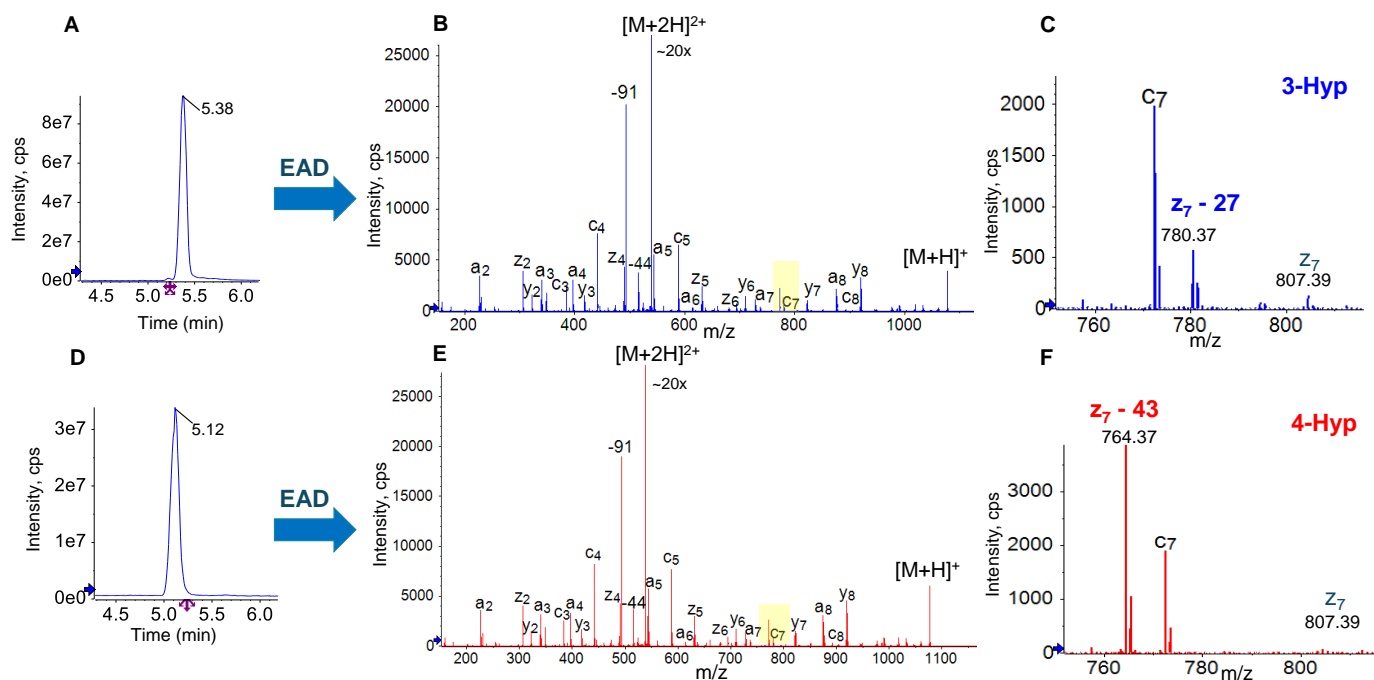


Figure 2. Differentiation of bradykinin peptides RPXGSPSPFR (X = 3-Hyp or 4-Hyp) by EAD. The 2 Hyp isomers were eluting at similar RTs of 5.38 min (A) and 5.12 min (D) under the chromatographic condition employed in this work. EAD led to similar high-quality MS/MS spectra of the 3-Hyp (B) and 4-Hyp isomers (E). The zoomed-in views of m/z ~750-810 Da showed the diagnostic ions of z_7 -27 (m/z 780.37) and z_7 -43 (m/z 764.37) generated from the 3-Hyp (C) and 4-Hyp (F) peptides, respectively. The peaks labeled with -44 and -91 in B and E are the neutral-loss products from the side chain cleavages of the Arg and Phe residues, respectively. In B and E, not all fragments were labeled for spectral clarity.

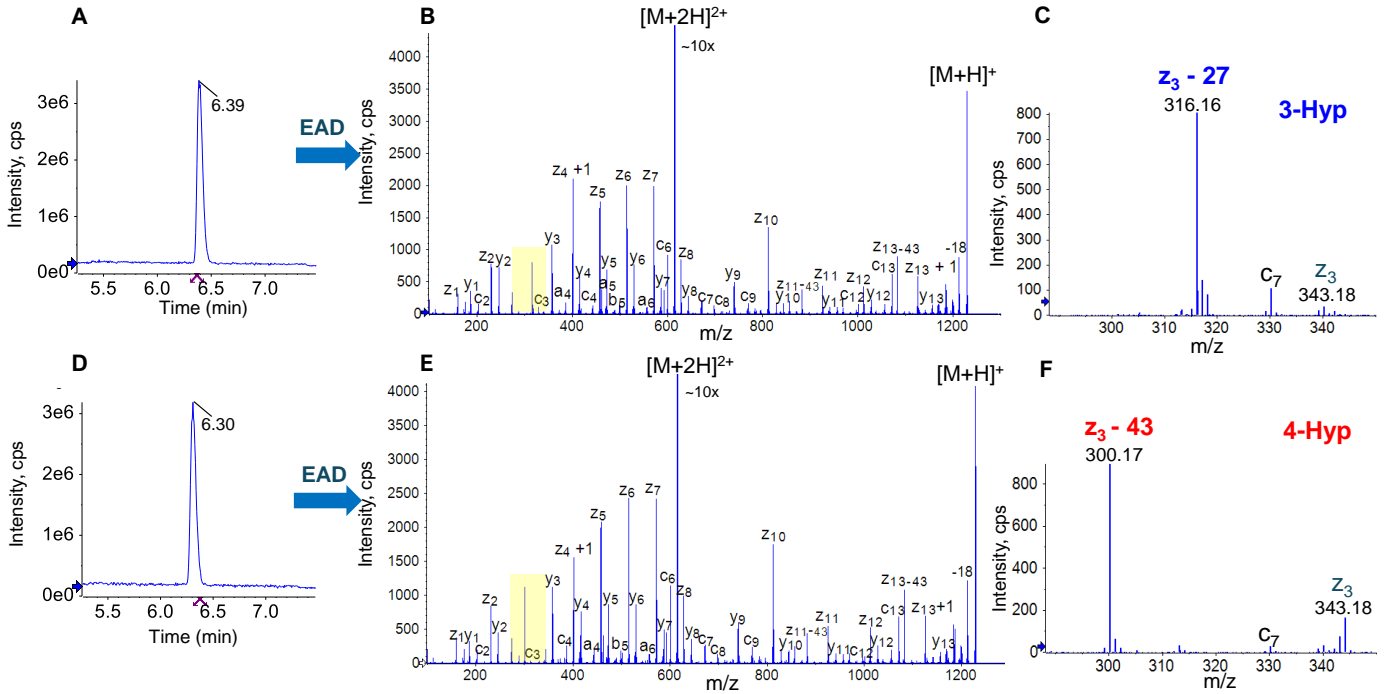


Figure 3. Differentiation of Hyp-modified G4P linker peptides SLSLSPGGGGXAR (X = 3- or 4-Hyp) using EAD. 2 Hyp-containing linker peptides were eluting at similar RTs of 6.39 min (A) and 6.31 min (D). EAD spectra of 2 doubly charged precursors were featured with rich sequence ions for confident sequence confirmation (B and E). The fragments z_{11-43} and z_{13-43} were the signature fragments of 2 Leu residues near the N-terminus (B and E). The detection of diagnostic ions z_3-27 at m/z 316.16 (C) and z_3-43 at m/z 300.17 (F) enabled the confident differentiation of the 3- and 4-Hyp isomers of this linker peptide. Not all fragments were labeled in B and E for spectral clarity.

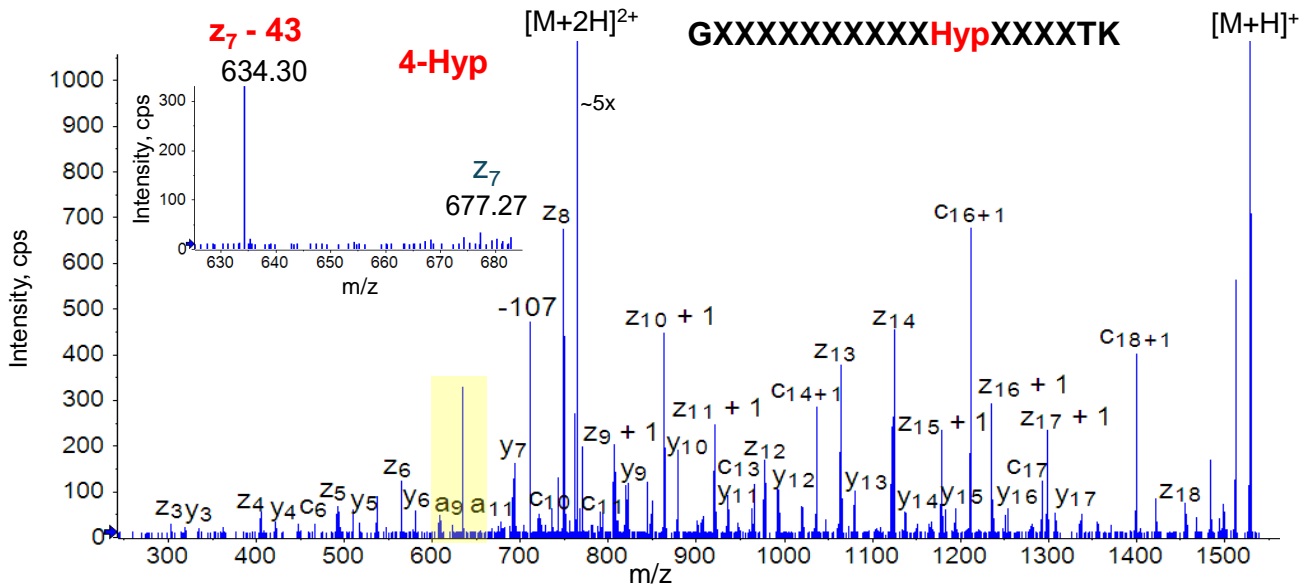


Figure 4. Identification and elucidation of a Hyp-containing peptide GXXXXXXXXXXHypXXXXTK from a reagent protein using an EAD DDA method. EAD produced rich sequence fragments of the Hyp-containing peptide from the reagent protein for its confident identification. The detection of signature z_7-43 confirmed the presence of 4-Hyp in this peptide (inset). The amino acid residues marked "X" were covered for proprietary reasons. The loss of 107 from the precursor corresponds to the side chain cleavage of a Trp residue.

Case Study: Elucidation of a Hyp isomer identified in a real-world reagent protein sample

E. coli expressed reagent proteins have been approved as an oncology therapeutic. However, this version of reagent protein has a short *in vivo* half-life and toxic side effects when administered at high doses. Hence, there is considerable interest in developing CHO-expressed reagent protein. Recent LC-MS analyses of real-world tryptic digests from several CHO-expressed proteins led to the unexpected identification of a Hyp-containing peptide (data not shown). EAD was leveraged to determine whether hydroxylation of the proline residue occurred in the 3rd or 4th position of the pyrrolidine ring.

Figure 4 shows the EAD spectrum of the doubly charged peptide GXXXXXXXXXXHypXXXXTK from the reagent protein (amino acid residues marked "X" were covered for proprietary reasons). EAD generated rich sequence fragments for confident identification of this Hyp-containing peptide in the reagent protein digest. The detection of a diagnostic ion $z-43$ (loss of C_2H_3O) at m/z 634.3 confirmed the presence of 4-Hyp instead of 3-Hyp in this reagent protein peptide (inset in Figure 4).

In summary, the data presented in this technical note demonstrates the power of EAD for confident and rapid elucidation of Hyp-containing peptides that are challenging to characterize using traditional chromatographic or fragmentation techniques. The ability of EAD for confident isomer differentiation was successfully leveraged to determine a Hyp isomer identified in real-world samples from the biotherapeutic industry.

Conclusions

- EAD provided confident identification and clear differentiation of 3- and 4-Hyp isomers that are challenging to elucidate using traditional fragmentation techniques.
- EAD produced the signature $z-27$ and $z-43$ fragments for 3- and 4-Hyp isomers, respectively, enabling unambiguous differentiation of these 2 isomers.
- Streamlined EAD methods were employed to differentiate synthetic peptides containing 3- and 4-Hyp residues confidently.
- An unexpected Hyp-containing peptide identified in the tryptic digests of a reagent protein was the 4-Hyp isomer based on the detection of a signature $z-43$ fragment using EAD.
- EAD workflows described in this technical note can be leveraged to identify and differentiate Hyp-containing peptides from protein therapeutics rapidly.

References

1. Arnik Shah *et al.* (2024) Characterization of charge variants, including post-translational modifications and proteoforms, of bispecific antigen-binding protein by cation-exchange chromatography coupled to native mass spectrometry. [Talanta. 266\(1\):125062.](#)
2. Oksana Tyshchuk *et al.* (2019) Characterization and prediction of positional 4-hydroxyproline and sulfotyrosine, two post-translational modifications that can occur at substantial levels in CHO cells-expressed biotherapeutics. [MAbs.11\(7\):1219-1232.](#)
3. Fengfei Ma *et al.* (2018) Large-scale differentiation and site specific discrimination of hydroxyproline isomers by electron transfer/higher-energy collision dissociation (ETHcD) mass spectrometry. [Anal. Chem. 90\(9\):5857-5864.](#)
4. Takashi Baba *et al.* (2021) Dissociation of Biomolecules by an Intense Low-Energy Electron Beam in a High Sensitivity Time-of-Flight Mass Spectrometer. [J. Am. Soc. Mass Spectrom. 32\(8\):1964-1975.](#)
5. Comprehensive differentiation isomers from forced degradation by electron activation dissociation (EAD). SCIEX technical note, [RUO-MKT-02-14730-A.](#)
6. An evaluation of single injection platform method for advanced characterization of protein therapeutics using electron activation dissociation (EAD). [SCIEX technical note, RUO-MKT-02-13965-A.](#)
7. Mary Krause and Erinc Sahin. (2019) Chemical and physical instabilities in manufacturing and storage of therapeutic proteins. [Current Opinion in Biotechnology. 60:159-167.](#)
8. Anshu Kuriakose *et al.* (2016) Immunogenicity of Biotherapeutics: Causes and Association with Posttranslational Modifications. [J. Immunol. Res. 2016:1298473.](#)
9. Anna Robotham and John Kelly (2020) LC-MS characterization of antibody-based therapeutics: recent highlights and future prospects. [Approaches to the Purification, Analysis and Characterization of Antibody-Based Therapeutics.](#) Chapter 1:1-33.

The SCIEX clinical diagnostic portfolio is For In Vitro Diagnostic Use. Rx Only. Product(s) not available in all countries. For information on availability, please contact your local sales representative or refer to <https://sciex.com/diagnostics>. All other products are For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd. or their respective owners in the United States and/or certain other countries (see www.sciex.com/trademarks).

© 2023 DH Tech. Dev. Pte. Ltd. MKT-30610-A



Headquarters

500 Old Connecticut Path | Framingham, MA 01701
USA
Phone 508-383-7700
sciex.com

International Sales

For our office locations please all the division
headquarters or refer to our website at
sciex.com/offices