

Comprehensive mapping of disulfide linkages of adalimumab using electron activated dissociation (EAD) and collision induced dissociation (CID)

Mona Hamada¹, Zoe Zhang² and Amandine Boudreau¹

¹SCIEX, Canada; ²SCIEX, USA

This technical note describes a comprehensive mapping of expected and scrambled disulfide linkages and free thiols in adalimumab using EAD and CID. Data were acquired on the ZenoTOF 7600 system with SCIEX OS software, and automated data interpretation was performed using Biologics Explorer software.

Disulfide linkages are recognized by regulatory agencies such as the FDA, EMA and ICH as critical quality attributes (CQAs) of protein biotherapeutics.^{1,2} Reduction and scrambling of the disulfide linkages can occur during manufacturing and storage, disrupting the proper folding of biotherapeutics and consequently impacting their functions.^{1,2} Therefore, accurate mapping of disulfide linkages and quantitative assessment of their variants are crucial to ensure the safety, efficacy, and stability of the drug product.

CID is the most widely adopted fragmentation mode in LC-MS bottom-up peptide mapping analysis.³ However, CID may not fragment disulfide linkages, and their presence can further hinder effective peptide backbone fragmentation, resulting in limited sequence information.^{2,3} In contrast, EAD effectively cleaves the disulfide linkages, further empowering its efficient peptide backbone fragmentation.^{2,3} EAD is an exceptionally powerful tool

that allows the comprehensive mapping of the cysteine residues and the peptides involved in these complex linkages.³

The data presented in this technical note shows the in-depth characterization of the complex disulfide linkage structure of adalimumab, including the identification of impurities such as free thiols and scrambled disulfide peptides.

Key features of disulfide linkage analysis using Biologics Explorer software

- **Confident identification:** High-quality MS/MS data obtained by EAD and CID fragmentation allows solid sequence confirmation and characterization of expected and scrambled disulfide linkages and free thiols.
- **Accurate localization:** Post-translational modifications (PTM), including amino acid isomerization, can be unambiguously localized on disulfide-linked peptides in the same run.
- **Powerful, streamlined solution:** EAD combined with superior data processing algorithms in Biologics Explorer software provides a complete solution for challenging disulfide linkages analysis.

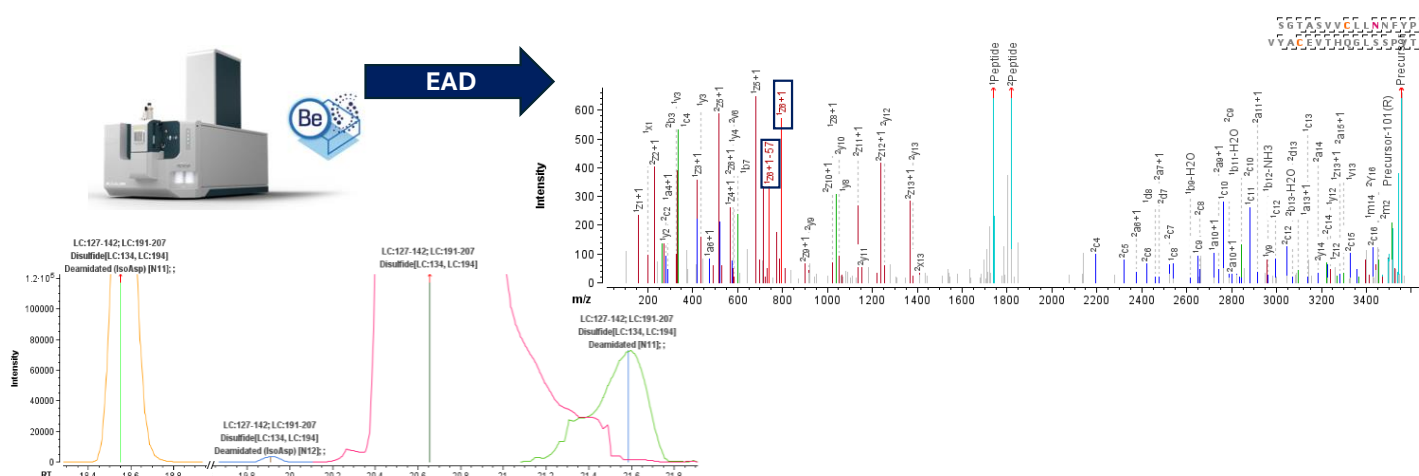


Figure 1. Characterization of disulfide linkages and PTMs on the same peptide using EAD with the ZenoTOF 7600 system. An intrachain disulfide-linked peptide from the light chain was chromatographically resolved from its deamidated forms (left panel). The highly sensitive EAD method offered by the ZenoTOF 7600 system led to rich fragmentation spectra of the disulfide-linked peptides (right panel). The detection of diagnostic fragments produced by EAD confirms the isoAsp formation at positions N11 (Asn137) and N12 (Asn138) following deamidation.

Introduction

Adalimumab is a recombinant human IgG1 monoclonal antibody that prevents the interaction of tumor necrosis factor (TNF) with TNF receptors. IgG1 antibodies are stabilized by 16 disulfide linkages (12 intrachain and 4 interchain) formed between all 32 Cys residues.^{1,4} Both intra- and inter-chain disulfide bonds are critical for establishing the tertiary and quaternary structure of biologics.⁴ Normally, Cys residues pair with their correct partner residues. However, under stress conditions such as heat, radicals, improper pH or agitation, unexpected linkages can be formed via free thiols, resulting in potential detrimental effects on the physicochemical and biological properties of biologics.¹

Methods

Sample preparation: Adalimumab was denaturated with 7.2M guanidine hydrochloride in 50mM Tris buffer (pH 7.0). Free cysteine was capped with 5mM iodoacetamide. Protein digestion was performed with trypsin/Lys-C enzyme at 30°C for 16 hours at pH 7. The reaction was stopped with 1% formic acid.

Chromatography: 4 µg (CID) or 8 µg (EAD) of the trypsin/Lys-C digest were separated with a CSH C18 column (2.1 × 100 mm, 1.7 µm, 130 Å, Waters) using an ExionLC AD system. The mobile phase A consisted of water with 0.1% formic acid, while the mobile phase B was 0.1% formic acid in acetonitrile. A gradient profile was used at a flow rate of 300 µL/min (Table 1). The column temperature was maintained at 50°C.

Table 1. LC gradient for peptide separation.

Time (min)	Mobile phase A (%)	Mobile phase B (%)
Initial	99	1
5	99	1
6	90	10
25	75	25
65	60	40
70	40	60
70.5	10	90
74	10	90
74.1	99	1
75	99	1

Mass spectrometry: Data were acquired on the ZenoTOF 7600 system in positive ionization mode with a DDA method using SCIEX OS software. EAD and CID MS/MS data were acquired in separate injections. Detailed method parameters are described in Table 2.

Table 2. DDA parameters using CID and EAD.

Parameter	MS	MS/MS
Scan mode	TOF-MS	IDA dependent
Gas 1		50
Gas 2		50
Curtain gas		35
Source temperature °C		450
Ion spray voltage		5500
Declustering potential		80
CAD gas		7
Charge state		2-10
Q1 resolution		Low
Time bins to sum		8
Start mass	200 m/z	100
Stop mass	2000 m/z	3000
Zeno trap	N/A*	ON
Zeno threshold	N/A*	100,000 cps
Maximum candidate ions	10	N/A*
Intensity threshold exceeds	125	N/A*
Accumulation time	0.25 s	0.09 s
CE	12 V	12 (EAD) Dynamic (CID)
Electron beam current (EAD)	N/A*	5500 nA
Electron KE (EAD)	N/A*	7 eV
Reaction time (EAD)	N/A*	10 ms
EAD RF	N/A*	100 Da

*N/A: Not applicable

Data processing: EAD and CID DDA data were analyzed in Biologics Explorer software version 5.0 using the *Pepmap_Extended_Be5.0* workflow. The following fixed disulfide linkages were specified in the first peptide mapping activity node: [HC:22, HC:96], [HC:148, HC:204], [HC:224, LC:214], [HC:230, HC:230], [HC:233, HC:233], [HC:265, HC:325], [HC:371, HC:429], [LC:23, LC:88] and [LC:134, LC:194]. The second peptide mapping activity node was set for the de-novo mapping of unexpected disulfide linkages. For all other parameters, the default workflow settings were used.

Identification of expected disulfide linkages

The ZenoTOF 7600 system introduces an alternative fragmentation mechanism, EAD. In addition, the Zeno trap improves the MS/MS sensitivity up to 7-fold and further enhances the detection of fragment ions generated by EAD or CID. Consequently, rich MS/MS spectra are generated for the analysis of disulfide-linked peptides, augmenting data confidence. Combined with automated data processing using Biologics Explorer software, a streamlined solution is provided for complex disulfide linkages analysis.

The characterization results obtained from EAD and CID were compared. Both techniques successfully identified all expected disulfide linkages in adalimumab with a mass error below 8 ppm. Overall, EAD demonstrated efficient disulfide linkage cleavage during the fragmentation, leading to a higher coverage of the peptide backbone in the MS/MS spectra, especially for the longer peptides.

Figure 2 shows an example of EAD and CID MS/MS spectra for peptide $\text{SLRLSCAASGFTFDDYAMHWVR}=\text{AEDTAVYYCAK}$ ($z=+5$, 753.55 m/z). The peptide, with a molecular weight of 3,762 Da, contains a disulfide bond as shown. In EAD analysis, the disulfide-linked peptide was identified in 3 different charge states ($z=+4$, $+5$, $+6$) with a high score of 328. Notably, EAD generates dominant fragments resulting from the cleavage of the disulfide bond, leading to efficient fragmentation of the peptide backbones. Consequently, it results in a higher amino acid sequence coverage (73%, 84% and 81% for the 3 charge states of the linked peptide, respectively). On the other hand, using CID, the disulfide-linked peptide was detected in charge state $z=+5$ with a lower score of 270. Due to the inability of CID to fragment the disulfide linkage, a lower amino acid sequence coverage (60%) was observed for the linked peptide.

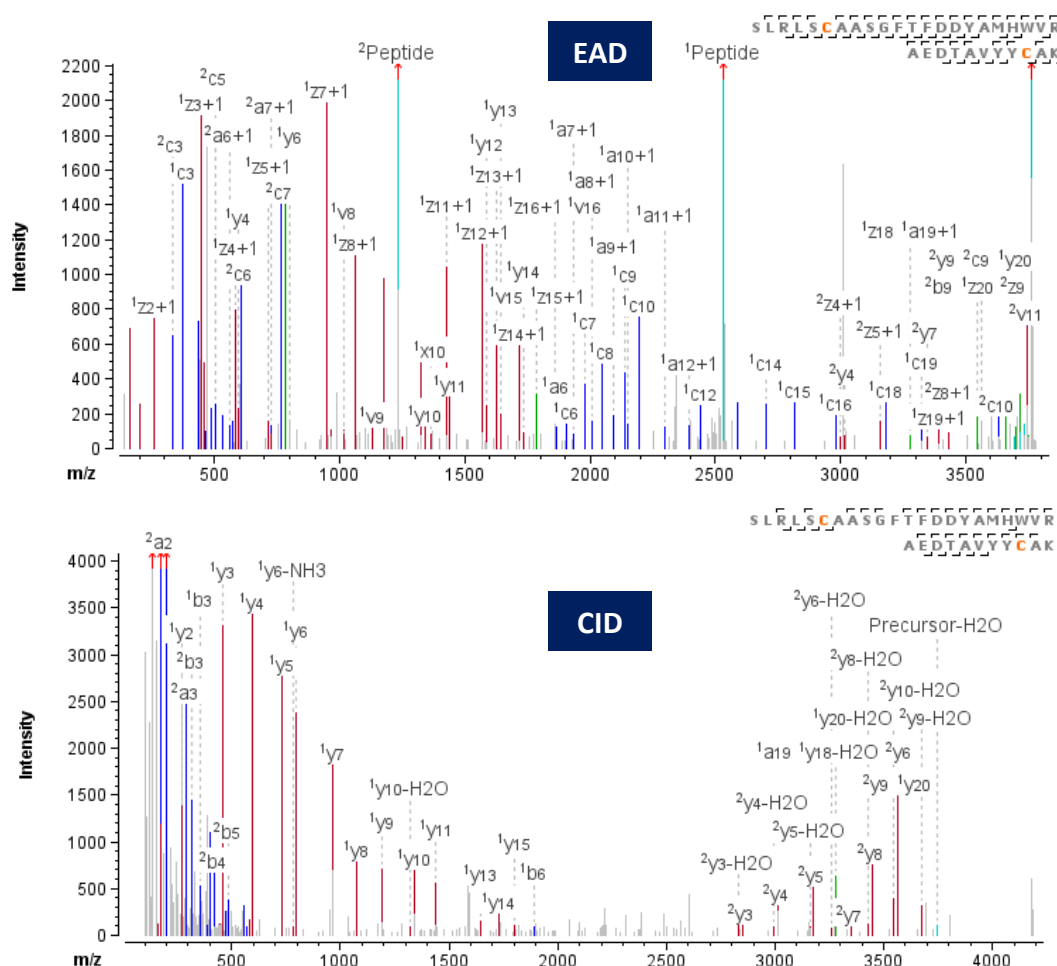


Figure 2. EAD (top) and CID (bottom) MS/MS spectra of the disulfide-linked peptide $\text{SLRLSCAASGFTFDDYAMHWVR}=\text{AEDTAVYYCAK}$, Cys22=Cys96, an intrachain disulfide bond from the heavy chain.

PTM characterization on disulfide-linked peptides

EAD significantly enhances the characterization and localization of post-translational modifications (PTMs) on disulfide-linked peptides, enabling comprehensive characterization of challenging PTMs such as isomerization. Figure 1 shows the chromatographic separation of SGTASVVCLLN^FYPR=VYACEVTHQGLSPVTK, an intrachain disulfide-linked peptide formed via Cys134 and Cys194 of the light chain, along with its deamidated forms. Both Asn137 and Asn138 are possible locations for deamidation. Further, isomerization of Asp to isoAsp can happen after the deamidation.

The localization and differentiation of isomers can be easily confirmed owing to the effective backbone fragmentation by EAD and its ability to generate signature fragment ions from isomers. As shown in Figure 3, the presence of the signature fragment ion pair z^*_6 and z^*_6-57 confirms the deamidation at Asn 137 and the formation of isoAsp, with a mass error of less than 2 ppm.

On the other hand, despite the very low abundance of the peptide observed at 19.9 min (Figure 1), EAD successfully generated the signature fragment ion pair z^*_5 and z^*_5-57 (with a mass error <7 ppm), further confirming the formation of isoAsp at Asn138 (Figure 3).

Identification of complex disulfide linkages

The most complex disulfide linkage of adalimumab lies in the hinge region, where the heavy chain is stabilized by 2 interchain disulfide bonds at Cys230-Cys230 and Cys233-Cys233. As such, non-reduced digest is anticipated to yield peptides connected by double or triple disulfide bonds.


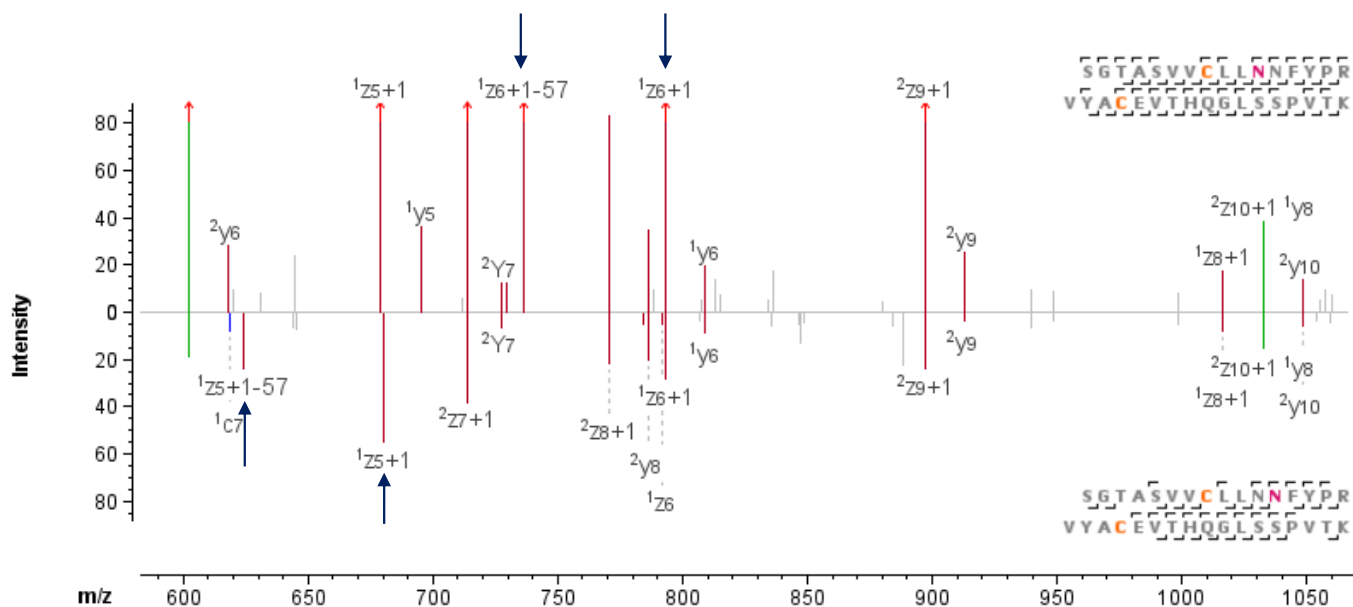


Figure 3. EAD MS/MS spectra of the disulfide linked peptide SGTASVVCLLN^FYPR=VYACEVTHQGLSPVTK, Cys134-Cys194 of the light chain. Asp and isoAsp can be localized and differentiated between Asn137 (top) or Asn138 (bottom) due to the signature fragment ion pair generation by EAD.

Figure 4 displays the EAD and CID MS/MS spectra of 3 peptides linked by 3 disulfide bonds, producing a large peptide complex ($z=+8$, m/z 838.163, mass error=1.72 ppm). Notable differences were observed between the 2 fragmentation techniques for this challenging peptide. EAD exhibited a preference for cleaving the disulfide linkages, generating fragments corresponding to the individual peptides. However, these peptides were absent in the CID spectrum due to its incapacity to fragment the disulfide linkages effectively. In addition, the disulfide linkage between peptide1 (Cys224, HC) and peptide3 (Cys214, LC) was confirmed in the EAD spectrum by 2 fragment ions (1c2 and 1c3, with mass errors of -1.9 and -9.4 ppm, respectively).

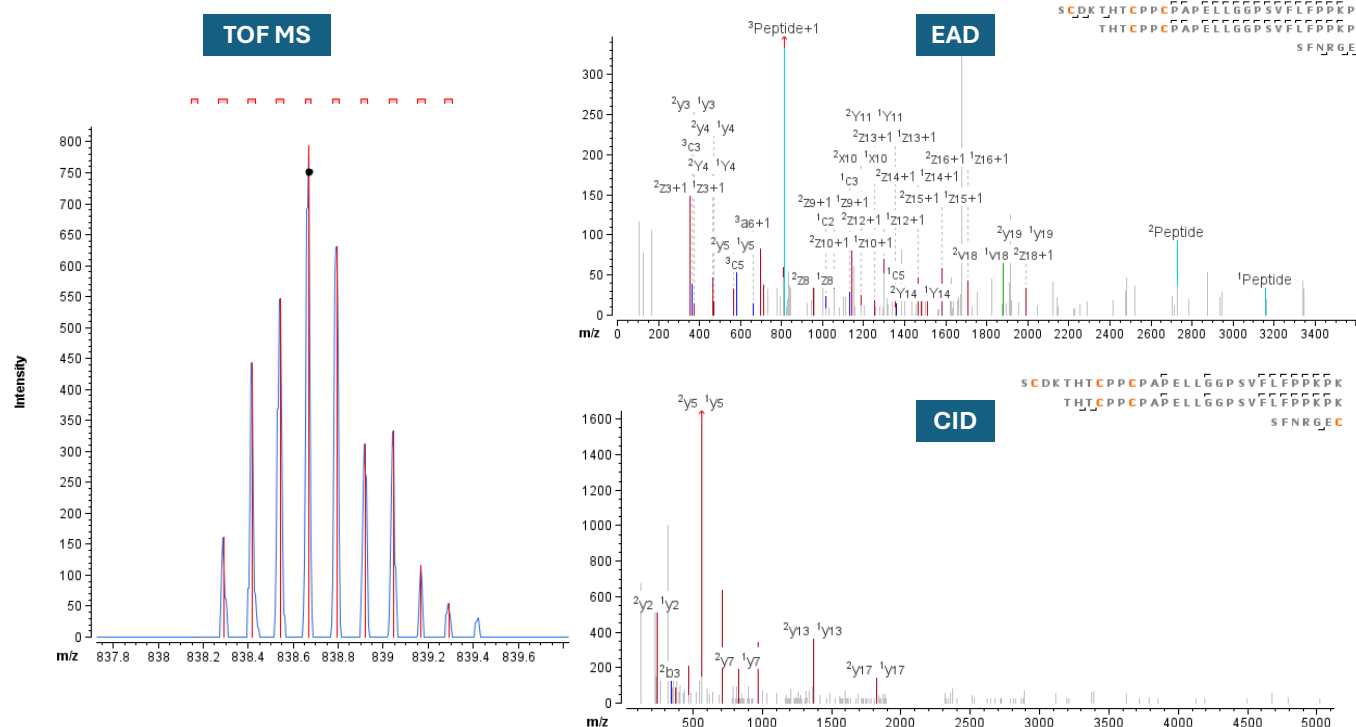

Scrambled disulfide linkages

Qualitative and quantitative assessment of scrambled disulfide linkages is a critical quality attribute of biologics. As such, several studies have investigated the impact of various sample preparation or storage conditions on the degree of scrambling of Cys residues in different IgG biotherapeutics.^{1,2,5} For example, Cys214 stabilizes the higher-order structure of adalimumab via the formation of an inter-chain disulfide linkage with Cys224 of the heavy chain^{4,5}, and heating adalimumab at 70 °C for 30 min resulted in 100% scrambling at Cys214 residue of the light chain⁵. In general, interchain linkages are more susceptible to reduction, incomplete formation and shuffling than intrachain linkages.¹

Figure 5 shows the peptide chromatogram, EAD and CID MS/MS spectra of VYACEVTHQGLSSPVTK=SFNRRGEC (z=+5, 526.45 m/z), where a scrambled disulfide linkage is formed between Cys194 and Cys214 of the light chains. Compared with the expected disulfide linkage at Cys224:HC-Cys214:LC, this scrambled-linked

peptide is of low abundance, representing only 0.4% of disulfide linkages at Cys214 (Figure 5, left). Once again, EAD demonstrated high peptide backbone amino acid sequence coverage of 84% compared to CID coverage of 55%. This was achieved by generating fragment ions corresponding to the individual peptides after breaking the disulfide linkage.

EAD generated the signature fragment for Leu, z⁺7-43 (w₇, mass error=5.6 ppm, Figure 5), which further confirmed the presence of Leu residue and differentiated it from its isomer Ile. This additional level of information obtained through single injection of an EAD method enabled the confirmation of the identity of the connected peptides. This eliminates the possibility of misassignment, especially in the case of low-abundance scrambled disulfide linkages, thus minimizing the instrument time.

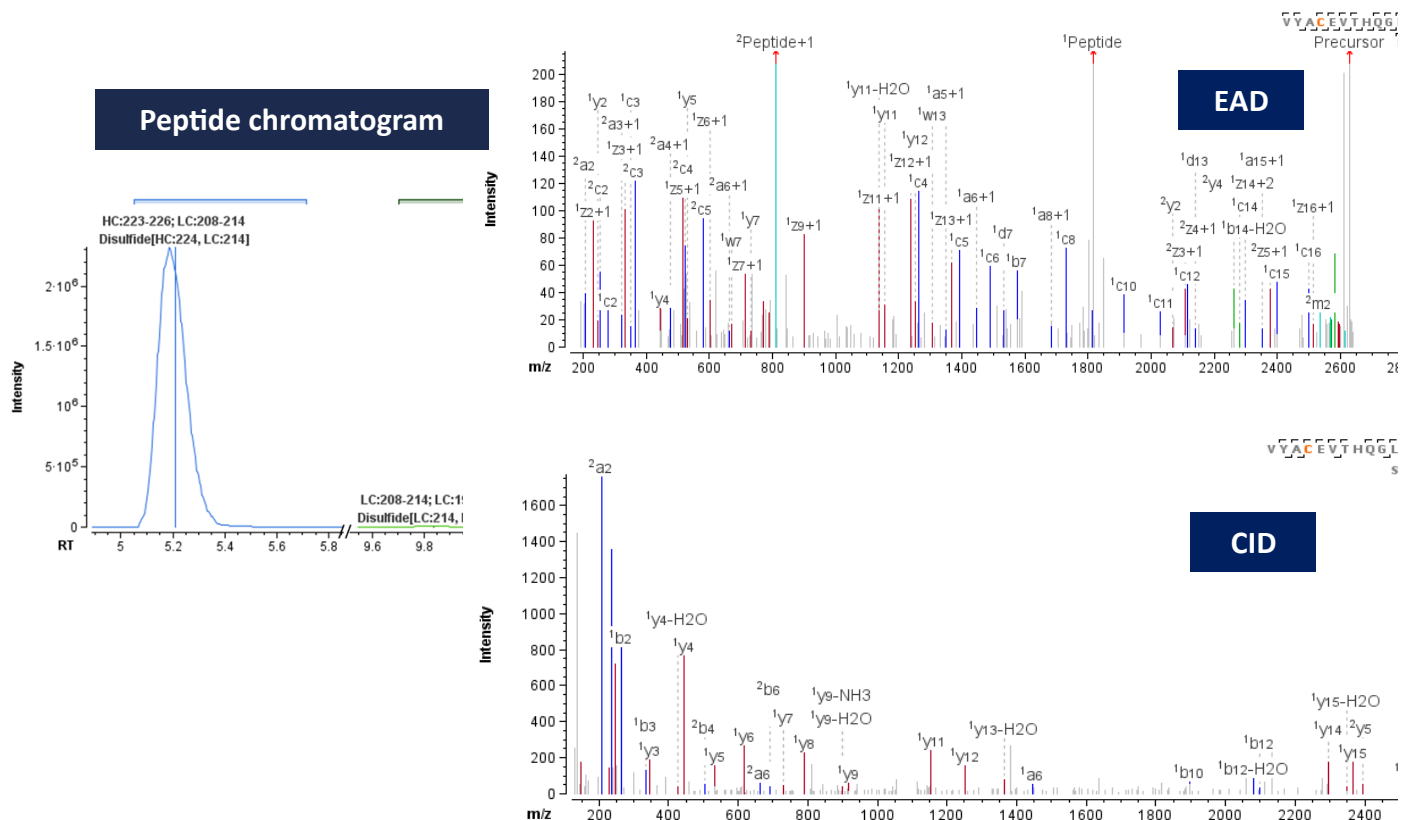


Figure 5. EAD (top right) and CID (bottom right) MS/MS spectra of the disulfide-linked peptide VYACEVTHQGLSSPVTK=SFNRRGEC, Cys194=Cys214, a scrambled disulfide linkage from the light chain that is chromatographically resolved from its native form (left).

Free Cys have been reported as a cause of forming non-native disulfide bonds during manufacturing², which can impact the protein's 3-dimensional structure. Therefore, it is necessary to identify and monitor existing free thiols to avoid potential risks to drug safety. To accurately detect free thiols, all free Cys residues were capped with iodoacetamide before enzymatic digestion, leading to a +57 Da mass increase per alkylation. Low levels of free thiols were detected at different Cys residues of adalimumab. Figure 6 illustrates an example from peptide SGTASVVCLLNNFYPR $z=+3$, where carbamidomethylation was detected at Cys134. Both EAD and CID methods provided highly informative spectra with sequence coverage of 93% and 73%, respectively, thereby confirming the existence of a free thiol. Furthermore, the EAD spectrum offers additional capability to confirm the presence of 2 Leu residues through the detection of z^*_{7-43} (w_7) and z^*_{8-43} (w_8) signature fragments with mass error = 0.14 and -1.52 ppm, respectively.

- Confident sequence and disulfide linkage confirmation of long disulfide-linked peptides were achieved with EAD
- For peptides that contain 2 or more disulfide bonds, EAD generated rich fragments allowing reliable confirmation of individual peptides involved in the disulfides
- EAD spectra demonstrate superior fragmentation coverage on disulfide-linked peptides compared to CID
- Free thiols were detected at low levels with high confidence using both EAD and CID
- EAD produces signature ions to facilitate isomer differentiation, like Asp and isoAsp, Leu and Ile
- Biologics Explorer software facilitates automatic, state-of-the-art data processing for both routine and advanced characterization of complex linkages in biotherapeutics and standard monoclonal antibodies in a reproducible manner

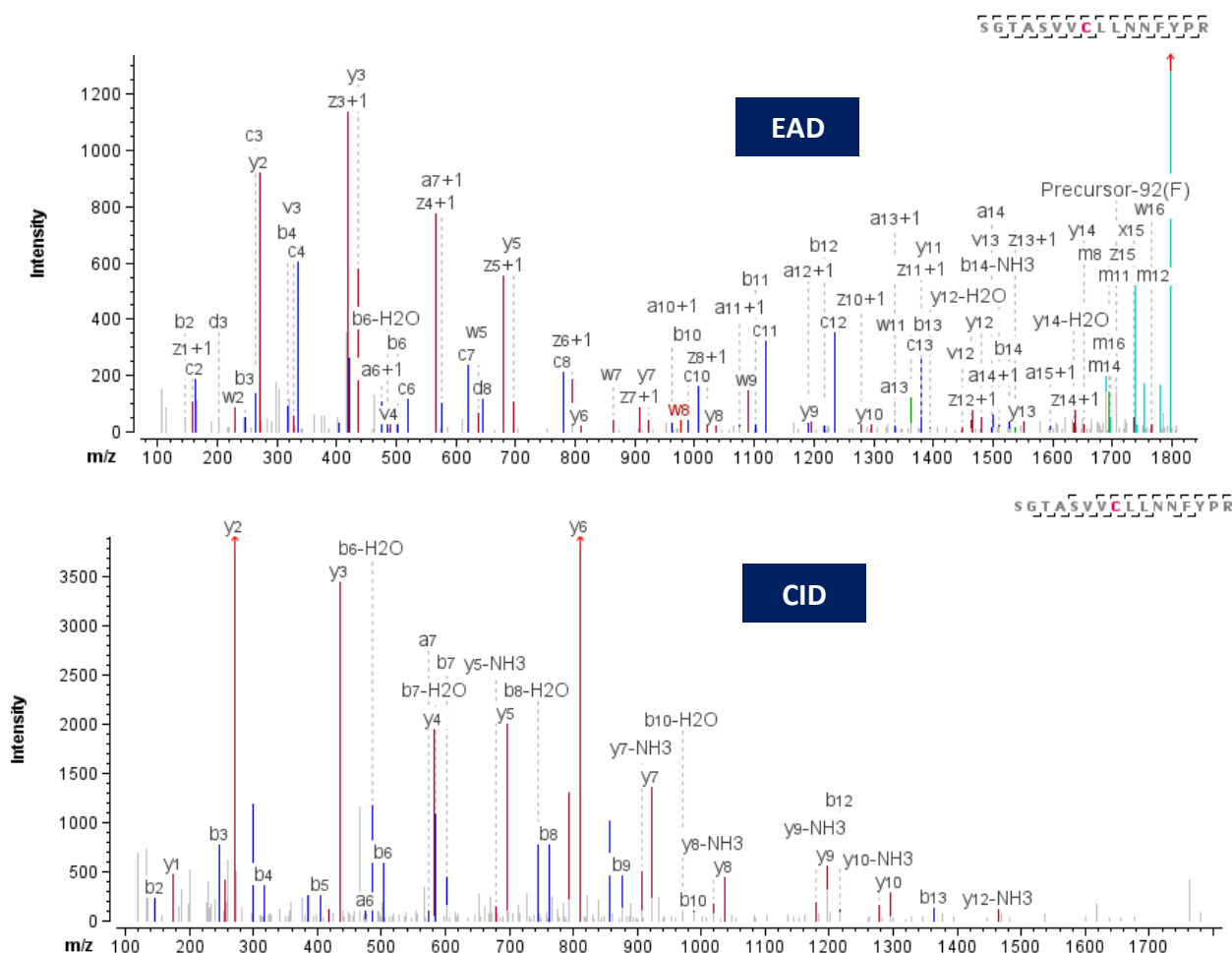


Figure 6. EAD (Top) and CID (bottom) MS/MS spectra of the peptide SGTASVVCLLNNFYPR, where a free thiol was capped using iodoacetamide at Cys 134.

References

1. Coghlan J et al. (2022) Streamlining the characterization of disulfide bond shuffling and protein degradation in IgG1 biopharmaceuticals under native and stressed conditions, *Front. Bioeng. Biotechnol.* 10, 862456.
2. Lakbub J et al. (2018) Recent mass spectrometry-based techniques and considerations for disulfide bond characterization in proteins, *Anal. Bioanal. Chem.* 410, 2467-2484.
3. Huang L et al. (2019) Complete mapping of disulfide linkages for etanercept products by multi-enzyme digestion coupled with LC-MS/MS using multi-fragmentations including CID and ETD, *J Food Drug Anal.* 27.2, 531-541.
4. Kwon J et al. (2021) Physicochemical and biological similarity assessment of LBAL, a biosimilar to adalimumab reference product (Humira®). *Anim. Cells Syst.* 25.3, 182-194.
5. Resemann A et al. (2018) Rapid, automated characterization of disulfide bond scrambling and IgG2 isoform determination, *MAbs.* 10.8,1200-1213.

The SCIEX clinical diagnostic portfolio is For In Vitro Diagnostic Use. Rx Only. Product(s) not available in all countries. For information on availability, please contact your local sales representative or refer to <https://sciex.com/diagnostics>. All other products are For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd. or their respective owners in the United States and/or certain other countries (see www.sciex.com/trademarks).

© 2024 DH Tech. Dev. Pte. Ltd. MKT-32122-A



Headquarters

500 Old Connecticut Path | Framingham, MA 01701
USA
Phone 508-383-7700
sciex.com

International Sales

For our office locations please all the division
headquarters or refer to our website at
sciex.com/offices