



Automated solid-phase extraction system for the analysis of PFAS in drinking water

Craig M. Butt¹, Lily Sanchez², Prem Parmar², Sam Lodge³

¹ SCIEX, USA; ² Orange County Water District, Fountain Valley, CA;

³ Phenomenex, USA

This technical note demonstrates the use of the PromoChrom automated solid-phase extraction (SPE) system for the analysis of PFAS in drinking water following EPA Method 533. Using the SCIEX QTRAP 6500+ system, negligible background contamination was observed as well as excellent method performance in spikes into reagent water and raw untreated groundwater samples. Milli-Q water blank spikes at 2, 40 and 70 ng/L showed mean recovery ranging from 96% to 115% with the mean precision <10% coefficient of variation [CV]. Sample matrix and duplicate spikes at 2 and 40 ng/L showed mean recovery between 95% and 122% with <13%CV [representative chromatograms shown in **Figure 1**]. Analysis of raw untreated groundwater samples showed the detection of 5 out of the 6 EPA-regulated PFAS compounds.

Key benefits of the PromoChrom automated SPE system for the analysis of PFAS in drinking water

- **Automated sample extraction.** Reduced sample preparation time through the fast, reproducible extraction and clean-up of drinking water samples
- **Low system background contamination.** Quality control blanks predominantly showed non-detectable PFAS levels
- **Excellent method performance in quality control spikes.** Method performance was evaluated with spikes into Milli-Q water and raw untreated groundwater samples
- **Real-world water samples.** 10 raw untreated groundwater samples analyzed to demonstrate method applicability

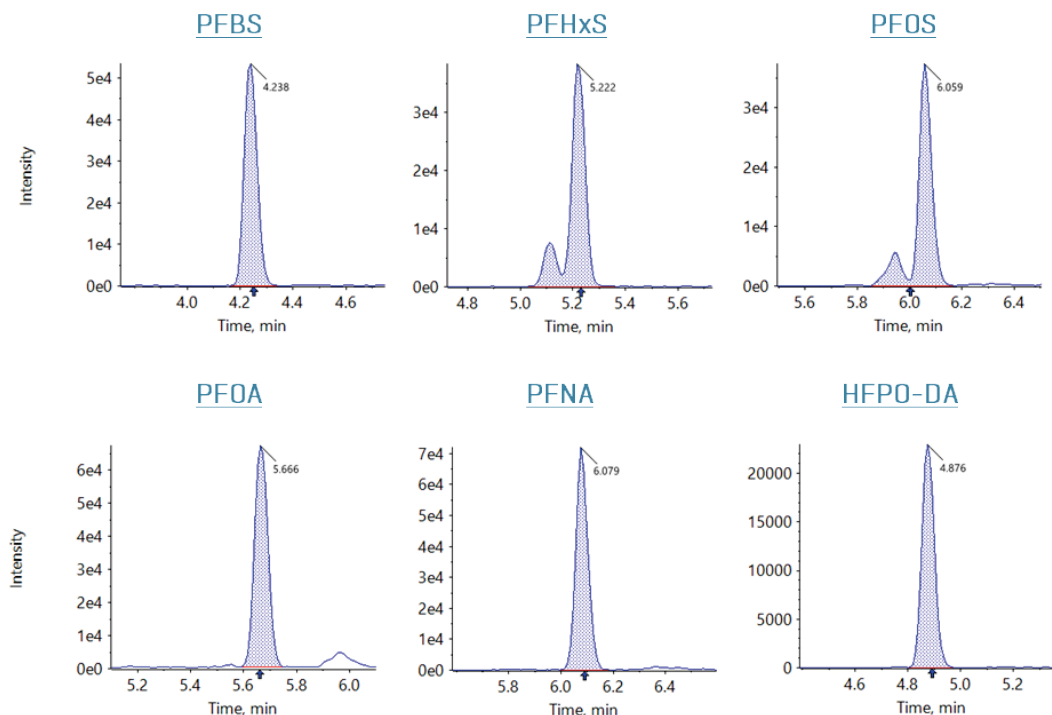


Figure 1. Representative extracted ion chromatograms (XICs) in a 2 ng/L laboratory fortified sample matrix (LFSM) sample for the 6 EPA-regulated PFAS compounds. Raw, untreated groundwater was used for the LFSM sample.

Introduction

The contamination of PFAS in drinking water is well-known, widespread and global in nature.¹⁻³ PFAS drinking water levels are formally regulated, or under proposed regulations, in several countries and regions, including the United States⁴ and the European Union.⁵ In the United States, the Environmental Protection Agency (EPA) has promulgated several methods for PFAS analysis in drinking water [EPA 537, 537.1, 533] as well as non-potable water samples such as surface water, groundwater and wastewater [EPA 1633, 1633A]. The EPA methods utilize solid-phase extraction (SPE) for sample concentration and clean-up. Traditionally, the SPE procedure has been manually performed with vacuum manifolds, which is time-intensive and occasionally error-prone. However, in recent years automated SPE systems have become commercially available with the aim of increasing throughput, reducing manual effort and improving reproducibility. These systems can extract multiple samples in parallel, following an EPA-compliant method, and can be modified to reduce fluoropolymer components, reducing background PFAS levels. In this technical note, a method is described for PFAS analysis in drinking water following EPA Method 533, using the PromoChrom SPE-03 automated SPE system and the SCIEX QTRAP 6500+ system.

Methods

Samples and reagents: Calibration standards were purchased from Phenova [Golden, Colorado] and Absolute Standards [Hamden, CT]. Stable-isotope labelled standards were purchased from Wellington Laboratories [Guelph, Ontario].

Quality control (QC) samples: The method performance was evaluated using 3 types of QCs, laboratory reagent blank [LRB], laboratory fortified blank [LFB] and, laboratory fortified sample matrix [LFSM] and duplicate [LFSMD] samples. LRB samples [n=38] consisted of Milli-Q water spiked with the isotope dilution analogues [IDAs] only. LFBs were comprised of Milli-Q water spiked with the target analytes and IDAs. LFB spikes were performed at 2 ng/L [n=38], 40 ng/L [n=38] and 70 ng/L [n=37]. LFSM and LFSMD samples were from untreated groundwater samples spiked with the target analytes at 2 ng/L [n=13] and 40 ng/L [n=23], as well as the IDA mix. All QC samples were

extracted and instrumentally analyzed using the procedures described below.

Real-world groundwater samples: 10 raw untreated groundwater samples collected from the influent of a PFAS treatment facility were analyzed between February 2024 and January 2025. These samples represent source water that had not yet undergone treatment and do not represent the final treated drinking water delivered to the community. The purpose of analyzing the raw groundwater samples was to demonstrate the applicability of the PromoChrom extraction and instrumental analysis methods.

Sample preparation: All samples were extracted during May-September 2024 using the Promochrom SPE-03 automated SPE system with the MOD-004 [sample bottle rinsing] and MOD-005 [minimal Teflon option] modifications. The automated SPE procedure is outlined in **Table 3**. The sample volume was 250 mL, and the SPE cartridge was the Phenomenex Strata SBD-L [500 mg/5 mL, [P/N: 8B-S014-HCH](#)]. QC samples were collected between May and September 2024 and were comprised of 38 laboratory reagent blanks [LRB], 38 laboratory fortified blanks [LFB] at 2, 40 and 70 ng/L, 13 laboratory fortified sample matrix [LFSM] and duplicate [LFSMD] samples at 2 ng/L, and 23 LFSM and LFSMD samples at 40 ng/L.

LC chromatography: Chromatographic separation was performed using an Agilent 1200 LC system. The analytical column was a Phenomenex Gemini C18 column [3 μ m, 100 x 2 mm, P/N: [00D-4439-B0](#)] and the delay column was a Phenomenex Luna C18[2] column [5 μ m, 30 x 2 mm, P/N: [00A-4252-B0](#)]. Mobile phase A was water with 20mM ammonium acetate, and mobile phase B was methanol. The gradient conditions used are presented in **Table 1**. The flow rate was 430 μ L/min, the injection volume was 5 μ L, and the column oven was 40°C.

Table 1: Chromatographic gradient for the analysis of PFAS in drinking water.

Time [min]	Mobile phase A [%]	Mobile phase B [%]
0.0	95	5
0.1	45	55
5.4	1	99
9.0	1	99
9.1	95	5
11	95	5

Mass spectrometry: Samples were analyzed using the [QTRAP 6500+ system](#) with electrospray ionization operated in negative mode. Data was acquired using multiple reaction monitoring (MRM) modes using optimized source and gas conditions ([Table 2](#)) and compound-specific parameters ([Table 4 in the Appendix](#)).

Data processing: Data acquisition was performed using Analyst (version 1.7.2) and processed using MultiQuant (version 3.0.3).

Table 2: Source and gas parameters for the analysis of PFAS in drinking water using the QTRAP 6500+ system.

Parameter	Value
Polarity	Negative
Ion source gas 1	40 psi
Ion source gas 2	40 psi
Curtain gas	35 psi
Source temperature	300°C
Ion spray voltage	-4500 V
CAD gas	10 psi

Table 3: Automated Promochrom extraction procedure for EPA Method 533.

Step	Action	Inlet	Flow [mL/min]	Volume [mL]	Time [min]
1	Elute W2	Methanol	5	5	
2	Wait				1
3	Elute W2	Methanol	3	5	
4	Wait				2
5	Elute W1	0.1M Phosphate buffer	5	5	
6	Wait				2
7	Elute W1	0.1M Phosphate buffer	5	5	
8	Wait				1
9	Elute W1	0.1M Phosphate buffer, then Milli-Q water	3	5	
10	Wait				2
11	Add sample W1	Sample	5	285	
12	Rinse W1	QC water	10	7.5	
13	Rinse W1	QC water	10	7.5	
14	Add sample W1	Sample	10	4.5	
15	Elute W2	Methanol	10	1	
16	Air-purge1	Air	10	5	
17	Blow N ₂				5 @ 2.5 L/min
18	Rinse 2	98:2 (w/w) methanol/ammonium hydroxide	5	5	
19	Wait				2
20	Rinse 2	98:2 (w/w) methanol/ammonium hydroxide	5	5	
21	Wait				2
22	Collect 1	Sample	5	4.5	
23	Air-purge1	Air	10	10	

W1 = aqueous waste, W2 = organic waste

System background contamination: Laboratory reagent blanks (LRB)

The extensive series of LRB samples (n=38) evaluated the contamination originating from the PromoChrom system during the sample preparation protocol. None of the targeted PFAS analytes were detected except for one sample that showed low 6:2 FTS levels [2.2 ppt]. **Figure 2** shows representative XICs for the 6 EPA regulated PFAS compounds in the LRBs. These results demonstrate the ability of the PromoChrom automated SPE system to generate extracts with very low background PFAS levels. This is critical for achieving the low parts-per-trillion [ppt] reporting levels necessary to meet the 2024 EPA drinking water maximum contaminant levels [MCLs]: 4 ppt for PFOA and PFOS, 10 ppt for PFNA, PFHxS and HFPO-DA.

Method performance in quality control spikes: Fortified blank and matrix samples

Method performance was evaluated in two different types of QC samples: LFBs and LFSMs/LFSMDs. LFBs consisted of spiked Milli-Q water samples processed through the PromoChrom system and instrumentally analyzed. The LFBs were prepared at three spiking levels: 2 ng/L (n=38), 40 ng/L (n=38) and 70 ng/L (n=37). The EPA 533 criteria specify that the LFB recovery must be within 50-150% at spiking levels near the 2 ng/L method reporting limit [MRL] and within 70-130% for the higher spike levels. Considering all targeted PFAS analytes, the mean LFB recovery ranged from 99.5% to 115% for the 2 ng/L spikes, 95.7% to 105% for the 40 ng/L spikes and 95.6 to 103% for the 70 ng/L spikes. **Figure 2** shows representative XICs for the 2 ng/L LFB spikes. **Figure 3** shows the mean analyte recovery for each spiking level and associated %CV bars as well as the $\pm 30\%$ accuracy tolerance lines [red dashed line]. The mean precision was <10%CV for all analytes across the 3 spiking levels.

The LFSM and LFSMD samples consisted of spiked untreated groundwater samples processed through the PromoChrom system and instrumentally analyzed. The corresponding unspiked matrix samples were also analyzed to correct for background PFAS levels present. The LFSM and LFSMD samples were prepared at 2 ng/L (n=13) and 40 ng/L (n=23). Similar to the LFB samples, the EPA 533 criteria specify 50-150% recovery at the 2 ng/L MRL and 70-130% recovery for higher spiking concentrations. **Table 4** shows the mean analyte recovery and precision [%CV] for the LFSM and LFSMD samples. The mean recovery range in the 2 ng/L spikes was 98-122% and 97-118% for the LFSM and LFSMD samples. For the 40 ng/L spikes, the mean recovery range was 96-110% and 95-109% in LFSM and LFSMD samples. The mean precision was <13%CV for all analytes in the 2 spiking levels. The mean relative percent difference [RPD] for the duplicate samples was 7% and 6% in the 2 ng/L and 40 ng/L spikes, respectively. All RPDs were within the EPA 533 criteria of <50% for the MRL spike and <30% for higher-level spikes.

Overall, the results from both the fortified reagent blank [LFBs] and sample matrix and duplicate [LFSMs/LFSMDs] samples demonstrate the ability of the PromoChrom automated SPE system and QTRAP 6500+ system to generate excellent quantitative data for PFAS in drinking water. Excellent recovery and precision were observed at both the MRL and higher levels.

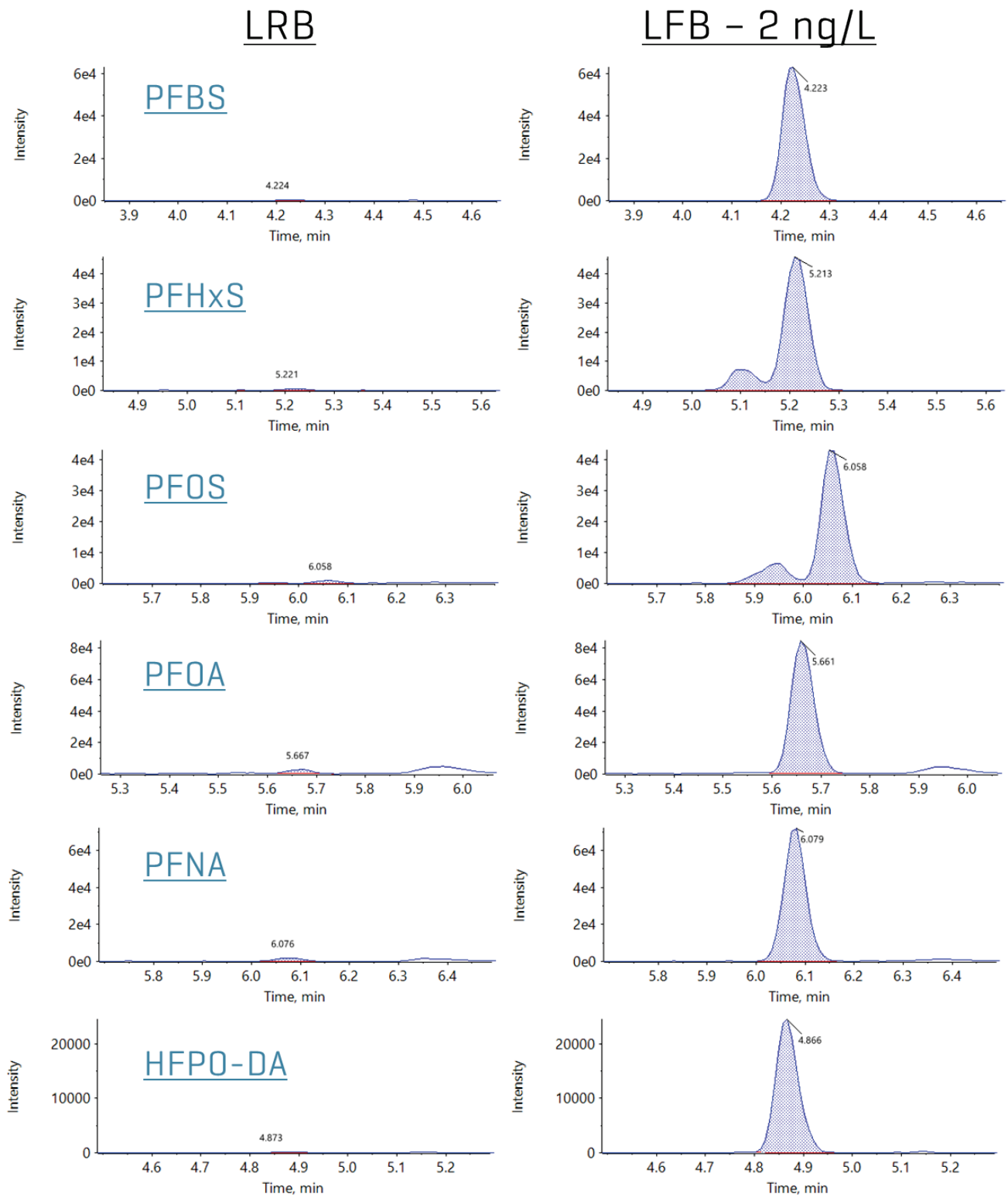


Figure 2. Extracted ion chromatograms (XICs) in representative laboratory reagent blank and laboratory fortified blank [2 ng/L spike level] samples for the six EPA regulated PFAS compounds [PFBS, PFHxS, PFOS, PFOA, PFNA and HFPO-DA].

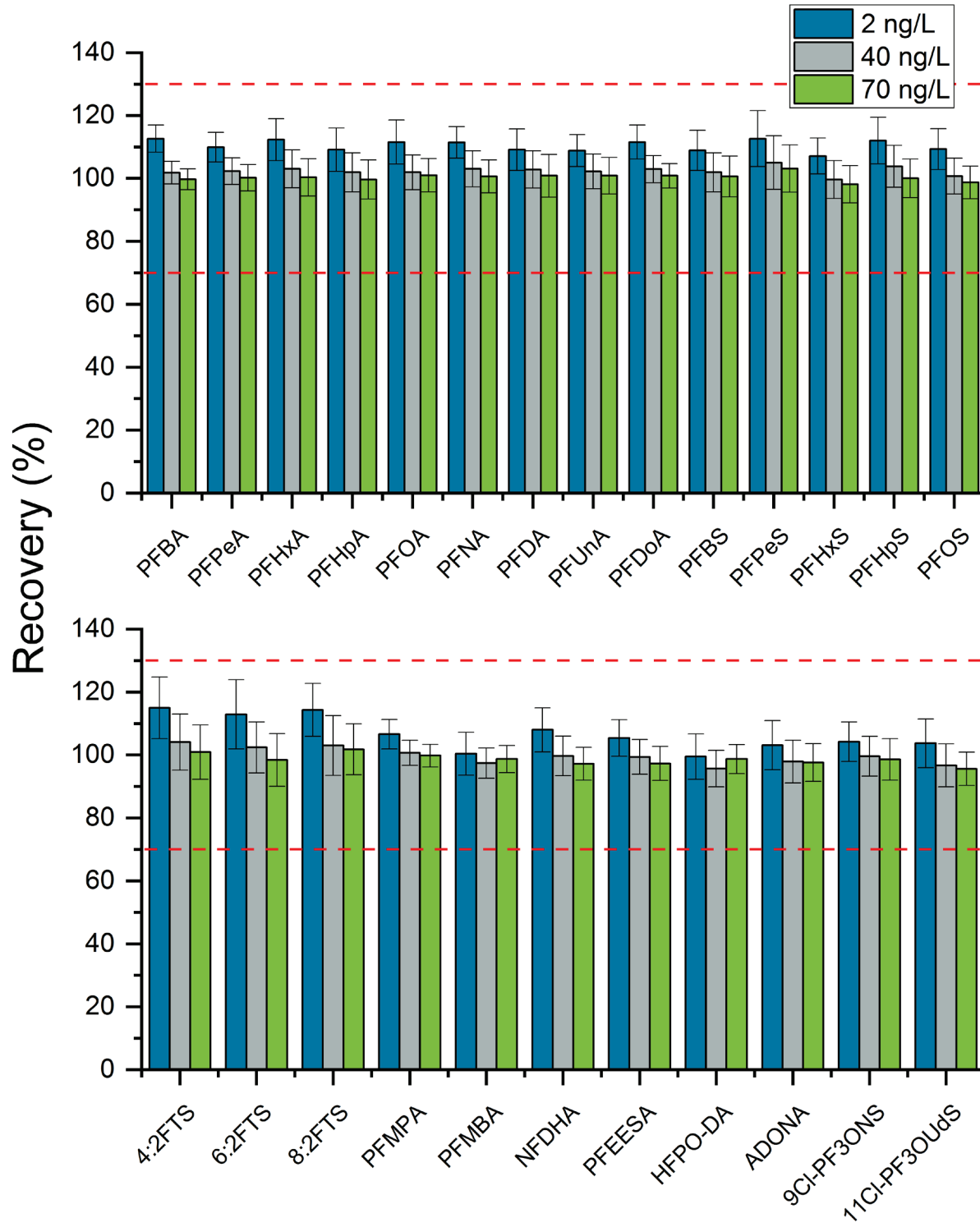


Figure 3. Mean PFAS recovery in laboratory fortified blank samples (n=38 for 2 ng/L and 40 ng/L spikes, n=37 for 70 ng/L spike). Error bars represent the coefficient of variation percentage [%CV]

Table 4: Mean recovery [%] and precision [%CV] in the laboratory fortified sample matrix (LFSM) and duplicates (LFSMD) for 2 ng/L (n=13) and 40 ng/L (n=23) spikes.

	2 ng/L - LFSM		2 ng/L - LFSMD		40 ng/L - LFSM		40 ng/L - LFSMD	
	Recovery [%]	%CV	Recovery [%]	%CV	Recovery [%]	%CV	Recovery [%]	%CV
PFBA	111	8.7	112	8.2	102	3.7	101	3.5
PFPeA	111	13	113	7.2	103	4.5	103	6.0
PFHxA	117	7.9	116	7.2	103	5.8	101	6.7
PFHpA	112	8.6	111	7.0	102	6.3	102	4.8
PFOA	114	7.7	111	8.6	102	6.2	102	7.9
PFNA	117	6.9	115	6.7	100	6.6	103	5.1
PFDA	114	7.1	115	6.8	103	6.6	102	5.9
PFUnA	115	6.4	113	7.1	101	5.5	102	6.8
PFDoA	114	9.3	112	7.2	104	5.2	103	5.3
PFBS	111	6.8	111	6.3	101	6.2	101	6.5
PFPeS	116	7.4	112	10	104	10	107	7.0
PFHxS	110	9.4	106	7.7	98.0	5.7	99.7	7.9
PFHpS	117	10	116	9.4	104	6.4	102	7.0
PFOS	113	9.5	113	7.2	101	6.9	99.9	6.5
4:2FTS	112	10	118	8.7	102	8.9	105	6.9
6:2FTS	122	8.6	115	12	102	10	98.5	9.6
8:2FTS	117	9.9	115	11	106	8.0	105	9.9
PFMPA	110	6.4	112	6.0	110	12	109	11
PFMBA	101	8.5	102	10	96.9	5.3	95.6	5.2
NFDHA	110	6.4	112	6.1	103	8.2	101	7.4
PFEESA	106	8.3	109	7.3	96.0	7.8	95.2	6.8
HFPO-DA	98.2	7.7	97.0	7.0	99.6	9.9	98.3	9.3
ADONA	108	5.8	107	6.5	99.9	7.7	100	7.2
9Cl-PF3ONS	108	9.1	104	8.4	97.3	6.9	97.7	5.3
11-Cl-PF3OUdS	105	12	107	6.3	98.5	5.4	105	6.9

Real-world samples

The PromoChrom automated SPE sample preparation method was applied to raw untreated groundwater samples collected from February 2024 to January 2025 and analyzed using the 6500+ system. Concentrations were lowest for PFNA [range: 2.1-3.3 ng/L] and highest for PFOS [range: 54.7-73.7 ng/L] while HFPO-DA levels were <MRL. See **Table 5** for the full data set. The samples were collected prior to treatment, and after undergoing treatment, all samples were non-detect for all PFAS analytes.

Conclusions

- Automated sample preparation using the PromoChrom SPE system produced drinking water extracts with low background contamination, satisfactory to meet EPA maximum contaminant levels [MCLs] for PFAS
- Excellent PFAS recovery and precision in QC spike samples, meeting data quality requirements in EPA Method 533 using the SCIEX 6500+ system

References

1. Ackerman Grunfeld, D.; Gilbert, D.; Hou, J. *et al.* Underestimated burden of per- and polyfluoroalkyl

substances in global surface waters and groundwaters. *Nat. Geosci.* **2024**, *17*, 340-346. DOI: [10.1038/s41561-024-01402-8](https://doi.org/10.1038/s41561-024-01402-8).

2. Smalling, K.L.; Romanok, K.M.; Bradley, P.M. *et al.* Per- and polyfluoroalkyl substances (PFAS) in United States tapwater: Comparison of underserved private-well and public-supply exposures and associated health implications. *Environ. Int.* **2023**, *178*, 108033. DOI: [j.envint.2023.108033](https://doi.org/10.1016/j.envint.2023.108033).
3. European Environment Agency. *PFAS pollution in European waters*. December 9, 2024. <https://www.eea.europa.eu/en/analysis/publications/pfas-pollution-in-european-waters> [accessed 2025-02-10].
4. United States Environmental Protection Agency. *Per- and polyfluoroalkyl substances (PFAS): Final PFAS national primary drinking water regulation*. <https://www.epa.gov/sdwa/and-polyfluoroalkyl-substances-pfas> [accessed 2025-02-10].
5. Directive [EU] 2020/2184 of the European Parliament and of the Council of 16 December 2020 on the quality of water intended for human consumption. <http://data.europa.eu/eli/dir/2020/2184/oj> [accessed: 2025-02-10].

Table 5. PFAS concentrations for the 6 EPA-regulated compounds in raw untreated groundwater samples collected between February 2024 and January 2025.

Date	PFBS (ng/L)	PFHxS (ng/L)	PFOS (ng/L)	PFOA (ng/L)	PFNA (ng/L)	HFPO-DA (ng/L)
02/21/24	7.6	18.8	59.5	21.6	2.1	< MRL
02/20/24	8.2	18.9	62.5	21.0	2.2	< MRL
02/23/24	8.1	18.6	54.7	21.7	2.1	< MRL
03/27/24	11.2	20.9	61.2	24.4	2.8	< MRL
03/26/24	9.1	19.9	59.7	23.6	2.5	< MRL
04/10/24	10.7	21.6	63.2	26.2	2.9	< MRL
05/20/24	10.4	21.3	65.4	29.2	3.0	< MRL
07/10/24	9.2	18.0	73.7	23.0	2.9	< MRL
10/03/24	11.0	20.7	68.1	27.0	2.9	< MRL
01/06/25	11.2	19.9	68.4	27.1	3.3	< MRL

Appendix

Table 4: Compound-specific MRM parameters for the analysis of PFAS in drinking water by EPA Method 533 using the QTRAP 6500+ system.

	Precursor ion [Da]	Fragment ion [Da]	DP [V]	CE [V]	CXP [V]
PFBA	213.0	168.9	-10	-14	-7
PFPeA	262.9	218.9	-5	-12	-21
PFHxA	312.9	268.9	-5	-14	-23
PFHpA	362.9	318.8	-15	-14	-21
PFOA	413.0	368.9	-15	-14	-27
PFNA	462.9	418.9	-5	-16	-19
PFDA	512.8	468.9	-5	-16	-27
PFUnA	562.9	518.9	-5	-16	-7
PFDoA	612.9	568.8	-15	-18	-29
PFBS	298.9	79.9	-5	-68	-11
PFPeS	348.8	79.9	-85	-74	-9
PFHxS	398.8	79.9	-60	-86	-13
PFHpS	448.9	79.9	-30	-100	-9
PFOS	498.8	79.9	-85	-108	-39
4:2FTS	326.9	306.9	-55	-28	-34
6:2FTS	426.9	406.8	-55	-32	-19
8:2FTS	526.9	506.8	-60	-40	-31
PFMPA	229.0	84.9	-5	-20	-11
PFMBA	278.9	84.9	-5	-22	-9
NFDHA	294.9	200.9	-5	-14	-17
PFEESA	314.9	134.9	-30	-30	-13
HFPO-DA	284.7	168.9	-15	-12	-17
ADONA	376.9	250.9	-5	-14	-17
9CI-PF3ONS	530.8	350.8	-20	-36	-23
11-CI-PF3OUdS	630.8	450.8	-5	-42	-25

The SCIEX clinical diagnostic portfolio is For In Vitro Diagnostic Use. Rx Only. Product(s) not available in all countries. For information on availability, please contact your local sales representative or refer to <https://sciex.com/diagnostics>. All other products are For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd. or their respective owners in the United States and/or certain other countries [see www.sciex.com/trademarks].

© 2025 DH Tech. Dev. Pte. Ltd. MKT-34209-A



Headquarters

500 Old Connecticut Path | Framingham, MA 01701
USA
Phone 508-383-7700
sciex.com

International Sales

For our office locations please call the division headquarters or refer to our website at sciex.com/offices