

A sensitive method for the quantitation of bimatoprost in human plasma

Sensitive quantitative performance on the SCIEX 7500 system with improved front-end technology

Lakshmanan Deenadayalan,¹ Sashank Pillai,¹ Rahul Baghla² and Eshani Nandita²

¹SCIEX, India; ²SCIEX, USA

This technical note demonstrates a sensitive method for the quantitation of an ocular drug, bimatoprost, in human plasma. The method employs a simple liquid-liquid extraction sample preparation to separate bimatoprost from human plasma. A lower limit of quantitation (LLOQ) of 0.5 pg/mL was achieved with %CV <3% (Figure 1).

Eye diseases, next to cardiovascular and respiratory diseases, are the most common among the general population.¹ Aging is one of the major risk factors for glaucoma, which leads to optic nerve cell damage and ultimately causes blindness.² Glaucoma develops when fluids cannot drain from the eye, leading to increased pressure. Common treatments include laser surgery or treatment with prostaglandin analogs and topical drugs (eye drops) such as bimatoprost and latanoprost.

Bimatoprost is a synthetic prostamide that is similar in structure to prostaglandin F₂α (PGF₂α). It functions by substituting an electrochemically neutral amide group for the carboxylic acid group in PGF₂α. Glaucoma and other eye conditions caused by high blood pressure can be effectively treated with bimatoprost. The usual dosage is one drop of 0.03% bimatoprost. It predominantly resides in plasma and has an 88% plasma protein binding rate. Bimatoprost has an estimated C_{max} of 60-80 pg/mL and the maximum plasma concentration is often reached in less than 10 minutes.⁵ Therefore, it is important to have a sensitive bioanalytical method for the quantitation of bimatoprost across a wide range of concentrations in biological matrices to support pharmacokinetic studies.⁴

The bioanalytical method described in this technical note uses 400 μL of human plasma and a simple liquid-liquid extraction procedure to detect low levels of bimatoprost.

Key features for the analysis of bimatoprost using the SCIEX 7500 system

- **Sensitive and reproducible of quantitation of low-dose, high-potency drug modalities:** Achieve a 0.50 pg/mL LLOQ for bimatoprost in human plasma with %CV <3%
- **Low plasma consumption and simple sample preparation:** Use a low sample volume of 400 μL of human plasma paired with a simple liquid-liquid extraction methodology to quantify bimatoprost
- **Enhanced sensitivity unlocked:** Improved front-end technology with the D Jet ion guide, OptiFlow Pro ion source and E Lens probe for enhanced ion generation, capture and transmission, enable users to reach their desired quantitative sensitivity
- **Streamlined data management:** Data acquisition and processing are integrated into SCIEX OS software, a 21 CFR Part 11 compliance-ready platform

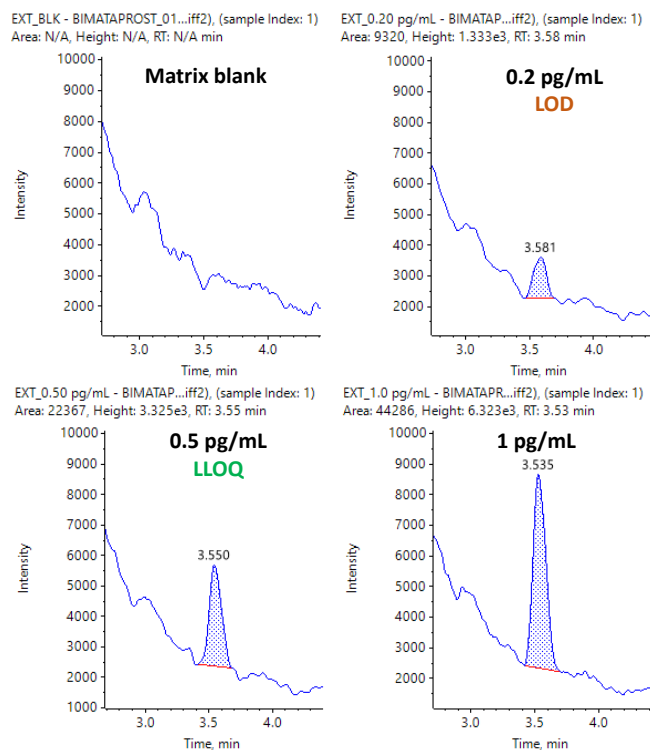


Figure 1. Representative extracted ion chromatograms (XICs) for bimatoprost in human plasma. The XICs of the matrix blank (top left), limit of detection (LOD, 0.2 pg/mL, top right), LLOQ (0.5 pg/mL, bottom left) and 1 pg/mL (bottom right) are shown.

Methods

Spiked sample preparation: Bimatoprost was spiked into 400 μ L of human plasma at concentrations ranging from 0.2 to 500 pg/mL. A 100 μ L aliquot of 0.1N sodium hydroxide in water was added to the sample and vortexed for 2 minutes. A 4 mL aliquot of 80:20 (v/v), ethyl acetate/n-hexane was added to the samples, vortexed for 10 minutes and centrifuged at 1204 rcf for 5 minutes. The supernatant was collected and dried under a nitrogen stream at 40°C. The dried samples were reconstituted using 150 μ L of 30:70 (v/v), acetonitrile/5mM ammonium formate in water. A 20 μ L sample injection was used for analysis.

Chromatography: An ExionLC system with a Phenomenex [Kinetex XB-C18 column \(2.1 x 100 mm, 1.7 \$\mu\$ m, 100 Å\)](#) was used for chromatographic separation. The LC column was operated at 40°C. Mobile phase A was 5mM ammonium formate in water and mobile phase B was acetonitrile. Table 1 summarizes the LC gradient conditions used.

Table 1. LC gradient.

Time (min)	Flow (mL/min)	Mobile phase A (%)	Mobile phase B (%)
0.0	0.6	70	30
4.5	0.6	70	30
4.6	0.6	20	80
5.5	0.6	20	80
5.6	0.6	70	30
7.0	0.6	70	30

Mass spectrometry: Samples were analyzed using the [SCIEX 7500 system](#) equipped with the OptiFlow Pro ion source. The system was controlled by SCIEX OS software. The optimized source and MS parameters are listed in Table 2.

Data processing: Data processing was performed using SCIEX OS software, version 3.1.6. Peaks were automatically integrated using the MQ4 algorithm with a weighting of $1/x^2$.

Table 2. Optimized MS parameters.

Name	Q1/Q3 (m/z)	Q0D (V)	CE (V)	CXP (V)
Bimatoprost	398.2/362.2	25	14	12
Bimatoprost d_5	403.3/367.2	10	14	12
Source parameters	Value	Source parameters	Value	
Curtain gas	40 psi	CAD gas	7	
Ion source gas 1	30 psi	Ion spray voltage	1500 V	
Ion source gas 2	85 psi	Temperature	350°C	

Quantitative performance

A calibration curve was analyzed for concentrations ranging from 0.2 to 500 pg/mL. To evaluate reproducibility, each calibration standard was analyzed in triplicate.

Bimatoprost in human plasma was quantified at an LLOQ of 0.5 pg/mL with an LOD of 0.2 pg/mL. No interferences were observed in the matrix blank (Figure 1). Linearity was achieved across concentrations ranging from 0.2 to 800 pg/mL with a correlation of determination (r^2) of 0.994 (Figure 2). A linear dynamic range (LDR) of 3 orders of magnitude was achieved.

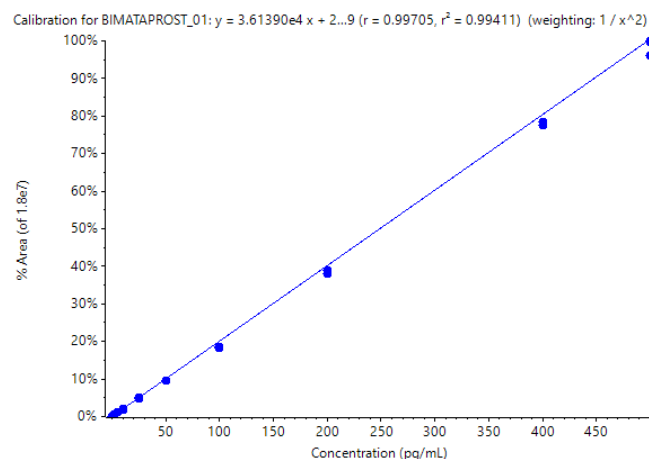


Figure 2. Calibration curve for the quantitation of bimatoprost in human plasma.

Analytical performance was evaluated based on the requirement that the accuracy of the calculated mean should be between 80% and 120% at the LLOQ and between 85% and 115% at higher concentrations. The %CV of the calculated mean concentration should be below 20% at the LLOQ and below 15% at all higher concentrations.³

The assay accuracy was within $\pm 10\%$ of the nominal concentration and the %CV was $< 6\%$ (Table 3), including at

the LOD. The calculated accuracy and %CV values were within the acceptance criteria at each concentration level (Table 3).

Table 3. Summary of the quantitative performance. Reproducibility and accuracy results were determined from the calibration curve across 3 replicates.

Concentration (pg/mL)	CV (%)	Accuracy (%)
0.2	5.14	93.3
0.5	2.97	110
1	2.45	109
2.5	1.46	108
5	0.62	106
10	1.99	101
25	2.23	96
50	0.71	97
100	2.10	97.9
200	0.64	95.1
400	2.85	94.8
500	1.71	92

Bimatoprost is a synthetic prostamide that has a similar structure to PGF₂α (Figure 3). Given its similarity to a lipid structure, these compounds can be difficult to separate from other phospholipids present in the human plasma. However, bimatoprost was successfully extracted from human plasma using a simple liquid-liquid extraction method that led to a recovery of >90% (Table 4).

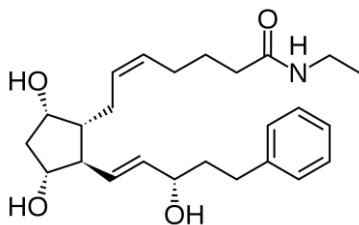


Figure 3. Chemical structure of bimatoprost.

Table 4. Recovery results using 5 pg/mL and 200 pg/mL of bimatoprost. Each concentration was evaluated in triplicate.

Concentration (pg/mL)	Response in post-spiked samples (Mean area ± SD)	Response in pre-spiked samples (Mean area ± SD)	% Recovery
5	224176±4153	208404±5326	93.0
200	7640368±56560	6953091±71700	91.0

Carryover was evaluated by injecting a blank sample before and after the upper limit of quantitation (ULOQ) at 500 pg/mL. Figure 4 shows no carryover was observed in the XICs run before and after the ULOQ at the retention time of the analyte (3.55 minute).

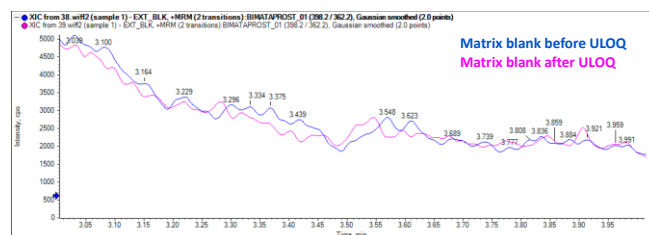


Figure 4. XICs of bimatoprost from matrix blank injection before (blue) and after (pink) the ULOQ (500 pg/mL).

Compliance-ready SCIEX OS software

SCIEX OS software is a closed system and requires records and signatures to be stored electronically, meeting the regulations outlined by 21 CFR Part 11. SCIEX OS software can open raw data files from any visible storage location within a closed network by using designated processing workstations. Figure 5 illustrates the features of SCIEX OS software used for monitoring the audit trail, acquiring and processing data and configuring user access.

The audit trail feature enables users to audit critical user actions and locks in data integrity. The Central Administrator Console (CAC) feature allows users to centralize acquisition and processing using a single platform to maximize efficiency for multi-instrument laboratories, independent of compliance standards. The configuration module allows users to assign roles and access as the administrator, method developer, analyst and reviewer.

Easily search and filter for specific high-risk events in audit trail viewer. Built-in data integrity features allow you to tailor each functionality specifically to meet compliance needs and data security requirements.

Filter Audit Trail dialog box with search criteria: Sample Name, Reason, and time range (From/To).

Users can manage groups, role definitions, workstations and projects across all systems using the SCIEX OS software Central Administrator Console (CAC). It supports all regulated and non-regulated compliance standards.

Central Administrator Console (CAC) interface showing 'Manage Workgroups and Assignments' and a 'Confirm Change Events' dialog box.

Assign users and access to administrator, method developer, analyst and reviewer roles under the audit trail module. Easily customize the role and specify level of access.

Configuration interface showing 'User Roles and Permission Categories' table and a 'Duplicate a User Role' dialog box.

Permission	Administrator	Method Developer	Analyst	Reviewer
Batch				
Submit unlocked methods	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Open	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Save as	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Submit	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Save	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Save ion reference table	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Add data sub-folders	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Configure Decision Rules	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Configuration				
General tab				
General: change regional setting	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
General: full screen mode	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
LIMS communication tab	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Audit maps tab	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Queue tab	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Queue: instrument idle time	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Queue: max. number of acquired samples	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
Queue: other queue settings	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>

Figure 5. Features of SCIEX OS software for monitoring user access and evaluating the audit trail. The audit trail view allows users to filter for high-risk events easily and enables data integrity features to meet compliance requirements. The software features a Central Administrator Console (CAC) to manage users and groups, role definitions, workstations and projects across all systems. The CAC feature supports both regulated and non-regulated compliance standards. The configuration module enables users to quickly set up roles and levels of access for the administrator, method developer, analyst and reviewer levels.

Conclusions

- An LLOQ of 0.50 pg/mL was achieved to quantify bimatoprost in human plasma with an LOD of 0.2 pg/mL
- Linearity was achieved at concentrations ranging from 0.2 pg/mL to 500 pg/mL with an r^2 of 0.994
- Desired LLOQ (0.5 pg/mL) for bimatoprost was achieved with 400 μ L of human plasma
- The method demonstrated accurate and highly reproducible quantitative performance (%CV <6%) for all concentration levels, including the LOD
- Sensitivity was achieved on the SCIEX 7500 system with an improved front-end technology for better ion generation, capture and transmission
- SCIEX OS software is compliance-ready to support 21 CFR Part 11 and integrates with a triple quadrupole mass spectrometer to support data acquisition, processing and management on a single platform

References

1. Trends in development and quality assessment of pharmaceutical formulations - F2 α analogues in the glaucoma treatment. [European Journal of Pharmaceutical Sciences, 180, Jan 2023, 106315.](#)
2. The number of people with glaucoma worldwide in 2010 and 2020. [Br. J. Ophthalmol. 90, 262–267.](#)
3. Bioanalytical Method Validation, [May 2018.](#)
4. Clinical pharmacology and biopharmaceutics review(s), [211911Orig1s000.](#)
5. Assessment report for Lumigan, 7 Jan 2010, [Doc. Ref:EMA/105752/2010](#)

The SCIEX clinical diagnostic portfolio is For In Vitro Diagnostic Use. Rx Only. Product(s) not available in all countries. For information on availability, please contact your local sales representative or refer to <https://sciex.com/diagnostics>. All other products are For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd. or their respective owners in the United States and/or certain other countries (see www.sciex.com/trademarks).

© 2023 DH Tech. Dev. Pte. Ltd. MKT-28441-A .



Headquarters
500 Old Connecticut Path | Framingham, MA 01701 USA
Phone 508-383-7700
sciex.com

International Sales
For our office locations please call the division headquarters or refer to our website at sciex.com/offices