

# Sensitive bioanalysis of galactosyl sphingosine (GalSPH) and glucosyl sphingosine (GluSPH) in cerebral spinal fluid

## Using the SCIEX 7500 system

Jack Steed<sup>1</sup>, Jessica Smith<sup>1</sup>, Freddy Oostebring<sup>2</sup>, Jianru Stahl-Zeng<sup>3</sup>  
<sup>1</sup>SCIEX, UK; <sup>2</sup>Ardena Bioanalysis Assen, The Netherlands; <sup>3</sup>SCIEX, Germany

GalSPH and GluSPH are useful biomarkers for several diseases such as Krabbe and Gaucher disease as well as being a potential marker for carriers of heterozygous mutated GBA1 gene, which increases the risk of Parkinson's.<sup>1,2</sup> Therefore, it is important to be able to accurately monitor these compounds within the brain. Cerebral spinal fluid (CSF) is analyzed for GalSPH and GluSPH due to being the closest matrix available to monitor changes in the brain. However, the volume available is extremely limited as typically several biomarker assays need to be performed on a single sample. Therefore, it is important to have as much sensitivity as possible to ensure that lower volumes of CSF can be used.

Currently, this analysis is performed by Ardena Bioanalysis Assen using the SCIEX Triple Quad 6500+ system but, due to the increased sensitivity available with the SCIEX 7500 system, a comparison study to evaluate impact was performed. The objective was to assess whether the increase in sensitivity allows for lower CSF volumes to be analyzed.




The analysis was performed using HILIC chromatography to provide separation between the two isomers (Figure 1), with the triple quadrupole mass spectrometers being operated in positive mode (electrospray ionization) using MRM acquisition.

## Methods

**Sample preparation:** Standards for galactosyl sphingosine (GalSPH) and glucosyl sphingosine (GluSPH) and their deuterated internal standards, GluSPH-D<sub>5</sub> and GalSPH-D<sub>5</sub>, were obtained from Avanti Polar Lipids. Extracted standards and cerebral spinal fluid (CSF) samples were prepared by Ardena Bioanalysis. Standards were prepared in surrogate CSF across a range of concentrations suitable for analyzing endogenous levels of the CSF samples, with the deuterated internal standards held add a constant concentration. The samples were then split to be analyzed on both the SCIEX Triple Quad 6500+ and SCIEX 7500 systems.

**LC-MS analysis:** Samples were analyzed using both the QTRAP 6500+ system and the SCIEX 7500 system. Separation of Gal- and Glu-SPH was achieved using HILIC chromatography. MRM analysis was performed using electrospray ionization in positive polarity.

**Data processing:** Datasets were processed using both SCIEX OS software and Analyst software.

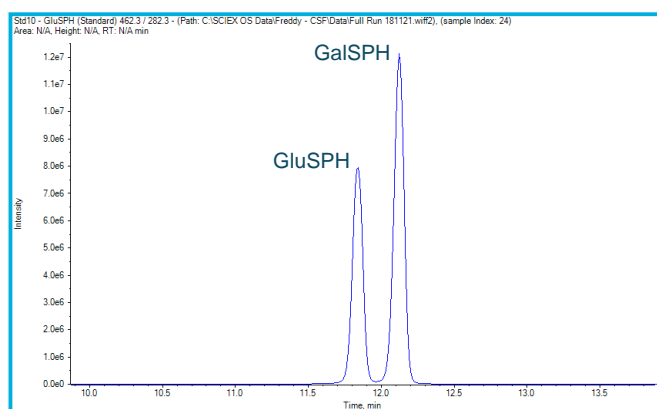


**Figure 1. Extracted ion chromatograms (XIC) of GluSPH and GalSPH.** Using HILIC chromatography, good separation, and peak shape of the two isomers was achieved. Both compounds are analyzed using the same MRM transition and so chromatographic separation is paramount.

## Results

To compare the sensitivities between the two systems, a set of standards and CSF samples were prepared by Ardena Bioanalysis Assen before being divided and analyzed on both MS systems. Figure 2 highlights the calibration curves obtained for GluSPH and GalSPH on the SCIEX 7500 system across a concentration range between 0.001 – 1.000 pmol/mL. Both provided an  $r$  value  $>0.99$ , with accuracies between 80 – 120% at each level (Table 1), providing good linearity across the 3 orders of dynamic range studied. A similar concentration curve was generated on the QTRAP 6500+ system (data not shown).

The %RSD and calculated concentrations for the bottom four surrogate CSF standards are also detailed in Table 2 highlighting that the %RSD values are below acceptable criteria of 10%.

The sensitivity difference between the two systems was next assessed by using peak-to-peak S/N with similar noise regions, across the concentration curves. A summary is shown in Table 3 for GluSPH and Table 4 for GalSPH. The S/N increase was found to be between 2.4 and 6.7x, with improvements seen across the concentration range analyzed.

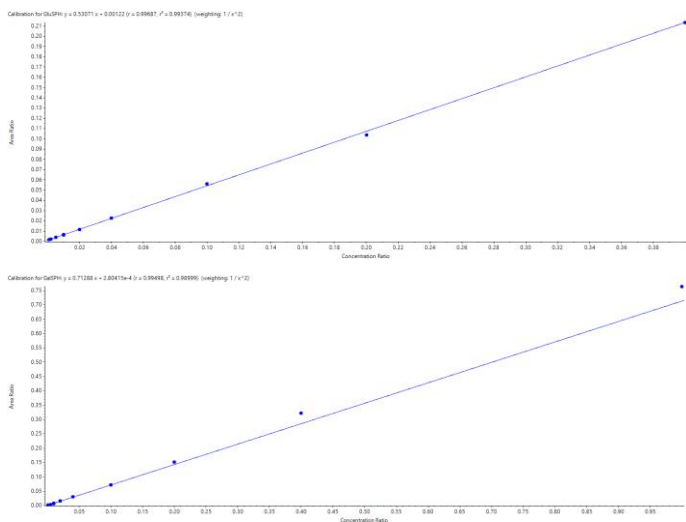

**Figure 2, Calibration curves using the SCIEX 7500 system.**

Calibration curves for both GluSPH (top) and GalSPH (bottom) were generated in surrogate CSF. Both have an  $r$  value  $> 0.99$  and use a  $1/X^2$  weighting, with a range between 0.001 – 1.000 pmol/mL. Results were obtained using internal standard correction with the deuterated forms of the analytes (GluSPH- $D_5$  and GalSPH- $D_5$ ).

**Table 1. The % accuracy values for GalSPH and GluSPH on the SCIEX 7500 system.** The accuracy values calculated from the calibration curve generated in surrogate CSF were within 80-120% for all concentrations analyzed.

Concentration (pmol/mL)	% Accuracy (GluSPH)	% Accuracy (GalSPH)
0.001	101.23	103.21
0.002	96.08	99.03
0.005	103.61	86.66
0.010	99.92	97.04
0.020	100.19	110.23
0.040	101.06	105.15
0.100	103.81	100.55
0.200	97.13	106.42
0.400	100.49	112.73
1.000	94.81	107.08

**Table 2. The %RSD and average calculated concentrations for GalSPH and GluSPH on the SCIEX 7500 system.** The peak area %RSD was calculated for the bottom 4 concentrations ( $n=3$ ) and was below acceptable criteria of 10% at each level assessed. The average calculated concentration (pmol/mL) was derived based on the calibration curve analyzed to show similarity to the theoretical values.

Concentration (pmol/mL)	%RSD of area (GluSPH)	Average calculated concentration (GluSPH, pmol/mL)	%RSD of area (GalSPH)	Average calculated concentration (GalSPH, pmol/mL)
0.0010	8.75	0.0010	4.07	0.0010
0.0020	8.97	0.0019	1.48	0.0020
0.0050	8.08	0.0052	3.14	0.0043
0.0100	4.13	0.0100	1.73	0.0097

**Table 3. Signal/Noise improvement for GluSPH.** The peak-to-peak S/N for GluSPH was measured at selected concentrations across the calibration curves for both the SCIEX 7500 system and the QTRAP 6500+ system. The peak-to-peak S/N improvement is between 2.4 and 6.5x.

Concentration (pmol/mL)	SCIEX 6500+ system - S/N (GluSPH)	SCIEX 7500 system - S/N (GluSPH)	S/N improvement
1.000	6233	21850	3.5x
0.020	119	284	2.4x
0.001	10	65	6.5x

**Table 4. Signal/Noise improvement for GalSPH.** The peak-to-peak S/N for GalSPH was measured at selected concentrations across the calibration curves for both the SCIEX 7500 system and the QTRAP 6500+ system. As can be seen the peak-to-peak S/N improvement is between 3.2 and 6.7x.

Concentration (pmol/mL)	SCIEX 6500+ system - S/N (GalSPH)	SCIEX 7500 system – S/N (GalSPH)	S/N improvement
1.000	6283	33324	5.3x
0.020	122	388	3.2x
0.001	6	40	6.7x

## Conclusions

- Sensitivity for two glycosphingolipids, galactosylsphingosine (GalSPH) and glucosylsphingosine (GluSPH) were evaluated in cerebral spinal fluid (CSF)
- Two MS systems were evaluated, the QTRAP 6500+ system was compared to the SCIEX 7500 system
- Improvements in S/N of between 2.4 and 6.7 x were achieved on the SCIEX 7500 system, as compared to the QTRAP 6500+ system
- Good accuracy and linearity were observed across the range evaluated on the SCIEX 7500 system.
- Improved S/N at similar concentrations of the standards in matrix suggests that lower volumes of CSF could be used in this assay on the SCIEX 7500 system, this hypothesis will be evaluated next.
- Based on the results, it is concluded that the objective to increase the sensitivity is met. Based on the improved S/N seen on the SCIEX 7500 system when compared to similar analysis on the SCIEX Triple Quad 6500+ System it may be possible to reduce the required assay volume of CSF for this method when using the SCIEX 7500 system.

## References

1. Rohini Sidhu et al. (2018). A HILIC-MS/MS Method for Simultaneous Quantification of the Lysosomal Disease Markers Galactosylsphingosine and Glucosylsphingosine in Mouse Serum. *Biomedical Chromatography*. **32(7)**. e4235.
2. Matthew Surface et al. (2021). Plasma Glucosylsphingosine in GBA1 Mutation Carriers with and without Parkinson's Disease. *Mov Disord*. **10**. 1-6.

The SCIEX clinical diagnostic portfolio is For In Vitro Diagnostic Use. Rx Only. Product(s) not available in all countries. For information on availability, please contact your local sales representative or refer to <https://sciex.com/diagnostics>. All other products are For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd. or their respective owners in the United States and/or certain other countries (see [www.sciex.com/trademarks](http://www.sciex.com/trademarks)).

© 2022 DH Tech. Dev. Pte. Ltd. RUO-MKT-02-14364-A.



**Headquarters**  
500 Old Connecticut Path | Framingham, MA 01701 USA  
Phone 508-383-7700  
[sciex.com](http://sciex.com)

**International Sales**  
For our office locations please call the division headquarters or refer to our website at [sciex.com/offices](http://sciex.com/offices)