

# Chemical Labeling Strategy and LC-MS/MS Analysis of Amine-containing Neurotoxin in Shark's Fins

See Chung Yip and Jason Neo  
AB SCIEX, Singapore

## Introduction


$\beta$ -Methylamino-L-alanine (BMAA), is a neurotoxin<sup>1</sup> and a non-proteinogenic amino acid. BMAA has been identified and quantified in cycad seeds and cyanobacteria by Rosen<sup>2</sup> et al. using liquid chromatography tandem mass spectrometry without pre-derivatisation. Kushnir et al.<sup>3</sup> has also developed a quantitative analysis of BMAA with iTRAQ<sup>®</sup> reagent as derivative in biological samples and plant extracts. No BMAA was detected in human serum samples and cerebrospinal samples. However, 50 mg/kg (ppm) of BMAA was found in the extract of the cycad seed.

Recently, a high concentration ranging from 144 to 1836 ng/mg of BMAA was discovered in shark fins by a group of University of Miami scientists.<sup>4</sup> In Mondo et al.'s analysis, BMAA was pre-labeled with a fluorescent tag 6-aminoquinolyl-N-hydroxysuccinimidyl carbamate (AQC) before injecting into LC-MS/MS.

The main challenges of determining BMAA are that the neurotoxin is highly polar and it provides low sensitivity for MRM when analyzed in the free form. In this paper, a simple labeling of BMAA using mTRAQ<sup>®</sup>  $\Delta 0$  reagent for detection with HPLC and AB SCIEX 4000 QTRAP<sup>®</sup> system is reported for shark's fin.

The mTRAQ<sup>®</sup>  $\Delta 0$  is an amine labeling reagent useful for performing relative quantitation experiments of targeted proteins, peptides and post-translational modifications by LC-MS/MS.

Here, mTRAQ<sup>®</sup> was used for labeling of BMAA prior LC-MS/MS detection using Multiple Reaction Monitoring (MRM).

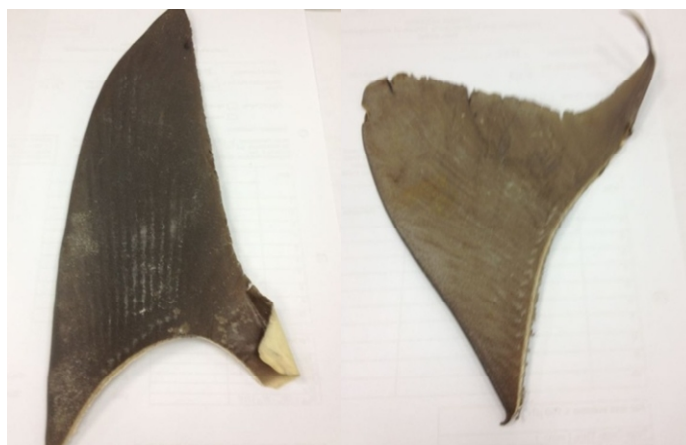

## Experimental

### Sample Preparation

The sample preparation protocol was adopted from the Mondo et al.<sup>4</sup> About 1 g of dried sample was weighed and digested with 6N HCl for 18 hours.

### Sample Labeling

5  $\mu$ L of the filtered sample was mixed with 15  $\mu$ L dissolution buffer (pH 8.5, 0.5 M triethylammonium bicarbonate). 50  $\mu$ L of isopropanol was added to the lyophilized mTRAQ<sup>®</sup> tube and vortexed. The mTRAQ<sup>®</sup> solution was then added to the 5  $\mu$ L sample and vortexed. The mixture was incubated at room temperature for 1 hour and then injected into LC-MS/MS.

### LC


A Shimadzu Prominence LC system with a Waters Atlantis HILIC column (100 x 2.1 mm, 3  $\mu$ m) at 40°C with a gradient of eluent A: 20 mM ammonium acetate + 0.1% formic acid in water and eluent B: 0.1% formic acid in acetonitrile was used at a flow rate of 300  $\mu$ L/min. The injection volume was set to 10  $\mu$ L.

## MS/MS

An AB SCIEX 4000 QTRAP<sup>®</sup> system with Turbo V<sup>™</sup> source and electrospray ionization (ESI) probe was used. 2 MRM transitions were monitored for BMAA; 399/228 was selected as quantifier and 399/172 was selected as qualifier transition.


## Results and Discussion

mTRAQ<sup>®</sup> reagent was used for labeling of BMAA. The proposed structure of the BMAA/mTRAQ<sup>®</sup> complex and the two monitored product ions are shown in Figure 1.

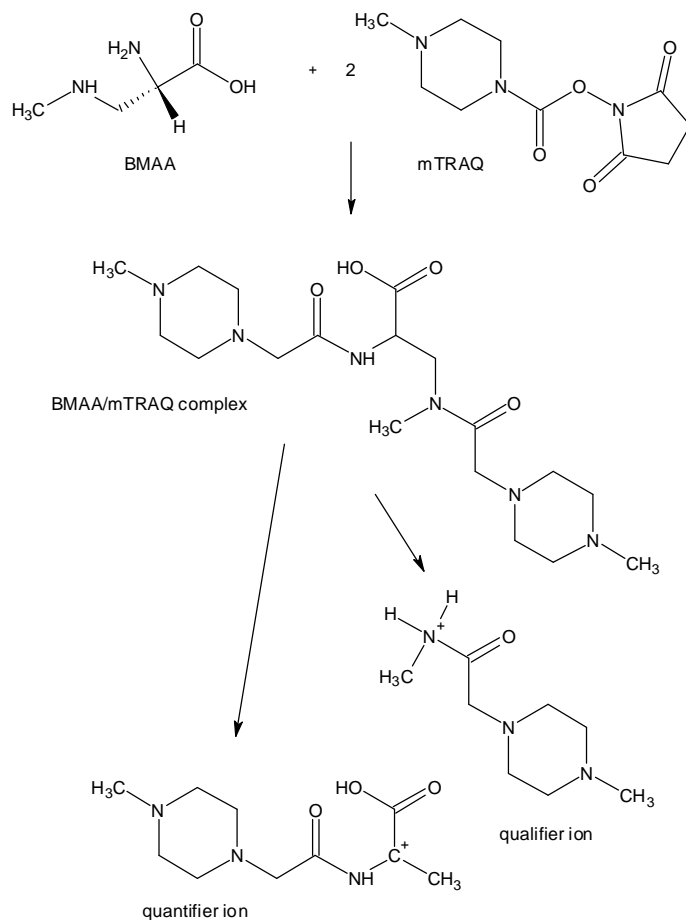


**Figure 1.** Proposed structure of the BMAA/mTRAQ<sup>®</sup> complex and the product ions monitored using MRM mode

An example chromatogram of a standard injection with a retention time of about 6.2 minutes is shown in Figure 2.

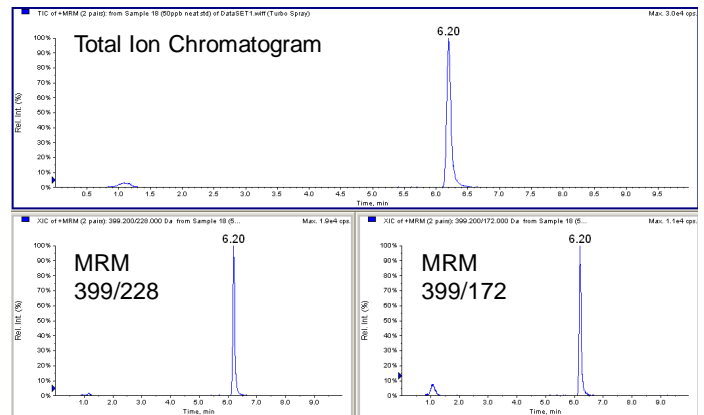

**Figure 2.** Detection of BMAA/mTRAQ<sup>®</sup> complex in neat standard of 50 ng/mL by LC-MS/MS

A 0.5, 1, 5, 10, 50, 100, 200, and 1000 mg/kg of BMAA was pre-spiked into the shark's fin, acid-digested, labeled with mTRAQ<sup>®</sup> and injected into the LC-MS/MS. A good linearity with r-value of greater 0.999 was observed (Figure 3). Repeatability was studied at 1, 5, 10, and 50 mg/kg over three consecutive days. The resulting coefficients of variation (%CV) are shown in Table 1.

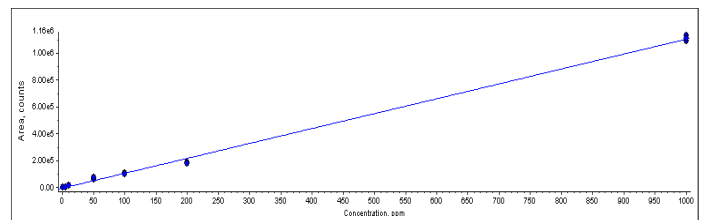


**Figure 3.** Calibration line from 0.5 to 1000 mg/kg BMAA spiked into shark's fin

**Table 1.** Repeatability of analysis on three consecutive days

%CV at	1 mg/kg	5 mg/kg	10 mg/kg	50 mg/kg
Day 1	19.7	14.5	3.2	9.5
Day 2	8.3	5.8	6.4	0.9
Day 3	10.9	7.8	6.4	1.2

The Limit of Quantitation (LOQ) was found to be 0.5 mg/kg BMAA in shark's fin with a signal to noise ratio of ~10 (Figure 4).

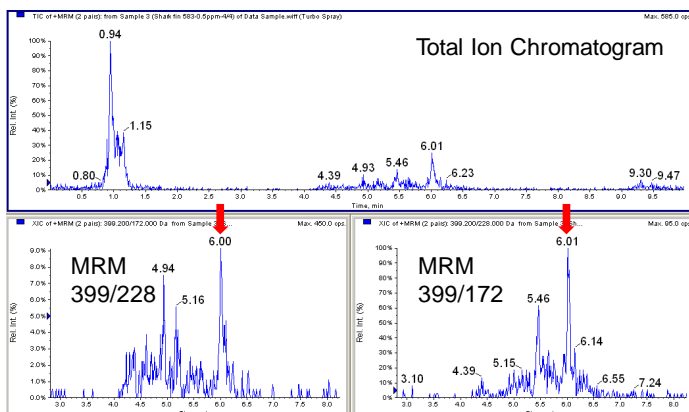

**Figure 4.** Detection of BMAA in 0.5 ppm spiked shark's fin after derivatization with mTRAQ<sup>®</sup> reagent

## Summary

A fast, simple and reliable labeling method with mTRAQ<sup>®</sup>  $\Delta$ 0 reagent, for the detection of BMAA in the shark's fin matrix, was developed for quantitative study on the AB SCIEX 4000 QTRAP<sup>®</sup> system. Multiple Reaction Monitoring (MRM) was used because of its high selectivity and sensitivity. The Limit of Quantitation (LOQ) of the method was found to be 0.5 mg/kg (ppm).

## Acknowledgements

The authors would like to thank the Veterinary Public Health Laboratory (VPHL) of the Agri-Food and Veterinary Authority of Singapore (AVA) for its collaboration in this study.

## References

- 1 <http://en.wikipedia.org/wiki/Beta-Methylamino-L-alanine>
- 2 Analyst 133 (2008) 1785-1789
- 3 Eur. J. Mass Spectrom 15 (2009) 439-443
- 4 Mar. Drugs 10 (2012) 509-520

**For Research Use Only. Not for use in diagnostic procedures.**

© 2012 AB SCIEX. The trademarks mentioned herein are the property of AB Sciex Pte. Ltd. or their respective owners. AB SCIEX™ is being used under license.

Publication number: 6640212-01

**Headquarters**  
500 Old Connecticut Path, Framingham, MA 01701 USA  
Phone 508-383-7700  
[www.sciex.com](http://www.sciex.com)

**International Sales**  
For our office locations please call the division  
headquarters or refer to our website at  
[www.absciex.com/offices](http://www.absciex.com/offices)